


Aug 15, 2018

Version 2

Sequential smFISH (Allen Institute) V.2

DOI

<https://dx.doi.org/10.17504/protocols.io.snnedde>

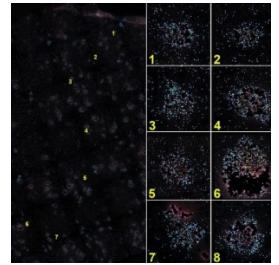

Jennie Close¹, Zoe Maltzer¹

¹Allen Institute for Brain Science

BICCN / BICAN



Stephanie Mok

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.snnedde>

Protocol Citation: Jennie Close, Zoe Maltzer 2018. Sequential smFISH (Allen Institute). **protocols.io**
<https://dx.doi.org/10.17504/protocols.io.snnedde>

License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: August 15, 2018

Last Modified: August 15, 2018

Protocol Integer ID: 14766

Keywords: smFISH, multiplexed single molecule fish protocol for use, multiplexed single molecule fish protocol, sequential smfish, mouse tissue, human tissue, protocol, nature protocol, tissue, fish

Abstract

We have developed a multiplexed single molecule FISH protocol for use at the Institute. This protocol was optimized on human tissue, but will work on mouse tissue as well. It was adapted from Lyubimova et. al., Nature Protocols, 2013.

Attachments



smFISH_Allen Institu...

16KB

Guidelines

Ensure that all reagents are in recombinant and RNase-free format, as we have noticed RNA degradation in solutions that contain enzymes derived from whole organisms.

We filter every solution with a 0.2um syringe filter prior to use. This reduces background spots and dust that interfere with imaging of diffraction limited spots.

For the SDS treatment after fixation and permeabilization, be gentle when dropping SDS onto the section, as well as during washes. This treatment is relatively harsh and the tissue must be treated somewhat delicately.

Materials

STEP MATERIALS

 4% PFA

 PBS

 PBS

 PBS

 2X SSC

 2X SSC

 65% formamide/2X SSC

 4% PFA

 PBS

 PBS

 PBS

 2X SSC

 2X SSC

 65% formamide/2X SSC

Protocol materials

 2X SSC

 PBS

 65% formamide/2X SSC

 4% PFA

Safety warnings

 Please refer to the SDS (Safety Data Sheet) for hazard information and safety warnings.
Avoid exposure to formamide, DAPI

Before start

Ensure all incubators and ovens are at the appropriate temperature prior to experiment.

Tissue and Sectioning

- 1 10-14 um cryosections are taken from fresh-frozen tissue, which are collected on poly-lysine-treated #1 coverslips at room temperature (RT). After 5-10 min at RT, sections are placed at 4°C until sectioning is complete. At that point, proceed immediately to fixation and permeabilization.

 00:05:00 RT

Fixation/Permeabilization

- 2 Post-fix sections for 15 min with 4% PFA @ 4 °C.

 4% PFA

 00:15:00 Post-fixing

 4 °C Post-fixing

- 3 Wash with PBS (1/3)

 PBS

- 4 Wash with PBS (2/3)

 PBS

- 5 Wash with PBS (3/3)

 PBS

- 6 Permeabilize with cold methanol at -20 °C for 10 min.

 -20 °C Permeabilizing

 00:10:00 Permeabilizing

- 7 Air dry for 30 min in fume hood (Stopping point: store coverslips at -80°C)

 00:30:00 Air drying

- 8 Optional: Treat sections with 8% SDS/PBS for 10 minutes, followed by 3 – 5 rinses with PBS or 2XSSC

 00:10:00

- 9 Add 2ml 2X SSC

 2X SSC

 2 µL 2X SSC



Hybridization

10 Pre-heat hyb oven to 37 °C

🔥 37 °C oven

11 Place sections in hyb buffer without probes.

12 Add 4 ul probe 400ul hyb buffer.

🧪 4 µL probe

🧪 400 µL hyb buffer

Note

Specific to 6-well plate format – if using perfusion chamber, this volume can be reduced.

13 Incubate at 37 C for 2H.

🔥 37 °C Incubation

🕒 02:00:00 Incubation

Wash

14 Add 2 ml wash buffer to each well.

🧪 2 µL wash buffer

15 Incubate at 37 C for 15 min.

🔥 37 °C Incubation

🕒 00:15:00 Incubation

16 Remove wash buffer.

17 Add 2 ml fresh wash buffer and incubate at 37 C for 15 min.

🧪 2 µL wash buffer

🔥 37 °C Incubation

🕒 00:15:00 Incubation

18 Replace wash buffer with fresh wash buffer + DAPI (final 5ug/mL) and incubate at 37 C for 15 min.



37 °C Incubation

00:15:00 Incubation

19 GLOX buffer step if performing antibody stain

20 Mount and image or store at 4 C in 2XSSC until imaging session

4 °C

2X SSC

Stripping

21 65% formamide/2X SSC, 10 min X 3, 30 C

65% formamide/2X SSC

00:10:00

30 °C

22 Wash in 2XSSC (1/3)

23 Wash in 2XSSC (2/3)

24 Wash in 2XSSC (3/3)

25 Following stripping, proceed to hybridization step.

[go to step #10](#) Hybridization step