

Jan 12, 2024

Short amplicons panels (Artic-like) for RSVA and RSVB

DOI

<https://dx.doi.org/10.17504/protocols.io.eq2lyj1nrlx9/v1>

Flora Donati¹, matthieu prot², Banujaa Jeyarajah¹, etienne.simon-loriere²

¹National Reference Center for viruses of respiratory infections, Institut Pasteur, Paris, France.;

²G5 Evolutionary Genomics of RNA Viruses, Institut Pasteur, Université Paris Cité, Paris, France.



etienne.simon-loriere

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.eq2lyj1nrlx9/v1>

Protocol Citation: Flora Donati, matthieu prot, Banujaa Jeyarajah, etienne.simon-loriere 2024. Short amplicons panels (Artic-like) for RSVA and RSVB. **protocols.io** <https://dx.doi.org/10.17504/protocols.io.eq2lyj1nrlx9/v1>

Manuscript citation:

Flora DONATI

Matthieu PROT

Banujaa JEYARAJAH

Etienne SIMON-LORIERE



License: This is an open access protocol distributed under the terms of the **[Creative Commons Attribution License](#)**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: November 07, 2023

Last Modified: January 12, 2024

Protocol Integer ID: 90547

Keywords: rsv sequencing, amplicon, panel, RSV, RSVA, RSVB, Respiratory syncytial virus, NGS, HTS, Illumina, ONT, Oxford Nanopore Technologies, viral genome, viral nucleic acid extract, rsva, generating cdna, short amplicons panel, rsvb, files for bioinformatic analysis, rsvb this sop, bioinformatic analysis, pcr, multiplex pcr, genome, cdna

Funders Acknowledgements:

Institut Pasteur

Abstract

This SOP describes the procedure for generating cDNA from respiratory syncytial virus A or B (RSVA or RSVB) viral nucleic acid extracts and subsequently producing ~400nt amplicons tiling the viral genome in a multiplex PCR.

The resulting products can be sequenced using short (Illumina) or long (Oxford Nanopore Technologies) reads approaches.

The panels and files for bioinformatic analysis are available at:

https://github.com/SimonLoriereLab/RSV_amplicons_panels

Guidelines

The panels have been optimized using RSV A or RSV B positive samples circulating in Europe in 2018-2022. Processing samples from previous years or from regions outside of Europe might require further adjustments.



Materials

Primers

Primer schemes for RSVA and RSVB have been designed and optimized to cover the diversity of viruses circulating in the Northern hemisphere (2018-2023). Processing samples from previous years or from regions outside of Europe might require further adjustments. See the file below for primer sequences and their ratio.

We work with typed samples, but it may be possible to combine the panels for A and B.

Multiple assays are available for typing:

(1) Todd A et al, J Virol Methods, 2021 - Rapid detection of human respiratory syncytial virus A and B by duplex real-time RT-PCR. doi:10.1016/j.jviromet.2021.114171

(2) Wang et al, J Virol Methods, 2019 - Duplex real-time RT-PCR assay for detection and subgroup-specific identification of human respiratory syncytial virus. doi: 10.1016/j.jviromet.2019.113676

Other panels, including a panel compatible with both RSVA and RSVB have been developed:

Dong et al, J Clin Virol, 2023 - A simplified, amplicon-based method for whole genome sequencing of human respiratory syncytial viruses. doi: 10.1016/j.jcv.2023.105423

Primers may be ordered from any oligonucleotide company. For primer preparation instructions see **Step 5** of the protocol.

Consumables

Filtered pipette tips

1.5 and 2mL microcentrifuge tubes

15 and 50 mL Falcon tubes

8-strip PCR tubes or PCR plates with caps or heat-sealing film

Reservoirs

Waste containers

Magnetic rack for PCR tubes or 96w plates

Molecular Biology

LunaScript RT SuperMix, M3010X, NEB - or equivalent

Q5 High Fidelity DNA polymerase, M0491L, NEB

dNTP Mix 10MM, 10319879, Thermo Fisher Scientific

Agencourt AMPure beads XP, A63880, Beckman Coulter

Nuclease Free Water, 10526945, Thermo Fisher Scientific

Qubit dsDNA HS Assay kit, Q32854, Thermo Fisher Scientific



Protocol materials

✕ Qubit® dsDNA HS Assay Kit **Thermo Fisher Scientific Catalog #Q32854**

✕ Agencourt AmPure XP beads **Catalog #A63880**

Safety warnings

! Wear appropriate personal protective equipment.

Before start

It is recommended to use to the forward flow principle to avoid contamination. All mastermixes should be made in a **mastermix cabinet** and aliquoted . Samples should be added in the **extraction and sample cabinet**. Amplicons should be purified in a **post-PCR cabinet**.

Mastermix, extraction and sample and **post-PCR cabinet** should be cleaned with appropriate disinfectant and UV-sterilized before and after use.



RNA preparation

1 Sample selection:

Viral RNA input from a clinical sample (recommended Ct value below 28).

It is recommended to dilute high viral load samples (below Ct 14-15) to reduce the likelihood of PCR inhibition.

2 Viral RNA extraction:

Viral RNA should be extracted with QIAamp Viral RNA extraction kit (Qiagen) or equivalent, following the manufacturer instructions.

Note

This step is also compatible with NucleoSpin 8x48 virus Core kit (Macherey-Nagel)

cDNA preparation

3 In a **mastermix cabinet**, mix the following components:

Component	Volume
LunaScript Master Mix	 2 µL
Template RNA	 8 µL

Mix by pipetting gently and spin down the tube.

Note

Always add a no-template RT-PCR control with nuclease-free water.

Most cDNA kits should work for this step (also tested with SuperScript III or IV (Invitrogen))

4 Incubate as follows:

 25 °C for  00:02:00

 55 °C for  00:20:00

23m





95 °C for 00:01:00

Hold at 4 °C

Then, place On ice .

Primer pool preparation

- 5 Resuspend lyophilized primers at a concentration of 100µM each.

Note

Primer schemes for RSVA and RSVB have been designed and optimized to cover the diversity of viruses circulating in the Northern hemisphere (2018-2023). They generate overlapping ~400bp amplicons.
See the file below for primer sequences and their ratio.

- 6 In a **mastermix cabinet**, generate working primer stocks by diluting primers at the adequate concentration in nuclease-free water (for example in a 96-well plate).
Note that some primers need to be added at a different concentration to help normalize sequencing coverage.

- 7 For each RSV subtype, prepare both primer pool stocks by adding 5 µL of each primer pair to a 1.5 mL Eppendorf tube labelled either RSV A Pool 1, RSV A Pool 2, RSV B Pool 1 or RSV B Pool 2.

Note

RSV A:

Pool 1 has **50** primers, Pool 2 has **50** primers.

By adding 5 µL , total volume should be 250 µL for Pool 1 and 250 µL for Pool 2.

RSV B:

Pool 1 has **52** primers, Pool 2 has **54** primers.

By adding 5 µL , total volume should be 260 µL for Pool 1 and 270 µL for Pool 2.

- 8 Dilute the primer pool stocks to 1:10 in nuclease-free water to generate working primer pool stocks.

Note

Make aliquots of both primer pools stocks in case of degradation or contamination. Store at  -20 °C .

Primers need to be used at a final concentration of  0.015 micromolar (μM) per primer.

Multiplex PCR

- 9 In a **mastermix cabinet**, set up the multiplex PCR reactions as follows:

>RSV A:

Component	Pool 1 or Pool 2
5X Q5 reaction buffer	 5 μL
10mM dNTP	 0.5 μL
Primer Pool 1 or 2 (working stock)	 1.88 μL
Nuclease-free Water	 14.9 μL
Q5 Polymerase	 0.25 μL
Total	 22.5 μL

>RSV B:

Component	Pool 1	Pool 2
5X Q5 reaction buffer	 5 μL	 5 μL
10mM dNTP	 0.5 μL	 0.5 μL
Primer Pool 1 or 2 (working stock)	 1.95 μL	 2.02 μL
Nuclease Free Water	 14.73 μL	 14.8 μL
Q5 Polymerase	 0.25 μL	 0.25 μL
Total	 22.5 μL	 22.5 μL



It is recommended to prepare a mastermix corresponding to the number of samples to be processed, always including a no template control per reaction.

Note: This highly multiplex reaction has been designed and tested with the NEB Q5 enzyme. Other types of enzymes may not be compatible.

- 10 In the **extraction and sample addition cabinet**, add  2.5 µL of cDNA to each tube and mix well by pipetting.

Note

Add a negative RT-PCR control with nuclease free water.

- 11 Spin down the tubes/strips/plates.

- 12 Set-up the following program on a thermal cycler:

5m 45s



Steps	Temperature	Time	Cycles
Heat Activation	 98 °C	 00:00:30	1
Denaturation	 98 °C	 00:00:15	35
Annealing	 65 °C	 00:05:00	35
Hold	 16 °C	infinite	1

PCR clean-up

- 13 *The following steps should be performed in a **post-PCR cabinet**.*

Combine the entire contents of Pool 1 and Pool 2 PCR reactions for each biological sample into a single tube.



14 Run  5 μL of each pooled sample on a 1% agarose gel.



Expected result

Expected amplicon size is ~400bp

15 Clean-up the amplicons using the following protocol.

15.1 **Vortex** SPRI beads (at room temperature) thoroughly. The solution should be homogenous if well resuspended.



 Agencourt AmPure XP beads **Catalog #A63880**

15.2 Add 1,8X of SPRI beads ( 45 μL if no PCR product was run on a gel) to the sample tube and mix gently by pipetting.

15.3 Spin down the tubes to collect all liquid at the bottom of the tube.

15.4 Incubate  00:05:00 to  00:10:00 at  Room temperature

15m

15.5 Place the tubes on an appropriate magnetic rack and incubate for  00:05:00 to  00:10:00 until the beads have pelleted and the supernatant is clear.

15m

15.6 Carefully remove and discard the supernatant, being careful not to touch the bead pellet.

15.7 Add  200 μL of **fresh**,  Room temperature 80% ethanol to the pellet while in the magnetic rack, and incubate for  00:01:00 .

1m

15.8 Carefully remove and discard ethanol, being careful not to touch the beads pellet.



- 15.9  [go to step #15.7](#) and repeat ethanol wash.
- 15.10 Spin down the tubes to collect all liquid at the bottom of the tube and carefully remove as much residual ethanol as possible using a P10 pipette. Do not touch the beads.
- 15.11 Incubate the tubes for  00:05:00 at  Room temperature with the lid open (drying step). 5m
- 15.12 Add  20 μL of nuclease-free water, remove the tubes from the magnetic rack and resuspend the pellet by gently pipetting. Incubate for  00:02:00 . 2m
- 15.13 Place the tubes back on the magnetic rack until the solution is clear and transfer the supernatant into clean tubes (e.g. 0.2mL 8-strip PCR). Make sure not to transfer any beads.

Quantification and normalisation

- 16 Quantify the DNA concentration using the Qubit High Sensitivity DNA kit (or equivalent) from 1 μL of each product.

 Qubit® dsDNA HS Assay Kit Thermo Fisher Scientific Catalog #Q32854

Expected result

Expected range = 10-100 ng/ μL DNA.

Library preparation

- 17 The PCR products can be used to prepare libraries for short or long reads sequencing.