



May 30, 2024

Version 2

Simultaneous detection of miRNA and mRNA at the single-cell level in plant tissues (v2) V.2

DOI

<https://dx.doi.org/10.17504/protocols.io.q26g714e8gwz/v2>

Chi-Chih Wu¹

¹Biodiversity Research Center, Academia Sinica, Taipei 11529, Taiwan



Chi-Chih Wu

IPMB

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.q26g714e8gwz/v2>

Protocol Citation: Chi-Chih Wu 2024. Simultaneous detection of miRNA and mRNA at the single-cell level in plant tissues (v2). protocols.io <https://dx.doi.org/10.17504/protocols.io.q26g714e8gwz/v2> Version created by [Chi-Chih Wu](#)

License: This is an open access protocol distributed under the terms of the [Creative Commons Attribution License](#), which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working



Created: May 30, 2024

Last Modified: May 30, 2024

Protocol Integer ID: 100911

Keywords: mirna molecules in specific cell, mirna molecule, transgenic rice lines with different gene expression level, mirna, transgenic rice line, cell level in plant tissue, plant tissue, lna probe, mrna, expression levels of mrna, gene expression, different gene expression level, single cell, locked nucleic acid, detection signal

Abstract

Detecting the simultaneous presence of a microRNA (miRNA) and a mRNA in a specific tissue can provide support for the prediction that the miRNA regulates the mRNA. We develop a method that uses sequence-specific miRNA-locked nucleic acid (LNA) and mRNA-LNA probes. Moreover, it augments the detection signal by rolling circle amplification, achieving a high signal-noise ratio at the single-cell level. Dot signals are counted for determining the expression levels of mRNA and miRNA molecules in specific cells. We show a high sequence specificity of our miRNA-LNA probe, revealing that it can discriminate single-base mismatches. Numerical quantification by our method is tested in transgenic rice lines with different gene expression levels.



1 Section permeabilization

	A
	1. The slides with sections are taken out the freezer and equilibrated to RT for 40 mins.
	2. Permeabilized in 20 ug/ml proteinase k for proper duration
	3. Quickly wash in DEPC-PBS
	4. Quickly dehydrate the slides in EtOH (50, 70, 99 %) and then air dry
	5. Mount the secure seal reaction chambers onto the slides.
	6. Add 1x DEPC-PBS-tween 0.05 % (Wash buffer) into the chambers to keep the slides wet until RT reaction ready

2 Mixture for miRNA hybridization (50 ul)

	A	B	C
		Stock	Final
	Formamide	100%	50%
	SSC	20x%	5x
	tRNA	10 mg/ml	0.5 ug/ul
	Denhardt's	50x	1x
	LNA probe A	10 uM	2-3 pmole
	DEPC-H ₂ O	-	

1. process

	A
	Hybridization below the predicted melting temperature of the probe, about 2 C, for an hour
	Wash with 0.1X SSC three times at the temperature set in the Step 1



	A
	Wash with 2X SSC at RT once
	Wahs wtih the wash buffer (PBS, 0.05% Tween-20) once

3 Mixture for mRNA cDNA synthesis (50 ul)

	Reagent	Stock	Final
	NEB Tag DNA ligase	40U/ul	0.5 U/ul
	Rnase H	5 U/ul	0.4 U/ul
	Ribolock Rnase inhibitor	40 U/ul	1 U/ul
	NEB Tag ligase buffer	10x	1x
	BSA	20 ug/ul	0.2 ug/ul
	KCl	1 M	0.05 M
	Formamide	100%	20%
	Pd_A	10 uM	0.1 uM
	Pd_B	10 uM	0.1 uM
	Pd_C	10 uM	0.1 uM
	DEPC-H2O		

Process

	1. Add ligation mixture in chambers, seal with adhesive film
	2. Incubate for 30 min at 37 C followed by 45 min at 48 C

	3. Wash 2x, 1x DEPC-PBS-Tween 20, 0.05%

4 Mixture for miRNA and mRNA padlock probe hybridization and ligation (50 ul)

	Reagent	Stock	Final
	NEB Tag DNA ligase	40U/ul	0.5 U/ul
	Rnase H	5 U/ul	0.4 U/ul
	Ribolock Rnase inhibitor	40 U/ul	1 U/ul
	NEB Tag ligase buffer	10x	1x
	BSA	20 ug/ul	0.2 ug/ul
	KCl	1 M	0.05 M
	Formamide	100%	20%
	Pd_A	10 uM	0.1 uM
	Pd_B	10 uM	0.1 uM
	Pd_C	10 uM	0.1 uM
	DEPC-H2O		

process

	A
	1. Add ligation mixture in chambers, seal with adhesive film
	2. Incubate for 30 min at 37 C followed by 45 min at 48 C
	3. Wash 2x, 1x DEPC-PBS-Tween 20, 0.05%



5 Rolling circle amplification (50 ul)

	Reagent	Stock	Final
	Phi 29 polymerase	10 U/ul	1 U/ul
	Ribolock Rnase inhibitor	40 U/ul	1 U/ul
	10 X phi 29 buffer	10 x	1 x
	dNTP	10 mM	0.25 mM
	BSA	20 ug/ul	0.2 ug/ul
	Glycerol	50%	5%
	DEPC-H ₂ O		

process

	A
	1. Add reaction mixture and seal chamber
	2. Incubate for over night at 37 C
	3. Wash 2x, DEPC-PBS-Tween 20, 0.05 %

6 Mixture for detection oligo hybridization (50 ul)

	A	B	C
	Reagent	Stock	Final
	Hyb mixture	4 x	2 x
	Detection oligo 1-FITC	1 uM	0.1 uM



	A	B	C
	Detection oligo 2-Cy3	1 uM	0.1 uM
	Detection oligo 3-Cy5	1 uM	0.1 uM
	DEPC-H2O		

process

	1. Add reaction mixture
	2. Incubate for 30 min at 37 C
	3. Wash 2x, DEPC-PBS-Tween 20, 0.05 %
	4. Dehydrated by EtOH 50, 70, 99 %; then air dry.
	5. Mount cover slips