

Oct 18, 2020

Single digested

DOI

dx.doi.org/10.17504/protocols.io.bni5mcg6

Jiaxin Li¹

¹South China University of Technology



Jiaxin Li

OPEN  ACCESS



DOI: dx.doi.org/10.17504/protocols.io.bni5mcg6

Protocol Citation: Jiaxin Li 2020. Single digested. **protocols.io** <https://dx.doi.org/10.17504/protocols.io.bni5mcg6>

License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: October 18, 2020

Last Modified: October 18, 2020

Protocol Integer ID: 43325



- 1 Add the following reagents to a tube.

Total	10ul
10 x Buffer	1ul
Template (500ng)	According to different template
Enzyme	0.5ul
ddWater	Add to 10ul

- 2 Pipette up and down thoroughly.
- 3 Put at 37°C for about 1.5 hours (according to different enzymes)