



Nov 05, 2023

Sleep-Wake Recording in Dan lab

DOI

<https://dx.doi.org/10.17504/protocols.io.x54v9pkd4g3e/v1>

Xiaolin Huang¹

¹University of California, Berkeley



Xiaolin Huang

The University of California

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.x54v9pkd4g3e/v1>

Protocol Citation: Xiaolin Huang 2023. Sleep-Wake Recording in Dan lab. **protocols.io**

<https://dx.doi.org/10.17504/protocols.io.x54v9pkd4g3e/v1>

License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: October 27, 2023



Last Modified: November 05, 2023

Protocol Integer ID: 89994

Keywords: wake recording in dan lab surgery, wake cycles in mice, wake recording, hour natural sleep, wake cycle, sleep, recording, dan lab surgery, mice, setups for recording

Funders Acknowledgements:

ASAP

Grant ID: 020551

Abstract

Surgery and Setups for Recording 24-Hour Natural Sleep/Wake Cycles in Mice.

Surgery

- 1 Mice were anesthetized with 1.5% isoflurane and placed in a stereotaxic frame.
- 2 Body temperature was maintained using a heating pad.
- 3 The skin was incised to expose the skull after asepsis and connective tissue was removed.
- 4 To implant electroencephalogram (EEG) and electromyogram (EMG) recording electrodes, two stainless steel screws were inserted into the skull 2.5 mm from the midline and 3 mm posterior to the bregma, and two EMG electrodes were also inserted into the neck muscles.
- 5 A reference screw for grounding was placed on top of the left cerebellum.
- 6 Insulated leads from the EEG and EMG electrodes were soldered to a pin header, which was secured to the skull using dental cement.

Recording

- 7 After a minimum of 7 days of undisturbed post-surgery recovery, behavioral experiments were carried out in home cages placed in sound-attenuating boxes.
- 8 EEG and EMG electrodes were connected to flexible recording cables via a mini-connector.
- 9 Recordings started after 1-2 days of habituation and continued for 3-4 days on a 12-h dark/12-h light cycle.
- 10 Mouse behavior was recorded using a video camera at 10 frames per second (fps), along with EEG and EMG recordings.



- 11 The signals were recorded with a TDT RZ10x/PZ5-32 or LR10-SK1 system.