

Feb 05, 2020 Version 1

Slide Preparation for Mounting Samples for TissueCyte 1000 Sectioning and Imaging V.1

DOI

dx.doi.org/10.17504/protocols.io.bbx3ipqn

Allen Institute for Brain Science¹

¹Allen Institute

BICCN / BICAN

Allen Institute for Brain S...



Dillan Brown

OPEN  ACCESS



DOI: dx.doi.org/10.17504/protocols.io.bbx3ipqn

Protocol Citation: Allen Institute for Brain Science 2020. Slide Preparation for Mounting Samples for TissueCyte 1000 Sectioning and Imaging. **protocols.io** <https://dx.doi.org/10.17504/protocols.io.bbx3ipqn>

License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: January 29, 2020

Last Modified: February 05, 2020

Protocol Integer ID: 32475

Keywords: TissueCyte, Sectioning, Mounting, Imaging,

Abstract

This protocol describes the preparation of 1×3 glass slides for mounting embedded mouse brain samples for subsequent imaging and sectioning on TissueCyte 1000.

Attachments



PF0260_TissueCyte_sl...

160KB

