

Oct 17, 2023

🌐 Soil extract for cocultures host-parasites media



Forked from [Soil extract for algal media](#)

DOI

<https://dx.doi.org/10.17504/protocols.io.rm7vzxkq2gx1/v1>



Roscoff Culture Collection¹, Estelle Bigeard², Laure Guillou²

¹Station Biologique, CNRS-Sorbonne Université, Roscoff France; ²Station Biologique de Roscoff, France

Ecology of Marine Plank...

Roscoff Culture Collection

1 more workspace



Estelle Bigeard

Station Biologique de Roscoff, UMR7144 CNRS Sorbonne Univers...

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.rm7vzxkq2gx1/v1>

Protocol Citation: Roscoff Culture Collection, Estelle Bigeard, Laure Guillou 2023. Soil extract for cocultures host-parasites media. [protocols.io https://dx.doi.org/10.17504/protocols.io.rm7vzxkq2gx1/v1](https://dx.doi.org/10.17504/protocols.io.rm7vzxkq2gx1/v1)

License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: September 18, 2023

Last Modified: October 17, 2023

Protocol Integer ID: 87926

Keywords: parasites media, soil, cocultures host

Funders Acknowledgements:

ANR AQUARADOX, PARALEX, EPHEMER

Grant ID: SU & CNRS

Abstract

How to prepare soil extract

Materials

Equipments

Autoclave

Laminar flow cabinet (or biosafety cabinet)

Vacuum pump

Magnetic hot plate stirrers

Materials

Glass fibre prefilters (Millipore, AP1507500)

Filtration unit Stericup GP Merck Millipore, pore size 0.22 µm (Millipore, S2GPU02RE)

Before start

Please refer to our general recommendations to grow cultures :

<https://www.protocols.io/private/A48906DC1374AD6281495CB86A8F092F>



Soil sampling

- 1 Soil is sampled in a pine wood, located in Brittany, France.

To find a correct sampling area in your country, try to find a location devoid of pesticides and pollutants. It should be also a place where you know that you can collect the soil again and later in order to always use the same soil, especially if this one worked.

Collect the soil in plastic boxes.

Back to the laboratory, plastic boxes are placed at 37°C for 12 to 24 hours to let the soil dry. The soil must be dry before use, otherwise the boiling step should be longer.

Preparation of the soil extract solution (25g/l)

- 2 Weigh 30g of dried soil and place it in a 4L glass erlenmeyer.

- 3 Add 1200ml of MilliQ water to the soil.
Close the erlenmeyer with a aluminium foil.

- 4 Heat at 100°C with pemanent mixing during 1 hour to obtain an efficient boiling.
Adjust this time depending on the volume (more volume, more time).
Let cool.

- 5 Filtrate on 0.2 µm, using a glass prefilter and a pump vacuum.
As the prefilter will be rapidly clogged, better to add the boiled soil 100mL per 100mL.
If this happens, replace the glass prefilter, then the filter.
Transfer the solution in one or several glass bottles.

- 6 Autoclave at 121°C for 20 minutes.
Place the bottle of the filtered-boiled soil under a safety cabinet, and let it cool.

- 7 The day after, when the medium is at room temperature, place it at 37°C for 24 to 48 hours.
Then autoclave again at 121°C for 20 minutes.

This step allows to reactivate spores (fungi etc) contained in soil which have not been destroyed by the first sterilization, and to kill them during the second one.

- 8 Add 50ml of soil extract per 1 L of medium.

