

Jul 23, 2020 Version 2

SOLUTION- 02 - Phosphate Buffered Saline (PBS) V.2

DOI

dx.doi.org/10.17504/protocols.io.biahkab6

Marco Cosentino¹, Elisa Storelli¹, Alessandra Luini¹, Massimiliano LM Legnaro¹, Emanuela Rasini¹, Marco Ferrari¹, Franca Marino¹

¹Center for Research in Medical Pharmacology, University of Insubria (Varese, Italy)



Elisa Storelli

Center for Research in Medical Pharmacology, University of I...

OPEN  ACCESS



DOI: dx.doi.org/10.17504/protocols.io.biahkab6

Document Citation: Marco Cosentino, Elisa Storelli, Alessandra Luini, Massimiliano LM Legnaro, Emanuela Rasini, Marco Ferrari, Franca Marino 2020. SOLUTION- 02 - Phosphate Buffered Saline (PBS). **protocols.io**

<https://dx.doi.org/10.17504/protocols.io.biahkab6>

License: This is an open access document distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Created: July 07, 2020

Last Modified: July 27, 2020

Document Integer ID: 38953

Abstract

This recipe is used in the following protocols:

- Separation and purification of human PBMC from FRESH BLOOD
- Separation and purification of human PBMC from BUFFY COAT
- Magnetic bead-based CD4+ T cell isolation from PBMCs with Dynabeads: CD4 Positive Isolation Kit
- Magnetic bead-based TREG-TEFF cell isolation from PBMC with Miltenyi CD4+CD25+ Regulatory T cell Isolation Kit
- Staining of human PBMC or ISOLATED SUBSETS with Cell Proliferation Dye-eFluor™ 670 (CPD-eFluor670) for cell proliferation evaluation by Flow Cytometry

Phosphate-buffered saline (PBS 1X)

NaCl - 8 g (137 mM)

KCl - 0.2 g (2.7 mM)

Na₂HPO₄ - 1.44 g (10 mM)

KH₂PO₄ - 0.24 g (1.8 mM)

To prepare 1 L of PBS 1X, dissolve the reagents listed above in 800 mL of ultrapure H₂O.

Adjust the pH to 7.4 with HCl or NaOH, and then add H₂O to 1 L.

NaCl code: S9625, Sigma

KCl code: P9541, Sigma

Na₂HPO₄ code: 1.06585, Sigma

KH₂PO₄ code: P0662, Sigma

Fonte: (Cold Spring Harbor Laboratory Press 2006)