

Jul 23, 2020

Version 2

SOLUTION- 04 - Wash solution (RPMI/FBS) for PBMC V.2

DOI

dx.doi.org/10.17504/protocols.io.biaikace

Marco Cosentino¹, Elisa Storelli¹, Alessandra Luini¹, Massimiliano LM Legnaro¹, Emanuela Rasini¹, Marco Ferrari¹, Franca Marino¹

¹Center for Research in Medical Pharmacology, University of Insubria (Varese, Italy)



Farmacologia Medica

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.biaikace>

Document Citation: Marco Cosentino, Elisa Storelli, Alessandra Luini, Massimiliano LM Legnaro, Emanuela Rasini, Marco Ferrari, Franca Marino 2020. SOLUTION- 04 - Wash solution (RPMI/FBS) for PBMC. **protocols.io**

<https://dx.doi.org/10.17504/protocols.io.biaikace>

License: This is an open access document distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Created: July 07, 2020



Last Modified: November 21, 2020

Document Integer ID: 38954

Keywords: pbmc, wash solution, rpmi, fb

Troubleshooting



Sterile RPMI/FBS 10% (v/v)

Prepare with sterile RPMI and FBS (Fetal Bovine Serum). Both can be found in Fridge 2 - (Room PSO3).

Prepare 50 ml of RPMI/FBS by adding 5 ml of FBS to 45 ml of RPMI, work under a laminar flow hood (Room PSO3).

Storage:  4 °C Fridge 2 - (Room PSO3).

RPMI: catalog number ECM0495L, Euroclone, Italy.

FBS: catalog number ECS0180L, Euroclone, Italy.