



Aug 03, 2020

SOLUTION- 10 - TEFF/TREG isolation buffer

DOI

dx.doi.org/10.17504/protocols.io.bjadkia6

Marco Cosentino¹, Elisa Storelli¹, Alessandra Luini¹, Massimiliano LM Legnaro¹, Emanuela Rasini¹, Marco Ferrari¹, Franca Marino¹

¹Center for Research in Medical Pharmacology, University of Insubria (Varese, Italy)



Farmacologia Medica

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.bjadkia6>

Document Citation: Marco Cosentino, Elisa Storelli, Alessandra Luini, Massimiliano LM Legnaro, Emanuela Rasini, Marco Ferrari, Franca Marino 2020. SOLUTION- 10 - TEFF/TREG isolation buffer. **protocols.io**

<https://dx.doi.org/10.17504/protocols.io.bjadkia6>

License: This is an open access document distributed under the terms of the **[Creative Commons Attribution License](#)**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Created: August 03, 2020

Last Modified: August 03, 2020

Document Integer ID: 39973

Keywords: isolation, treg, teff

Troubleshooting



This buffer is composed of PBS, 0.5% BSA and 2mM EDTA.

It must be **cold (2-8°C) and sterile filtered before use.**

BSA 1.250 g,

EDTA 0.2235 g

Dissolve the following powders in **240 mL PBS 1X**, adjust the pH to 7.2 with HCl or NaOH, and then add PBS 1X to 250 mL.

Storage: 🌡 4 °C (Fridge - 2 Room PS03)

BSA: catalog number A2153 , Sigma, Italy

EDTA: catalog number ED2SS, Sigma, Italy