



Nov 21, 2020

SOLUTION- 13 - Complete culture medium

DOI

<https://dx.doi.org/10.17504/protocols.io.bpxkmpkw>

Marco Cosentino¹, Elisa Storelli¹, Alessandra Luini¹, Massimiliano Legnaro¹, Emanuela Rasini¹, Marco Ferrari¹, Franca Marino¹

¹Center for Research in Medical Pharmacology, University of Insubria (Varese)



Farmacologia Medica

University of Insubria

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.bpxkmpkw>

Document Citation: Marco Cosentino, Elisa Storelli, Alessandra Luini, Massimiliano Legnaro, Emanuela Rasini, Marco Ferrari, Franca Marino 2020. SOLUTION- 13 - Complete culture medium. **protocols.io**

<https://dx.doi.org/10.17504/protocols.io.bpxkmpkw>

License: This is an open access document distributed under the terms of the [Creative Commons Attribution License](#), which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Created: November 21, 2020

Last Modified: November 21, 2020

Document Integer ID: 44748

Keywords: solution, culture



Prepare 50 mL culture medium (**RPMI/FBS 10%/ Pen-Strep 100 U/mL**) with:

44,5 mL **RPMI 1640** with phenol-red supplemented with 5 mL **of FBS** (Fetal Bovine Serum) and 0,5 mL **penicillin/streptomycin** (equivament of 100 U/mL).

Alle reagents can be found in Fridge 2 - (Room PSO3). Work under a laminar flow hood (Room PSO3).

Storage:  4 °C Fridge 2 - (Room PSO3).

RPMI: catalog number ECM0495L, Euroclone, Italy.

FBS: catalog number ECS0180L, Euroclone, Italy.

Penicillin/Streptomycin: Catalog number ECB3001D - 100 ml, Euroclone, Italy.