


Feb 24, 2020

Version 2

SOP Appendix for Spleen V.2

DOI

<https://dx.doi.org/10.17504/protocols.io.bcwhixb6>

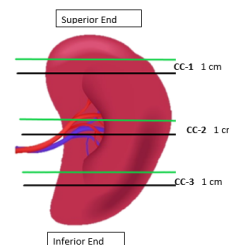

Francesca Farris¹, Marda Jorgensen¹

¹University of Florida

Human BioMolecular Atlas Program (HuBMAP) Method Development Community

Tech. support email: Jeff.spraggins@vanderbilt.edu



Marda Jorgensen

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.bcwhixb6>

Document Citation: Francesca Farris, Marda Jorgensen 2020. SOP Appendix for Spleen. **protocols.io**

<https://dx.doi.org/10.17504/protocols.io.bcwhixb6>

License: This is an open access document distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Created: February 24, 2020

Last Modified: October 03, 2020

Document Integer ID: 33449

Keywords: appendix for spleen sop, sop appendix for spleen, sop appendix, spleen sop, spleen, appendix, sop

Abstract


This is an appendix for spleen SOP.

Attachments



Appendix for SOP.doc...

115KB

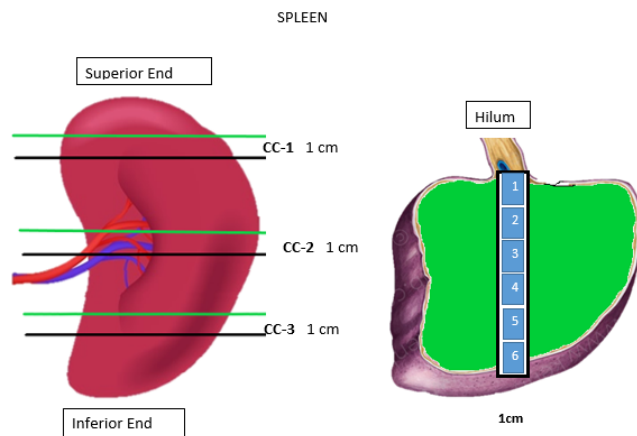

Distance between CC1 and CC2 _____ mm CC2 and CC3 _____ mm

	Block 1 (A)	Block 2 (B)	Block 3 (C)	Block 4 (D)	Block 5 (E)	Block 6 (F)
CC-1	upper (NBF-IMC)	upper (Fresh OCT)	10 X fresh tissue in medium	CLARITY PFA (megacassette)	upper (PFA)	
	lower (NBF)	lower (NBF)			lower (PFA OCT)	
CC-2	10 X fresh tissue in medium	upper (NBF-IMC)	upper (Fresh OCT)	upper CMC frozen	upper (PFA)	CLARITY PFA (megacassette)
		lower (NBF)	lower (NBF)	lower CMC frozen	lower (PFA OCT)	
CC-3	upper (NBF-IMC)	upper (Fresh OCT)	10 X fresh tissue in medium	CLARITY PFA (megacassette)	upper (PFA)	
	lower (NBF)	lower (NBF)			lower (PFA OCT)	

Photo document each step***Cut CCF regions. Orient slabs. Cut strips. Cut 1cm blocks. Bisect blocks as denoted for cassettes. Maintain up facing surface in cassettes. Cassettes will be embedded such that up surface is the sectioning surface.