

Feb 27, 2024

SOP for Flow cytometry after DSS-induced gut and brain injury

DOI

<https://dx.doi.org/10.17504/protocols.io.bp2l6x6zklqe/v1>

Malu Tansey¹

¹College of Medicine |University of Florida



Senthilkumar Karuppagounder

Duke University

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.bp2l6x6zklqe/v1>

Protocol Citation: Malu Tansey 2024. SOP for Flow cytometry after DSS-induced gut and brain injury. **protocols.io**
<https://dx.doi.org/10.17504/protocols.io.bp2l6x6zklqe/v1>

License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: January 28, 2024



Last Modified: May 31, 2024

Protocol Integer ID: 94272

Keywords: ASAPCRN, brain injury sop for flow cytometry, sop for flow cytometry, brain injury sop, flow cytometry, induced gut

Funders Acknowledgements:

Aligning Science Across Parkinson's (ASAP)

Grant ID: ASAP-020527

Abstract

SOP for Flow cytometry after DSS-induced gut and brain injury

Materials

LIVE/DEAD Fixable Blue stain (ThermoFisher, #L34962)

1% paraformaldehyde in a phosphate buffer

negative AbC Total Antibody Compensation Beads (ThermoFisher, #A10497)

Amine Reactive Compensation Beads (ThermoFisher, #A10346)

SPHERO Supra Rainbow particles (Spherotech, Inc., #SRCP-35-2A)



- 1 Following isolation of brain and circulating immune cells with magnetic beads, cells were transferred to a V-bottom plate for surface epitope labeling.
- 2 Samples were centrifuged at 300g for 5min at 4°C, washed once in PBS, re-pelleted, and resuspended in an antibody cocktail containing fluorophore-conjugated antibodies and LIVE/DEAD Fixable Blue stain (ThermoFisher, #L34962) in BD Horizon Brilliant Stain Buffer (BD Biosciences, #563794). 5m
- 3 Cells were incubated with antibodies for 20min at 4°C in the dark before three washes in FACS buffer containing 0.25mM EDTA, 0.01% NaN₃, and 0.1% BSA in PBS. 20m
- 4 Labeled cells were fixed in 1% paraformaldehyde in a phosphate buffer for 30min at 4°C in the dark. 30m
- 5 Cells were washed three times in FACS buffer and transferred to FACS tubes in FACS buffer for flow cytometry.
- 6 For compensation, 1uL of antibody was added to reactive and negative AbC Total Antibody Compensation Beads (ThermoFisher, #A10497) and 0.5uL of LIVE/DEAD Fixable Blue stain was added to ArC Amine Reactive Compensation Beads (ThermoFisher, #A10346).
- 7 Bead staining occurred at the same time as cell staining and was subjected to the same protocol described above without fixation (e.g., labeled beads remained in FACS buffer while cells were in fixative).
- 8 Samples were run on a BD FACSymphony A3, capturing either 100,000 single cells or the entire sample. Laser voltages were set with the assistance of SPHERO Supra Rainbow particles (Spherotech, Inc., #SRCP-35-2A) based on settings optimized during the development of this antibody panel. Compensation controls were calculated to ensure successful channel compensation prior to running samples. Raw files were analyzed using FlowJo (BD Biosciences) v.10.9.0 and gates were drawn with the assistance of fluorescence-minus-one controls.