

Jan 31, 2025

Version 1

SP3 protocol optimised for foodcrust protein extraction from archaeological cooking vessels V.1

 [PLOS One](#)

DOI

<https://dx.doi.org/10.17504/protocols.io.e6nvwb22wvmk/v1>

Virginia Harvey^{1,2}, Jessica Hendy²

¹University of Chester; ²University of York



Joannes Dekker

University of Turin

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

[Create free account](#)

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.e6nvwb22wvmk/v1>

External link: <https://doi.org/10.1371/journal.pone.0330195>

Protocol Citation: Virginia Harvey, Jessica Hendy 2025. SP3 protocol optimised for foodcrust protein extraction from archaeological cooking vessels. **protocols.io** <https://dx.doi.org/10.17504/protocols.io.e6nvwb22wvmk/v1>

Manuscript citation:

Dekker JAA, Hagan R, Collins M, Hendy J (2025) An unclean slate, discrepancies between food input and recovered protein signal from experimental foodcrusts. PLOS One 20(8). doi: [10.1371/journal.pone.0330195](https://doi.org/10.1371/journal.pone.0330195)

License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: January 24, 2025

Last Modified: January 31, 2025

Protocol Integer ID: 119014

Keywords: SP3, Palaeoproteomics, Organic residue analysis, Proteins, Pottery, Foodcrust, Archaeology, foodcrust protein extraction from archaeological cooking vessel, foodcrust protein extraction, archaeological cooking vessel, foodcrust, charred organic residue, extraction of protein, extraction, sp3 protocol, cooking vessel, sp3, protein, downstream contamination, contamination

Funders Acknowledgements:

Philip Leverhulme Prize

Grant ID: -

Abstract

A SP3-based protocol developed for the extraction of proteins from charred organic residues (foodcrust) found on cooking vessels. This protocol is to prepare samples for MALDI-ToF MS and LC-MS/MS analysis. We recommend starting with a batch size of 2-10 samples, and including flanking negative controls (blanks) to test for any downstream contamination.

Guidelines

Perform this protocol with the appropriate PPE and ideally in a clean lab setting.

Materials

Equipment

- Scale for laboratory use
- Magnetic tube rack
- Heat block/Thermoshaker
- Centrifuge
- Vortexer

Consumables

- Protein Low-Bind 1.5 mL tubes.
- Pipette tips: 1000, 200 and 10 μ L
- Pipettes
- C18 ZipTips or equivalent (e.g. C18 StageTips)

Reagents

- GuHCl (Guanidine hydrochloride), concentration: 6 M
- TCEP (Tris-(2-carboxyethyl)-phosphine), concentration: 100 mM
- CAA (Chloroacetamide), concentration: 100 mM
- Water, molecular grade
- Ethanol, 100% v/v
- ABC (Ammonium bicarbonate), concentration: 50 mM
- TFA (Trifluoroacetic acid), concentration: 5% v/v and 0.1% v/v
-  Sequencing Grade Modified Trypsin **Promega Catalog #V5111**
-  Sera-Mag Speedbead carboxylate-modified [E3] magnetic particles **Cytiva Life Sciences Catalog #65152105050250**
-  Sera-Mag™ SpeedBead Carboxylate-Modified [E7] Magnetic Particles **Cytiva Life Sciences Catalog #45152105050250**

Protocol materials

 Sequencing grade Trypsin **Promega Catalog #V5111**

 Sera-Mag Speedbead carboxylate-modified [E3] magnetic particles **Cytiva Life Sciences Catalog #65152105050250**

 Sera-Mag™ SpeedBead Carboxylate-Modified [E7] Magnetic Particles **Cytiva Life Sciences Catalog #45152105050250**

 Sequencing Grade Modified Trypsin **Promega Catalog #V5111**

Troubleshooting

Safety warnings

 Take care and adhere to all chemical safety data sheets.



Protein extraction

- 1 Prepare foodcrust samples by sampling into new Protein Low-Bind tubes. Suggested mass for each sample is 10-20 mg depending on sample availability.
- 2 Preheat a heat block to 65°C. 
- 3 Add 150 μ L 6 M GuHCl to each sample.
- 4 Briefly vortex to homogenise and centrifuge for 1 minute at 13k RPM. 
- 5 Incubate samples at 65°C for 1 hour on the heat block to extract the proteins from the foodcrust into the GuHCl. Then remove the samples from the heat block.  

Reduction & alkylation

- 6 Preheat a heat block to 99°C. 
- 7 Add 15 μ L of a solution containing 100 mM TCEP and 100 mM CAA. 
- 8 Briefly vortex to homogenise and centrifuge for 1 minute at 13k RPM. 
- 9 Incubate samples at 99°C for 10 minutes in a heat block, then remove the samples from the heat block and allow to cool down at room temperature for roughly 5 minutes.  

Bead clean up

- 10 Preheat a Thermomixer or equivalent heating and shaking device to 24°C or room temperature. 
- 11 Add 500 μ g of beads (i.e. 10 μ L of a 50 μ g/ μ L bead solution in ultrapure water) to each sample. Homogenise through gentle aspirating and dispensing with a pipette in the sample.  



Note: the ratio of protein to bead concentration should be 1:10 for optimal performance.

- 12 Add 175 μ L 100% ethanol to each sample and homogenise through pipetting.



Note: added ethanol volume should be equal to volume currently in the sample tube.

- 13 Incubate samples in a Thermomixer at 24°C at 1000 RPM for 5 minutes.
The proteins will now bind to the magnetic beads.



- 14 Place the samples on the magnetic separation rack and allow the beads to fully migrate for at least 2 minutes.

- 15 With the samples still on the magnetic rack, remove the supernatant.

Note: the supernatant can be discarded or saved to test for bead extraction efficiency.

- 16 Remove the samples from the magnetic rack and add 500 μ L of 80% ethanol. Resuspend the beads through gentle pipetting. Then place the beads back on the magnetic rack and allow the beads to migrate for at least 2 minutes.



- 17 With the beads on the magnetic rack, remove the supernatant and discard to waste.

- 18 Remove the samples from the magnetic rack and add 300 μ L of 80% ethanol. Resuspend the beads through gentle pipetting. Then place the beads back on the magnetic rack and allow the beads to migrate for at least 2 minutes.

- 19 With the beads on the magnetic rack, remove the supernatant and discard to waste.

- 20 Remove the samples from the magnetic rack and add 200 μ L of 80% ethanol. Resuspend the beads through gentle pipetting. Then place the beads back on the magnetic rack and allow the beads to migrate for at least 2 minutes.

- 21 With the beads on the magnetic rack, remove the supernatant and discard to waste. Ensure that as much supernatant as possible is removed.

- 22 Resuspend the beads in 100 μ L 50 mM ABC to elute the proteins from the beads and homogenise through gentle pipetting.

Separate aliquot for protein quantification (optional)

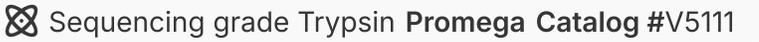
- 23 Place the samples on a thermomixer at 37°C and 750 RPM for 3 minutes.



- 24 Place the samples on the magnetic rack and allow the beads to migrate for at least 2 minutes. *
- 25 Remove 10 μL and store separately for protein quantification (i.e. Bradford assay or BCA).

Tryptic digestion

18h

- 26 Resuspend 20 μg of  with 100 μL of resuspension buffer to create a 0.2 $\mu\text{g}/\mu\text{L}$ solution. Add 1 μL of trypsin solution to each sample.  
- 27 Incubate the samples at 37°C and 750 RPM for 18 hours to digest. 18h
- 28 Remove the samples from the mixer and centrifuge at 13k RPM for 1 minute. 
- 29 Place the samples on the magnetic rack and allow the beads to migrate for at least 2 minutes. Transfer all the supernatant to new Protein Low-Bind tubes. The beads can now be discarded as all peptides are in the supernatant. 
- 30 Halt tryptic activity by acidifying with 10 μL 5% v/v TFA.  

Desalting clean up

- 31 Prime C18 ZipTips by aspirating 100 μL 50% ACN + 0.1% TFA. Discard solvent to waste. Repeat this step for a total of 2 primings.
- 32 Wash the ZipTip by aspirating 100 μL 0.1% TFA and discarding to waste. Repeat for a total of 2 washes.
- 33 Desalt samples by aspirating and dispensing 5-10 times in the sample. The peptides should now sit on the C18 filter of the ZipTip.
- 34 Wash the ZipTip by aspirating 100 μL 0.1% TFA and discarding to waste.

Repeat for a total of 2 washes.

- 35 Elute the peptides in a new Low-Bind tube by taking up 100 μL 50% ACN + 0.1% TFA and dispensing.
The peptides have now been eluted off the C18 filter into the solution.

Reagent preparation

- 36 Preparing a 100 mM TCEP and 100 CAA solution:
Weigh out 28.66 mg TCEP and 9.35 mg CAA in a darkened tube (or one wrapped with foil).
Add 1 mL ultrapure water and vortex for 1 minute to homogenise.

- 37 Preparing a 50 $\mu\text{g}/\mu\text{L}$ bead solution

- 37.1 Carefully invert the bottles of bead stock solutions and knock (!) them firmly on the counter to gain an equal distribution of the beads in the stock solution.

- 37.2 Pipette 500 μL of the

 Sera-Mag Speedbead carboxylate-modified [E3] magnetic particles **Cytiva Life Sciences Catalog #65152105050250**

into a tube and combine with 500 μL of

 Sera-Mag™ SpeedBead Carboxylate-Modified [E7] Magnetic Particles **Cytiva Life Sciences Catalog #45152105050250**

Place them on the magnetic rack and remove the supernatant.

- 37.3 Wash 3 times by removing the beads from the magnetic rack, resuspending in 1000 μL ultrapure water, homogenising through pipetting, placing back on the magnetic rack and removing the supernatant.

- 37.4 After the final wash resuspend the beads in 1000 μL ultrapure water and homogenise through gentle pipetting.

Protocol references

Developed from: Wilkin, S., Hagan, R., Hebestreit, S., Bleasdale, M., Nayak, A., Tang, L., Billings, T.N., Boivin, N. and Richter, K.K., 2021. *SP3 (Single-Pot, Solid-phase, Sample-Preparation) Protein Extraction for Dental Calculus* [online]. *protocols.io* <https://dx.doi.org/10.17504/protocols.io.bfgrijv6>

See also: Bleasdale, M., Richter, K.K., Janzen, A., Brown, S., Scott, A., Zech, J., Wilkin, S., Wang, K., Schiffels, S., Desideri, J. and Besse, M., 2021. Ancient proteins provide evidence of dairy consumption in eastern Africa. *Nature Communications*, 12(1), p.632.

Hughes, C. S., Foehr, S., Garfield, D. A., Furlong, E. E., Steinmetz, L. M., & Krijgsveld, J. (2014). Ultrasensitive proteome analysis using paramagnetic bead technology. *Molecular Systems Biology*, 10, 757. <https://doi.org/10.15252/msb.20145625>

Hughes, C. S., Moggridge, S., Müller, T., Sorensen, P. H., Morin, G. B., & Krijgsveld, J. (2019). Single-pot, solid-phase-enhanced sample preparation for proteomics experiments. *Nature Protocols*, 14(1), 68–85. <https://doi.org/10.1038/s41596-018-0082-x>

Cleland, T. P. (2018). Human Bone Paleoproteomics Utilizing the Single-Pot, Solid-Phase-Enhanced Sample Preparation Method to Maximize Detected Proteins and Reduce Humics. *Journal of Proteome Research*, 17(11), 3976–3983. <https://doi.org/10.1021/acs.jproteome.8b00637>

Acknowledgements

This work was supported by a Philip Leverhulme Prize awarded to Dr. Jessica Hendy. We thank Dr. Joannes Dekker for his assistance in typesetting this protocol.