

May 02, 2020

Version 1

SPARC_Duke_Grill_OT2-OD025340_VagusNerve_IHC_ChAT V.1

DOI

<https://dx.doi.org/10.17504/protocols.io.6hfhb3n>

J. Ashley Ezzell¹, Nicole A. Pelot¹, Kara A. Clissold¹, Warren M. Grill¹

¹Duke University



Nicole A Pelot

Duke University

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.6hfhb3n>

Protocol Citation: J. Ashley Ezzell, Nicole A. Pelot, Kara A. Clissold, Warren M. Grill 2020. SPARC_Duke_Grill_OT2-OD025340_VagusNerve_IHC_ChAT. protocols.io <https://dx.doi.org/10.17504/protocols.io.6hfhb3n>

License: This is an open access protocol distributed under the terms of the [Creative Commons Attribution License](#), which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: August 13, 2019



Last Modified: May 02, 2020

Protocol Integer ID: 26887

Keywords: Vagus nerve, peripheral nerve, immunohistochemistry, choline acetyltransferase, ChAT, vagal efferents, abdominal vagus nerve samples from rat, vagus nerve, abdominal vagus nerve sample, abdominal vagus nerve, od025340-vagusnerve-ihc-chat the protocol, od025340-vagusnerve-ihc-chat

Abstract

The protocol describes immunohistochemistry with anti-choline acetyltransferase, as it has been applied to cervical and abdominal vagus nerve samples from rats, pigs, and humans.

Materials

- Microscope slides with paraffin slices
- Xylene
- Ethanol
- Deionized water
- HIER Buffer L (Thermo, TA-135-HBL) for human samples or BioCare Borg Decloaker (BD1000MM) for rat/pig samples
- H₂O₂
- Tris buffer
- Tris Tween buffer
- DAKO Protein Block (X0909)
- Antibody Diluent OP Quanto (Thermo, TA-125-ADQ)
- Goat anti-choline acetyltransferase (Millipore, AB144P)
- Biotinylated SP-conjugated Affinipure donkey anti-goat IgG (H+L) (Jackson, 705-065-147)
- ABC Elite (Vector, PK-6100)
- DAB chromogen (Thermo, TA-125-QHDX)
- Harris hematoxylin (Thermo, 6765003)
- DPX mountant (Electron Microscopy Sciences, 13512)
- Microscope with color camera

Immunohistochemistry

- 1 Bake slides with sections of paraffin-embedded vagus nerve overnight at 50oC and then cool overnight.
- 2 Deparaffinize the slides and hydrate them to distilled water: xylene (2× 6 min), 100% ethanol (5 min), 95% ethanol (4 min), 70% ethanol (3 min), deionized water (2× 1 min).
- 3 Perform heat-induced epitope retrieval (HIER) at 120oC for 30 s (human) or at 95oC for 5 min (rat, pig) followed by 90oC for 10 s, using a buffer with pH 6.0 (Thermo, TA-135-HBL) for human samples or a buffer with pH 9.0 (BioCare, BD1000MM) for rat/pig samples.
- 4 Cool for 20 min at room temperature.
- 5 Rinse in deionized water (2× 2 min).
- 6 Block with 3% H₂O₂ diluted in deionized water for 10 min.
- 7 Rinse in deionized water (2× 2 min).
- 8 Rinse in Tris buffer (1× 2 min).
- 9 Block using DAKO Protein Block (X0909) for 10 min at room temperature.
- 10 Apply the primary antibody (goat anti-choline acetyltransferase, Millipore, AB144P) diluted in Thermo Antibody Diluent to a concentration of 1:100 for cervical & abdominal human samples, 1:50 for cervical pig samples, 1:100 for abdominal pig samples, and 1:200 for cervical & abdominal rat samples. Incubate overnight at 4oC.
- 11 Rinse in Tris Tween buffer (2× 2 min).
- 12 Rinse in Tris buffer (1× 2 min).

- 13 Apply the secondary antibody (biotinylated SP-conjugated Affinipure goat anti-rabbit IgG (H+L), Jackson, 111-065-144) diluted in Thermo Antibody Diluent to a concentration of 1:500, and incubate for 1 hour at room temperature.
- 14 Rinse in Tris Tween buffer (2× 2 min).
- 15 Rinse in Tris buffer (1× 2 min).
- 16 Apply ABC Elite (Vector, PK-6100) at a concentration of 1:50 for 30 min at room temperature.
- 17 Rinse in Tris Tween buffer (2× 2 min).
- 18 Rinse in Tris buffer (1× 2 min).
- 19 Apply DAB chromogen (Thermo, TA-125-QHDX) at room temperature for 2.5 min for cervical & abdominal human samples, 2 min for cervical pig samples, 3 min for abdominal pig samples, and 4 min for cervical & abdominal rat samples.
- 20 Rinse in deionized water (2× 2 min).
- 21 Counterstain using hematoxylin.
- 22 Dehydrate, clear, and coverslip using DPX mountant.

Microscopy

- 23 Each sample was imaged at 20x using a Nikon Ti2 microscope with a Photometrics Prime 95B-25MM camera (Nikon Instruments Inc.). We selected the best of four slices for each sample based on the quality of the slice (no tearing or fraying).