

Jul 11, 2019

SPARC Long-term exposure to intermittent hypoxia (or normoxia) using a custom in-cage computer controlled system

 In 1 collection

DOI

dx.doi.org/10.17504/protocols.io.2j3gcqn

Elisa Gonzalez-Rothi¹, Latoya Allen¹, Alec Simon¹, Yasin Seven¹, Marissa Ciesla¹, Gordon Mitchell¹

¹University of Florida

SPARC

Tech. support email: info@neuinfo.org



Elisa Gonzalez-Rothi

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.2j3gcqn>

Protocol Citation: Elisa Gonzalez-Rothi, Latoya Allen, Alec Simon, Yasin Seven, Marissa Ciesla, Gordon Mitchell 2019. SPARC Long-term exposure to intermittent hypoxia (or normoxia) using a custom in-cage computer controlled system. **protocols.io** <https://dx.doi.org/10.17504/protocols.io.2j3gcqn>

License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: May 02, 2019

Last Modified: July 11, 2019

Protocol Integer ID: 22875

Keywords: intermittent hypoxia, chronic intermittent hypoxia, acute intermittent hypoxia, IH, dAIH, CIH, oxygen, term exposure to intermittent hypoxia, term exposure to various intermittent hypoxia, various intermittent hypoxia, protocols in rat, normoxia, cage computer

Abstract

This protocol describes the procedures for conducting long-term exposure to various intermittent hypoxia (or normoxia) protocols in rats.

Troubleshooting

- 1 Rats are exposed to either long-term normoxia or IH treatments (various protocols) for up to 28 days
 - 1.1 Rats are housed in custom-designed PlexiGlas gas exposure cages (Volume~16L) with free access to food and water
 - 1.2 Cages are fitted with a custom cage top equipped to deliver protocol-specific gas (O₂ balanced with N₂) in specified cycle intervals via programmable mass-flow controller (Therapeutiq, Madison, WI).
 - 1.3 Cages are filled with low-dust emitting bedding (Tek-Fresh) and are changed 3 times per week
 - 1.4 Rats and gas exposure system are monitored at least 1x/hour during daily exposure to ensure system functionality and rat health/wellbeing
 - 1.5 Gas flow rate through cages is set to 16 L/min (~60 air changes per hour) to prevent CO₂ and ammonia buildup
 - 1.6 All gases are medical grade and are filtered for small microparticles and bacteria using a high purity compressed air filter (Parker Hannifin)
 - 1.7 Filtered air is humidified to reach a relative humidity of ~30-50% in the final gas mixture. The humidification system is comprised of a Nafion Fuel Cell Humidifier (Perma Pure, Lakewood, NJ) connected to the filtered air line and a water bath.
- 2 Exposure Protocols Include: Normoxia (Nx), daily acute intermittent hypoxia (dAIH), chronic intermittent hypoxia simulating mild sleep apnea (apnea-hypopnea equivalent = 6 episdoes/hour) (IH-5/5), or chronic intermittent hypoxia simulating moderate sleep apnea (apnea-hypopnea equivalent = 15 episdoes/hour) (IH-2/2)
- 3 Normoxia Exposure:
 - 3.1 Continuous gas flow of 21% O₂ for 8 hrs/day
- 4 Daily Acute Intermittent Hypoxia Exposure (dAIH)

- 4.1 10, 5-minute episodes of 10.5% O₂ alternating with 5 minute normoxic intervals (21% O₂) for a total duration of 1.5 hours/day (10 hypoxic episodes per day)
- 5 Chronic intermittent hypoxia simulating mild sleep apnea (IH-5/5): this protocol results in 6 hypoxic episodes per hour which is equivalent to the clinical criteria (apnea-hypopnea index, AHI) for diagnosing mild sleep apnea
- 5.1 5-minute episodes of 10.5% O₂ alternating with 5 minute normoxic intervals (21% O₂) for a total duration of 8 hours/day (6 hypoxic episodes per hour for 8 hours = 48 hypoxic episodes per day)
- 6 Chronic intermittent hypoxia simulating moderate sleep apnea (IH-2/2): this protocol results in 15 hypoxic episodes per hour which is equivalent to the clinical criteria (apnea-hypopnea index, AHI) for diagnosing moderate sleep apnea
- 6.1 2-minute episodes of 10.5% O₂ alternating with 2 minute normoxic intervals (21% O₂) for a total duration of 8 hours/day (15 hypoxic episodes per hour for 8 hours = 120 hypoxic episodes per day)