

Oct 26, 2023

Sputum Sample Processing for Single Cell Isolation and Live Recovery for Single Cell RNA Sequencing

DOI

<https://dx.doi.org/10.17504/protocols.io.j8nlko9k5v5r/v1>

Jason D Limberis¹, Alina Nalyvayko¹, Michele Tomasicchio², Shameem Jaumdally², anil pooran²,
Keertan Dheda², john.metcalfe¹

¹University of California, San Francisco; ²University of Cape Town, South Africa



Jason D Limberis

University of California, San Francisco

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.j8nlko9k5v5r/v1>

Protocol Citation: Jason D Limberis, Alina Nalyvayko, Michele Tomasicchio, Shameem Jaumdally, anil pooran, Keertan Dheda, john.metcalfe 2023. Sputum Sample Processing for Single Cell Isolation and Live Recovery for Single Cell RNA Sequencing.

protocols.io <https://dx.doi.org/10.17504/protocols.io.j8nlko9k5v5r/v1>

License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: September 11, 2023

Last Modified: October 26, 2023

Protocol Integer ID: 87649

Keywords: sputum, single cell isolation, single cell RNA sequencing, scRNAseq, infectious disease, clinical sputum samples for single live cell isolation, live recovery for single cell rna sequencing processing, single cell rna sequencing processing, single live cell isolation, clinical sputum sample, single cell, rna, cell, live recovery

Funders Acknowledgements:

UCB Catalyst

Grant ID: Chipotle

Abstract

Processing of clinical sputum samples for single live cell isolation for downstream applications.

Materials

✕ Zymo DNA/RNA Shield **Fisher Scientific Catalog #50-125-1706**

✕ Phosphate buffer solution (PBS)

✕ EDTA (0.5 M), pH 8.0 **Life Technologies Catalog #AM9260G**

✕ Bovine Serum Albumin (BSA) **Merck MilliporeSigma (Sigma-Aldrich) Catalog #A7906**

✕ Magnesium Chloride **Fisher Scientific Catalog #AC223210010**

✕ DMSO **Merck MilliporeSigma (Sigma-Aldrich) Catalog #472301**

✕ Fetal calf serum (FCS) **Gemini Bio-Products Catalog # 900-108**

✕ DNase I, Bovine Pancreas, >500 Kunitz U/mg **Gold Biotechnology Catalog #D-300**

✕ 100 g N-Acetyl-L-Cysteine **biorbyt Catalog #orb320385**

✕ Collagenase, Type IV, powder **Thermo Fisher Catalog #17104019**

Before start

Increasing the amount of sputum, by asking the patient to expectorate several time, will likely increase the final cell numbers.



Prepare Buffers

1 5mM EDTA in PBS

500ul 0.5M EDTA stock

49.5ml PBS. Store at 4C

10% BSA stock

Add 1g BSA to 10ml PBS, mix gently till dissolved

Filter sterilize & keep at 4C

50mM MgCl₂ solution

Add 238mg of MgCl₂ to 50ml of H₂O

10% DMSO/FCS freezing mixture

 5 mL DMSO

 45 mL fetal calf serum (FCS)

[M] 7.5 Mass Percent **NALC in PBS**

DNase

Aliquoted into 100ul aliquots at 224ku/ml and stored at -20C

Add 100ul to 40ml dissociation buffer to make a final conc of 0.56ku/ml.

Collagenase IV

Add  500 μ L to  40 mL dissociation buffer to make a final conc of 0.05mg/ml

Aliquoted  500 μ L at 4mg/ml and stored at -20C.

Make fresh for each use

1% BSA in PBS

 5 mL 10% BSA

 45 mL PBS

Dissociation buffer

 1.6 mL of [M] 50 millimolar (mM) MgCl₂

 100 μ L DNase

 500 μ L collagenase



🧪 37.8 mL PBS

Sputum collection

- 2 Collect an induced sputum, preferably

🧪 5 mL

 or more in volume, and process as below as soon as possible.
- 2.1 Weigh the sputum collection cup before and after the sputum collection.
- 2.2 Estimate the volume of sputum collected using a graduated container of the same size.
- 2.3 Record the viscosity (salivary, mucosalivary, purulent, mucopurulent)
- 2.4 Record the time of sputum collection

Bulk RNA sequencing

- 3 Record the time of sputum processing
- 4 Using a wide bore tip (or cut the bottom of a

🧪 1000 μ L

) draw up ~

🧪 200 μ L

 sputum. If the sample is too viscous for a pipette, use a syringe needle instead.
- 5 Add to a 1.5ml tube containing

🧪 500 μ L

 Zymo RNA Shield
- 6 Vortex tube for

🕒 00:00:10
- 7 Immediately transfer to

🌡️ -80 °C

10s

Make a single cell suspension (scRNAseq)



- 8 Add  40 mL dissociation buffer into the collection tube containing the sputum sample
- 9 Secure cap, manually invert the tube several times. Secure to a hula mixer (low setting) and incubate for  00:10:00 at  Room temperature . 20m

Note: Big clumps should dissolve after  00:10:00 . Some small fibers may will remain.
- 10 Add DTT to a final concentration of 1-2mM. Invert the tube several times and incubate for  00:05:00 5m
- 11 Pass the dissociated sample through a  100 μm strainer on top of a new  50 mL tube using an automated pipettor (do not pour).  350 rcf, 4°C, 00:10:00 10m
- 12 Wash the strainer with  2 mL of PBS
- 13 Spin tube in a sealed bucket at  250 rcf, 4°C, 00:10:00 10m
- 14 Remove  1000 μL supernatant and place in a cryovial for Luminex Cytokine assays. Add  140 μL of protease inhibitor and store at  -80 °C
- 15 Aspirate the remainder of the supernatant and discard.
- 16 Gently resuspend cell pellet in  1 mL EDTA/PBS using a wide bore  1000 μL tip
- 17 Count cells using Turks protocol followed by paraformaldehyde fixation to decontaminate

Long-term storage of cells (scRNAseq)

- 18 Pre-cool CellCool and cryovials at  4 °C for >  02:00:00 2h




- 19 Place cells On ice
- 20 Gently mix the cells
- 21 Centrifuge at 300 rcf, 4°C, 00:05:00 5m
- 22 Discard supernatant and resuspend cell pellets in pre-chilled (4 °C) 10% DMSO/FCS solution (add dropwise)
- 23 Dispense cell suspension in 500 µL aliquots into the pre-cooled cryovials and place into CoolCell at -80 °C


Expected results

24 Pilot samples

	A	B	C	D
Sample		S1	S2	S3
Viscosity		mucosal	mucosal	mucosal
Volume used (ml)		10	4	6
Total cells before freezing (10 ⁶)		2.2	0.9	2.9
Total cells recovered after thawing (2 weeks after frozen; 10 ⁶)		0.88	0.48	1.33
Percentage cell recovery post freezing (%)		40	53	46

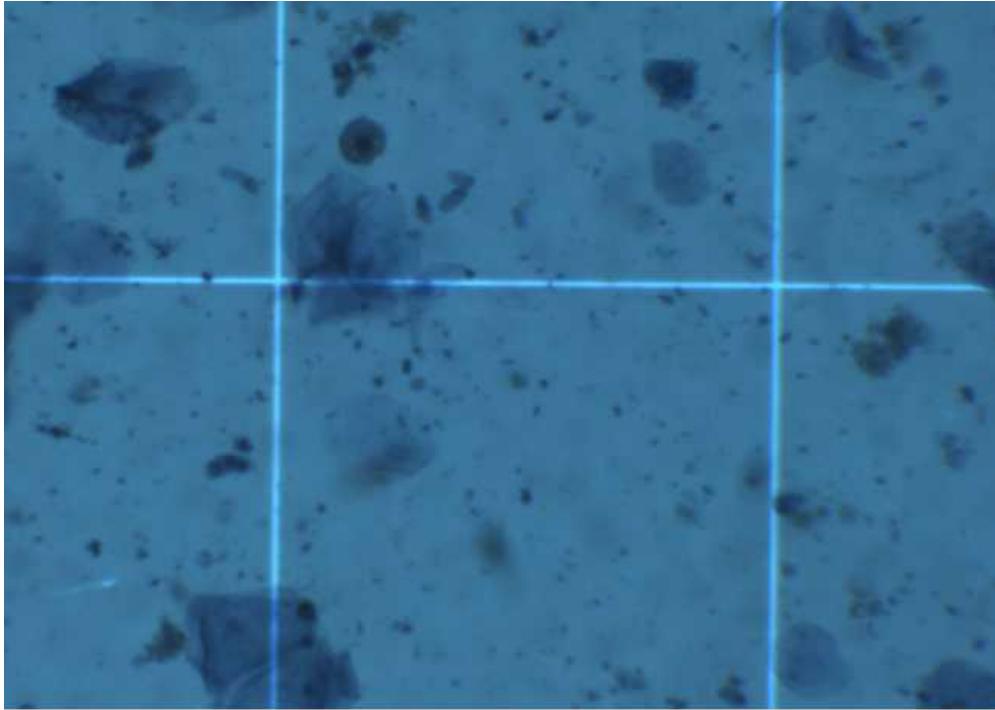

S2 400x

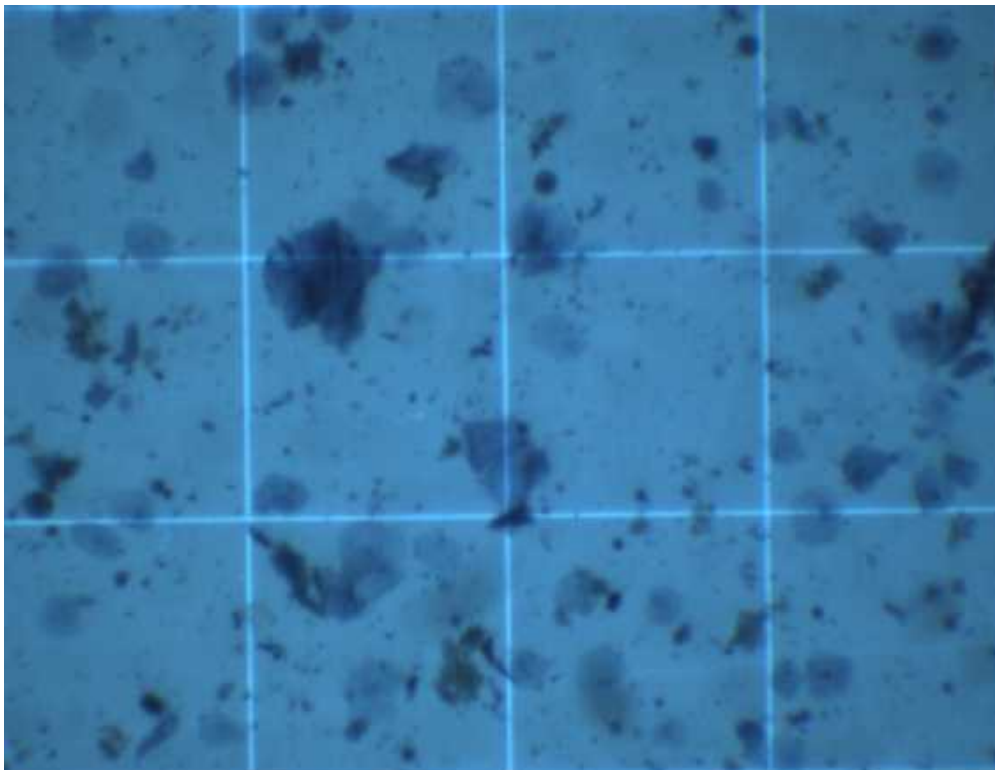

S3 100x

Clinical samples

	A	B	C	D	E	F	G	H	I	J	K	L	M	N
	Sample ID	Time to sputum processing (h. m)	weight of sputum (g)	volume of sputum (ml)	viscosity of sputum	Time to thaw (days)	Total cell count pre freeze (x10 ⁶)	Total cell count taken for ward pre freeze (x10 ⁶)	Total cell count post freeze (x10 ⁶)	Total cell count difference (x10 ⁶)	Live cell count	Dead cell count	Viability (%)	Viability of pre freeze total (%)
	CC 20	2.85	17.8	15	salivary	662	9.1	7	4.86	2.14	1.58	3.28	33	23
	CC 22	<6.00	2.1	2	mucoïd	658	1.3	1.3	0.60	0.71	0.15	0.44	26	12
	CC 23	2.10	2.9	3	mucoïd alivary	657	0.43	0.43	0.37	0.06	0.08	0.29	21	18
	CC 24	1.10	8.7	8	mucoïd alivary	657	1.1	1.1	0.81	0.30	0.19	0.62	23	17
	CC 25	4.90	0.7	0.9	mucoïd alivary	652	0.17	0.17	0.00	0.17	0.00	0.00	0	0
	CC 26	2.00	1.5	1.4	mucoïd alivary	652	1.13	1.13	0.73	0.40	0.17	0.57	23	15
	CC 27	2.20	3.2	3	mucoïd alivary	651	1.4	1.4	1.33	0.07	0.30	1.03	23	22
	CC 28	2.00	12	10	mucoïd	644	3.2	2	2.74	-0.74	0.29	2.45	11	15

	A	B	C	D	E	F	G	H	I	J	K	L	M	N
	CC 33	2.0 0	3	3	mu cos aliv ary	62 7	0.8	0.8	0.6 7	0.1 3	0.2 5	0.4 2	37	31
	CC 37	<6. 00	4.1	4	mu cos aliv ary	62 3	0.9 2	0.9 2	0.9 7	-0. 05	0.3 4	0.6 3	35	37