


Oct 14, 2019

## 🌐 Staining cells with IncuCyte Cytolight Rapid Dyes for flow cytometry or fluorescent microscopy

DOI

<https://dx.doi.org/10.17504/protocols.io.73nhqme>

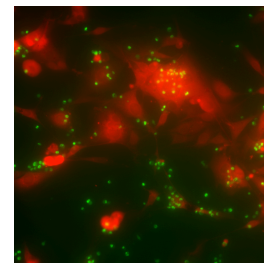

Elinor Gottschalk<sup>1</sup>, Bulent Arman Aksoy<sup>1</sup>, Pinar Aksoy<sup>1</sup>, Eric Czech<sup>1</sup>, Jeff Hammerbacher<sup>1</sup>

<sup>1</sup>Medical University of South Carolina

Hammer Lab

Tech. support phone: +18437924527 email: [arman@hammerlab.org](mailto:arman@hammerlab.org)



Elinor Gottschalk

### Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



**DOI:** <https://dx.doi.org/10.17504/protocols.io.73nhqme>

**Protocol Citation:** Elinor Gottschalk, Bulent Arman Aksoy, Pinar Aksoy, Eric Czech, Jeff Hammerbacher 2019. Staining cells with IncuCyte Cytolight Rapid Dyes for flow cytometry or fluorescent microscopy. **protocols.io**

<https://dx.doi.org/10.17504/protocols.io.73nhqme>

**License:** This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

**Protocol status:** Working

**We use this protocol and it's working**

**Created:** October 09, 2019

**Last Modified:** October 14, 2019

**Protocol Integer ID:** 28494

**Keywords:** cells with incucyte cytolight rapid dye, incucyte cytolight rapid dye, dye concentration for jurkat cell, staining cell, flow cytometry, dye concentration, melanoma cell, fluorescent microscopy this protocol, fluorescent microscopy, jurkat cell, labeling cell, cell

## Abstract

This protocol is for labeling cells with IncuCyte Cytolight Rapid dyes from Essen Bioscience. We optimized the dye concentration for Jurkat cells, pmel-1 T cells, and B16F10 melanoma cells.

## Guidelines

We optimized **IncuCyte® CytoLight Rapid Green Reagent for live cell cytoplasmic labeling** with Jurkat cells, pmel-1 T cells, and **IncuCyte® CytoLight Rapid Red Reagent** with B16F10 melanoma cells. The optimal concentrations (based on minimal decrease in cell viability and highest percentage of stained cells) were:

- Jurkat: 1  $\mu$ M (green)
- pmel-1 T cells: 0.11  $\mu$ M (green)
- B16F10 cells: 0.11  $\mu$ M (red)

We visualized the stained cells by flow cytometry or on the Keyence BZ-X710 fluorescence microscope equipped with GFP and Cy5 filters.

## Materials

### MATERIALS

 IncuCyte® CytoLight Rapid Green Reagent for live cell cytoplasmic labeling **Essen Biosciences Catalog #4705**

 1X PBS **VWR International (Avantor) Catalog #75800-986**

 IncuCyte® CytoLight Rapid Red Reagent for live cell cytoplasmic labeling **Essen Biosciences Catalog #4706**



- 1 Harvest cells and wash with 1X PBS
- 2 Centrifuge  350 x g , 5 minutes
  - 2.1 Resuspend cells in PBS at 1 million cells per ml
- 3 Add IncuCyte Dye
  - For pmel-1 T cells: we determined that 0.11  $\mu$ M of green dye was optimal
  - We made our working stock 11  $\mu$ M (so for every 1 ml of cell suspension, we add 10  $\mu$ l of working stock of dye)
  - For Jurkat cells, 1  $\mu$ M of green dye was optimal
  - For B16F10 cells, 0.11  $\mu$ M of red dye was optimal
- 3.1 Incubate the cells at  37 °C for  00:20:00 (we wrapped the tube in foil and placed it in a water bath).
  - invert the tube twice during incubation to mix the cells
- 3.2 Add a 6-fold volume of complete media (containing serum) to bind the excess dye.
- 3.3 Centrifuge  350 x g , 5 minutes
- 4 Resuspend cell in complete media at the desired concentration.
  - Two different cell types can be labeled with different dyes and then co-cultured and visualized
  - Stained T cells can be visualized migrating in collagen gel
- 5 Assay the cells via flowcytometry or fluorescence microscopy.  
Flow cytometry: FITC for green dye; APC for red dye  
Fluorescence microscopy: FITC filter for green dye; Cy5 filter for red dye