




Nov 23, 2022

🌐 Staining of Gfap, Iba1, and NeuN on PFA-fixed mouse brain sections

DOI

dx.doi.org/10.17504/protocols.io.4r3l27q5pg1y/v1

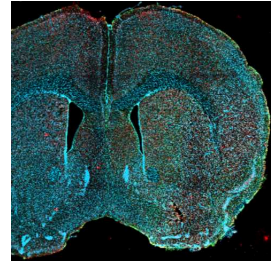

Daniel Manrique-Castano¹

¹Université Laval



Daniel Manrique-Castano

Université Laval

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.4r3l27q5pg1y/v1>

Protocol Citation: Daniel Manrique-Castano 2022. Staining of Gfap, Iba1, and NeuN on PFA-fixed mouse brain sections. **protocols.io** <https://dx.doi.org/10.17504/protocols.io.4r3l27q5pg1y/v1>

License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited



Protocol status: Working

Protocol established after several optimizing test

Created: November 23, 2022

Last Modified: November 23, 2022

Protocol Integer ID: 73191

Keywords: mouse brain section, fixed mouse brain section, neuron on pfa, gfap, mouse brain, protocol for gfap, pfa, staining protocol

Abstract

Staining protocol for Gfap, Iba1, and NeuN on PFA-fixed mouse brain sections.

Guidelines

Read the full protocol before starting the procedure.

Note that this protocol uses 3 hours (room temperature) incubation for primary antibodies.

Works with the listed antibodies and dilutions. For other references, preliminary tests are highly recommended.

Materials

Equipment

ImmEdge® Hydrophobic Barrier PAP Pen

NAME

Vector

BRAND

H-4000

SKU

<https://vectorlabs.com/products/histology/immedge-hydrophobic-barrier-pen>

LINK

Immunohistochemistry / Immunocytochemistry, Immunofluorescence, In situ hybridization SPECIFICATIONS

Permeabilization solution = [M] 0.3 % (v/v) Triton

⊗ Triton X-100 Merck MilliporeSigma (Sigma-Aldrich) Catalog #T8787-50ML / = [M] 0.1 % (v/v)

⊗ TWEEN 20 Merck MilliporeSigma (Sigma-Aldrich) Catalog #P7949 /

[M] 0.3 Molarity (M) ⊗ Glycine in **1x PBS**

Antibody buffer = [M] 0.1 % (v/v) ⊗ TWEEN 20 Merck MilliporeSigma (Sigma-Aldrich) Catalog #P7949 /

[M] 1 % (v/v) ⊗ Goat serum Merck MilliporeSigma (Sigma-Aldrich) Catalog #G9023

[M] 1 Mass / % volume

⊗ Bovine Serum Albumin (BSA) Merck MilliporeSigma (Sigma-Aldrich) Catalog #A7906

Primary antibodies:

⊗ GFAP Monoclonal Antibody (2.2B10) Thermo Fisher Scientific Catalog #13-0300

⊗ Anti Iba1 Rabbit antibody FUJIFILM Wako Pure Chemical Corporation Catalog #019-19741

⊗ Anti-NeuN Antibody Merck Millipore (EMD Millipore) Catalog #ABN91

Secondary antibodies:

⊗ Goat anti-Rat IgG (H L) Cross-Adsorbed Secondary Antibody Alexa Fluor 488 Thermo Fisher Scientific Catalog #A-11006



Cy3-AffiniPure Goat Anti-Rabbit IgG (H L) antibody **Jackson ImmunoResearch Laboratories, Inc. Catalog #111-165-003**



Goat anti-Chicken IgY (H L) Cross-Adsorbed Secondary Antibody Alexa Fluor Plus 647 **Thermo Fisher Scientific Catalog #A32933**



DAPI (46-Diamidino-2-Phenylindole Dilactate) **Invitrogen - Thermo Fisher Catalog #D3571**

Mounting media:



Fluoromount-G **Electron Microscopy Sciences Catalog #17984-25**



Protocol materials

- ✕ Anti Iba1 Rabbit antibody **FUJIFILM Wako Pure Chemical Corporation Catalog #019-19741**
- ✕ TWEEN 20 **Merck MilliporeSigma (Sigma-Aldrich) Catalog #P7949**
- ✕ Goat anti-Rat IgG (H L) Cross-Adsorbed Secondary Antibody Alexa Fluor 488 **Thermo Fisher Scientific Catalog #A-11006**
- ✕ Triton X-100 **Merck MilliporeSigma (Sigma-Aldrich) Catalog #T8787-50ML**
- ✕ TWEEN 20 **Merck MilliporeSigma (Sigma-Aldrich) Catalog #P7949**
- ✕ Goat anti-Chicken IgY (H L) Cross-Adsorbed Secondary Antibody Alexa Fluor Plus 647 **Thermo Fisher Scientific Catalog #A32933**
- ✕ Glycine
- ✕ GFAP Monoclonal Antibody (2.2B10) **Thermo Fisher Scientific Catalog #13-0300**
- ✕ Anti-NeuN Antibody **Merck Millipore (EMD Millipore) Catalog #ABN91**
- ✕ Cy3-AffiniPure Goat Anti-Rabbit IgG (H L) antibody **Jackson ImmunoResearch Laboratories, Inc. Catalog #111-165-003**
- ✕ Fluoromount-G **Electron Microscopy Sciences Catalog #17984-25**
- ✕ Goat serum **Merck MilliporeSigma (Sigma-Aldrich) Catalog #G9023**
- ✕ Bovine Serum Albumin (BSA) **Merck MilliporeSigma (Sigma-Aldrich) Catalog #A7906**
- ✕ DAPI (46-Diamidino-2-Phenylindole Dilactate) **Invitrogen - Thermo Fisher Catalog #D3571**
- ✕ Goat serum **Merck MilliporeSigma (Sigma-Aldrich) Catalog #G9023**
- ✕ Bovine Serum Albumin (BSA) **Merck MilliporeSigma (Sigma-Aldrich) Catalog #A7906**
- ✕ TWEEN 20 **Merck MilliporeSigma (Sigma-Aldrich) Catalog #P7949**
- ✕ TWEEN 20 **Merck MilliporeSigma (Sigma-Aldrich) Catalog #P7949**
- ✕ TWEEN 20 **Merck MilliporeSigma (Sigma-Aldrich) Catalog #P7949**
- ✕ Fluoromount-G **Electron Microscopy Sciences Catalog #17984-25**

Troubleshooting

Safety warnings

- ! DAPI is highly toxic. Handle it with care.



Tissue preparation and blocking

20m

- 1 Take out sections from -80 and place them in an incubator/plate for 00:20:00 at 37 °C . This step is performed to ensure tissue attachment to the crystal slides. 20m
- 2 Draw a hydrophobic barrier on each slide using **ImmEdge® Hydrophobic Barrier PAP Pen** and let dry for about 00:10:00 minutes. This will prevent buffers from escaping the tissue area. 10m

Note

There are other hydrophobic pens on the market. However, this one is recommended for its quality, consistency, and durability.

- 3 To initially rehydrate and permeabilize the tissue, place the slides in a slide jar containing the **permeabilization solution** with shaking for 00:30:00 30m

Note

The proposed permeabilization solution contains **glycine**, suitable to reduce PFA-related autofluorescence, especially in the 488 channel.

- 4 To prevent unspecific binding, incubate brain sections in **permeabilization solution** containing 5 % (v/v) 1h

Goat serum **Merck MilliporeSigma (Sigma-Aldrich) Catalog #G9023** and

1 Mass / % volume

Bovine Serum Albumin (BSA) **Merck MilliporeSigma (Sigma-Aldrich) Catalog #A7906**

for 01:00:00 at Room temperature .





Note

Please note that blocking serum must be chosen according to **secondary antibody species**. For the procedure depicted in this protocol, all secondary antibodies are raised in goat. However, donkey secondary antibodies will be also suitable.

Additionally, consider as well **avoiding BSA** when **primary antibodies** come from goat or sheep. BSA can generate unspecific background.

Antibody incubation

3h

- 5 When blocking is finished, decant the buffer (no washing is required) and incubate **primary antibodies** in **antibody buffer** for  03:00:00 at  Room temperature according to **table 1**.

3h



A	B	C	D	E
Antibody	Company	Reference	Specie	Dilution
Gfap	Invitrogen	13-0300	Rat	1:500
Iba1	Wako	019-19741	Rabbit	1:500
NeuN	Millipore	ABN91	Chicken	1:300




Table 1. Primary antibodies

Note

Please note that the antibody buffers **do not contain Triton X-100**. The reason is that this detergent tends to break the hydrophobic barrier, and the sections may not be adequately incubated. In addition, because sufficient permeabilization has already been performed previously, this detergent is not contemplated in this step.



Please note that the proposed antibody incubation lasts only **3 hours**. Although in most cases, staining protocols recommend overnight incubation for primary antibodies, the tests performed in our lab disclosed that, under the reported conditions, labeling for NeuN and Iba1 is superior at room temperature.

Also, the cited **Iba1 antibody** shows evident labeling and specificity superiority compared to others available in the market; therefore, highly recommended.

6 When primary antibody incubation is finished, wash the sections  00:05:00 5x with  0.05 % (v/v)  TWEEN 20 Merck MilliporeSigma (Sigma-Aldrich) Catalog #P7949 in **PBS**.

5m







7 Incubate **secondary antibodies** in  0.1 % (v/v)  TWEEN 20 Merck MilliporeSigma (Sigma-Aldrich) Catalog #P7949 for  01:00:00 at  Room temperature according to **table 2**.

1h



	A	B	C	D	E
	Antibody	Company	Reference	Channel	Dilution
	Goat anti-rat	Invitrogen	A11006	488	1:500
	Goat anti-rabbit	Jackson	111-165-003	Cy3	1:500
	Goat anti-chicken	Invitrogen	A32933	647	1:500
	Dapi	Invitrogen	D3571	405 (Dapi)	1:5000

Table 2. Secondary antibodies

8 When secondary antibody incubation is finished, wash the sections  00:05:00 3x with  0.05 % (v/v)  TWEEN 20 Merck MilliporeSigma (Sigma-Aldrich) Catalog #P7949 in **PBS**. Follow this with  00:05:00 2x washes with **PBS** to remove all detergent traces.

10m



9 Clean the remaining buffer on the slides using absorbent tissue and mount the sections with a drop of  Fluoromount-G Electron Microscopy Sciences Catalog #17984-25 .