


May 06, 2019

Standard iGEM Cell Measurement Protocol

 [ACS Synthetic Biology](#)

DOI

<https://dx.doi.org/10.17504/protocols.io.zgmf3u6>

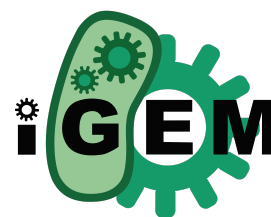

Richard Tennant¹, Paul Ruten¹

¹iGEM Measurement Committee

iGEM Measurement

Tech. support email: pauljruten@gmail.com



Paul Ruten

The University of Oxford

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.zgmf3u6>

External link: <https://2019.igem.org/Measurement>

Protocol Citation: Richard Tennant, Paul Ruten 2019. Standard iGEM Cell Measurement Protocol. **protocols.io**
<https://dx.doi.org/10.17504/protocols.io.zgmf3u6>

**Manuscript citation:**

Fedorec AJ, Robinson CM, Wen KY, Barnes CP, FlopR: An Open Source Software Package for Calibration and Normalization of Plate Reader and Flow Cytometry Data. ACS Synthetic Biology 9(9). doi: [10.1021/acssynbio.0c00296](https://doi.org/10.1021/acssynbio.0c00296)

License: This is an open access protocol distributed under the terms of the [Creative Commons Attribution License](#), which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

Created: March 24, 2019

Last Modified: August 07, 2019

Protocol Integer ID: 21741

Keywords: standard igem cell measurement protocol, igem interlab study, protocol, cell, standard, parts from the registry

Abstract

This is a protocol from the 2018 iGEM InterLab study which tests various 'standard' parts from the Registry.

Guidelines

Help with debugging:

- If you have measurements that are off scale ("OVERFLOW"), that data will not be usable. You need to adjust your settings so that the data will be in range and re-run your calibration.
- If your Abs600 measurements for your cell colonies are very close to that of your LB with Chloramphenicol, then your cells have probably not been transformed correctly or grown correctly.
- If your negative and positive control values are very close to each other, that probably means something has gone wrong in your protocol or measurement.

Materials

MATERIALS

✕ 1.5 mL Eppendorf tubes

✕ 96 well plate

✕ Falcon Tube (50 mL) **Fischer Scientific**

✕ Liquid LB medium

✕ Competent Cells

✕ Ice & ice bucket

✕ Chloramphenicol (25 mg/ml in EtOH)

STEP MATERIALS

✕ Chloramphenicol (25 mg/ml in EtOH)

Devices (from Distribution Kit, all in pSB1C3 backbone)

	Device	Part Number	Plate	Location
	Negative control	BBa_R0040	Kit Plate 7	Well 2D
	Positive control	BBa_J20270	Kit Plate 7	Well 2B
	Test Device 1	BBa_J364000	Kit Plate 7	Well 2F
	Test Device 2	BBa_J364001	Kit Plate 7	Well 2H
	Test Device 3	BBa_J364002	Kit Plate 7	Well 2J
	Test Device 4	BBa_J364007	Kit Plate 7	Well 2L
	Test Device	BBa_J36	Kit Plate	Well 2N

	ce 5	400 8	7	
	Test Devi ce 6	BBa_ J36 400 9	Kit Plate 7	Well 2P

The 96-well plate should preferably be black with a clear flat bottom.

Protocol materials

-  Liquid LB medium
-  1.5 mL Eppendorf tubes
-  Falcon Tube (50 mL) **Fischer Scientific**
-  Competent Cells
-  Ice & ice bucket
-  Chloramphenicol (25 mg/ml in EtOH)
-  96 well plate

Before start


Prior to performing this protocol you should complete the following calibration protocols:

- OD₆₀₀ Inter-equipment Conversion with LUDOX
- Particle Standard Curve with Microspheres
- Fluorescence Standard Curve with Fluorescein

Completion of the calibrations will ensure that you understand the measurement process and that you can take the cell measurements under the same conditions.

For all of these cell measurements, you must use the same plates and volumes that you used in your previous calibration protocols. You must also use the same settings (e.g., filters or excitation and emission wavelengths) that you used in your calibration measurements. If you do not use the same plates, volumes, and settings, the measurements will not be valid.

Read through this entire protocol carefully before you start your experiment and prepare any materials you may need. See the Guidelines tab on this page for debugging help.

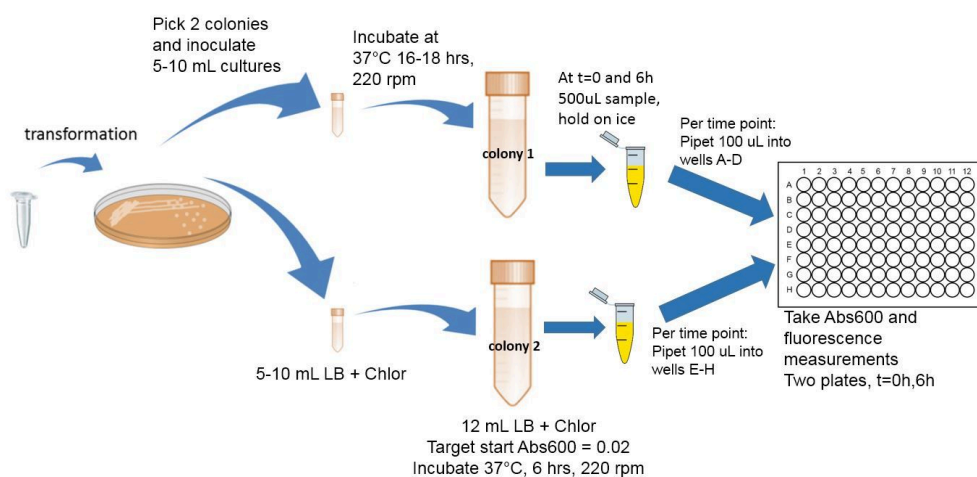



Day 1

- 1 Transform *Escherichia coli* DH5 α with these following plasmids (all in pSB1C3)

	Device	Part Number	Plate	Location
	Negative control	BBa_R0040	Kit Plate 7	Well 2D
	Positive control	BBa_I20270	Kit Plate 7	Well 2B
	Test Device 1	BBa_J364000	Kit Plate 7	Well 2F
	Test Device 2	BBa_J364001	Kit Plate 7	Well 2H
	Test Device 3	BBa_J364002	Kit Plate 7	Well 2J
	Test Device 4	BBa_J364007	Kit Plate 7	Well 2L
	Test Device 5	BBa_J364008	Kit Plate 7	Well 2N
	Test Device 6	BBa_J364009	Kit Plate 7	Well 2P



Note

Help Debugging Your Transformations:

- We STRONGLY recommend that you use the [iGEM protocol](#) to create your competent cells
- Once you have created your competent cells, we STRONGLY recommend that you measure the competency of your cells using the [Competent Cell Test Kit](#)
- Finally, we **strongly** recommend that you closely follow the [iGEM protocols](#) for resuspending DNA from the kit plates and performing the transformation
- Year after year, we have found that most teams are highly successful when they follow these protocols, even if alternative protocols are used within your lab. If you are having trouble transforming your test devices, please try the protocols above

Day 2

- 2 Pick 2 colonies from each of the transformation plates and inoculate in 5-10 mL LB medium + Chloramphenicol.

Grow the cells overnight (16-18 hours) at 37 °C and 220 rpm.



Liquid LB medium



Chloramphenicol (25 mg/ml in EtOH)

Day 3

- 3 Make a 1:10 dilution of each overnight culture in LB with Chloramphenicol (0.5 mL of culture into 4.5 mL of LB with Chloramphenicol)
- 4 Measure the Abs₆₀₀ of these 1:10 diluted cultures
- 5 Record the data in your notebook
- 6 Dilute the cultures further to a target Abs₆₀₀ of 0.02 in a final volume of 12 mL LB medium with Chloramphenicol in a 50 mL falcon tube

**Note**

Falcon tube should be amber, or covered with foil to block light

- 7 Take 500 μ L samples of the diluted cultures at 0 hours into 1.5 mL eppendorf tubes, prior to incubation. At each time point between 0 hours and 6 hours, you will take a sample from each of the 8 devices, two colonies per device, for a total of 16 eppendorf tubes with 500 μ L samples per time point, 32 samples total.

Place the samples on ice.

- 8 Incubate the remainder of the cultures at 37 °C and 220 rpm for 6 hours

- 9 Take 500 μ L samples of the cultures at 6 hours of incubation into 1.5 mL eppendorftubes. Place samples on ice.

Abs₆₀₀ and Fluorescence Measurement

- 10 At the end of the experiment, you should have two plates to read. You will have one plate for each time point: 0 and 6 hours. On each plate you will read both fluorescence and absorbance.

Samples should be laid out according to the plate diagram below.

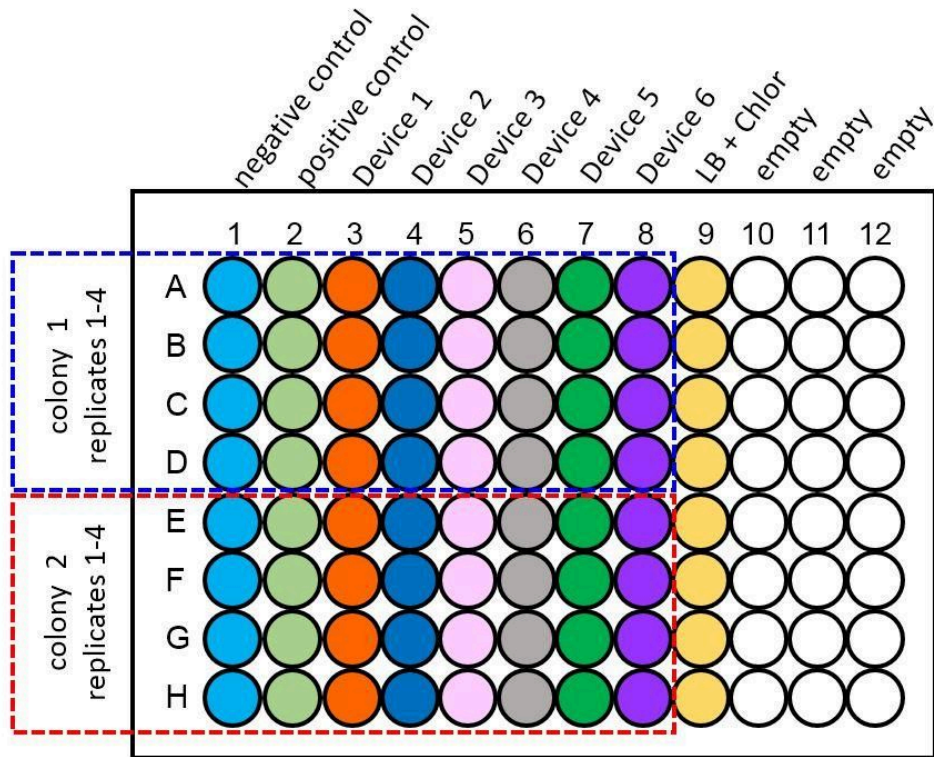


Plate layout for cell measurement calibration protocol

- 11 Pipette 100 μ L of each sample into each well. From 500 μ L samples in a 1.5 mL eppendorf tube:
 - 4 replicate samples of colony #1 should be pipetted into wells in rows A, B, C and D
 - Replicate samples of colony #2 should be pipetted into wells in rows E, F, G and H
- 12 Be sure to include 8 control wells containing 100 μ L each of only LB with Chloramphenicol on each plate in column 9, as shown in the diagram.

Set the instrument settings as those that gave the best results in your calibration curves (no measurements off scale). If necessary you can test more than one of the previously calibrated settings to get the best data (no measurements off scale).

Note

Instrument temperature should be set to room temperature (approximately 20-25 C) if your instrument has variable temperature settings

Congratulations!



13 You have now completed this calibration protocol