

Nov 06, 2024

Sucrose Gradient_Surface Isolation

DOI

<https://dx.doi.org/10.17504/protocols.io.rm7vz81m2vx1/v1>

Alexander Radaoui¹, hamiltonak^{1,2}, Karina L Konkrite¹

¹Children's Hospital of Philadelphia; ²University of Pennsylvania School of Medicine

Diskin Lab-CHOP

Tech. support phone: +12674253160 email: konkritek@email.chop.edu



Karina L Konkrite

Children's Hospital of Philadelphia

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.rm7vz81m2vx1/v1>

Protocol Citation: Alexander Radaoui, hamiltonak , Karina L Konkrite 2024. Sucrose Gradient_Surface Isolation. **protocols.io** <https://dx.doi.org/10.17504/protocols.io.rm7vz81m2vx1/v1>

License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited



Protocol status: Working

We use this protocol and it's working

Created: April 02, 2020

Last Modified: November 06, 2024

Protocol Integer ID: 35150

Keywords: sucrose gradient-surface isolation protocol, surface membrane of cell line, surface membrane, membrane, cell line, pdx tumor, patient tumor, surface

Funders Acknowledgements:

NIH

Grant ID: U54-CA232568

NIH

Grant ID: R01-CA204974

NIH

Grant ID: R01-CA237562

StandUp2Cancer- St Baldrick's Pediatric Dream Team

Grant ID: SU2C-AACR-DT1113

NIH

Grant ID: R03-CA230366

NIH

Grant ID: U01-CA199287

NIH

Grant ID: R35- CA220500

NIH

Grant ID: U01-CA263957

NIH

Grant ID: U01-CA263957

NIH

Grant ID: F31- CA225069

NIH

Grant ID: T32-CA009140

Disclaimer

This protocol needs prior approval by the users' institutional review board (IRB) or equivalent ethics committee(s).

Abstract

Protocol for extracting the surface membrane of cell lines or PDX tumors (or patient tumors) and preparing for mass spec.



Materials

MATERIALS

⊗ D-Sucrose **Fisher Scientific Catalog #BP220-1**

⊗ HEPES **Merck MilliporeSigma (Sigma-Aldrich) Catalog #H3375**

⊗ UltraPure™ 0.5M EDTA pH 8.0 **Invitrogen - Thermo Fisher Catalog #15575020**

⊗ Stericup-HV Sterile Vacuum Filtration System (0.45 um PVDF Membrane) **Merck MilliporeSigma (Sigma-Aldrich) Catalog #SCHVU11RE**

⊗ Sodium chloride 99 % crystalline powder PDV **VWR International (Avantor) Catalog #AAAA12313-0B**

⊗ Calcium chloride dihydrate **Merck MilliporeSigma (Sigma-Aldrich) Catalog #C3306**

⊗ Magnesium chloride solution BioUltra (2 M in H₂O) **Merck MilliporeSigma (Sigma-Aldrich) Catalog #68475**

⊗ Potassium Phosphate Dibasic **Merck MilliporeSigma (Sigma-Aldrich) Catalog #PX1570**

⊗ Trifluoroacetic Acid Optima™ LC/MS Grade **Fisher Scientific Catalog #A116-50**

⊗ Iodoacetamide **Merck MilliporeSigma (Sigma-Aldrich) Catalog #I1149**

⊗ Sodium Deoxycholate **Merck MilliporeSigma (Sigma-Aldrich) Catalog #D6750**

⊗ Dithiothreitol (DTT) 1M **Thermo Fisher Scientific Catalog #P2325**

⊗ Ammonium hydrogen carbonate ≥99.0% **Alfa Aesar Catalog #AA14249**

⊗ 13.2 mL Open-Top Thinwall Ultra-Clear Tube 14 × 89mm **Beckman Coulter Catalog #344059**

⊗ Amicon Ultra-0.5 Centrifugal Filter Unit (for up to 10 KDa) **Merck Millipore (EMD Millipore) Catalog #UFC501096**

⊗ Pierce™ BCA Protein Assay Kit **Thermo Fisher Scientific Catalog #23225**

⊗ Acetonitrile Optima™ LC/MS Grade **Fisher Scientific Catalog #A955-500**

⊗ 18 G x 1 1/2 in (Thin wall fill) **Becton Dickinson (BD) Catalog #305185**

⊗ 22 G x 1 1/2 in **Becton Dickinson (BD) Catalog #305156**

⊗ cOmplete™ Protease Inhibitor Cocktail **Roche Catalog #4693116001**

⊗ Sequencing Grade Modified Trypsin Frozen 5 × 20microg **Promega Catalog #PAV5113**

HOMOGENIZATION BUFFER (250 mM sucrose, 10 mM HEPES, 1 mM EDTA - for 1000 mL)

- 1) 85.58 g D-Sucrose
- 2) 2.38 g HEPES
- 3) 2 mL of 0.5 M EDTA
- 4) Bring to 900 mL with ddH₂O
- 5) Adjust the pH to 7.4 with either HCl or NaOH
- 6) Bring the final volume to 1000 mL with ddH₂O
- 7) Filter solution through a Durapore-Stericup 0.45 um filter

[Store solution at 4 °C]

HEPES BUFFER (115 mM NaCl, 1.2 mM CaCl₂, 1.2 mM MgCl₂, 2.4 mM K₂HPO₄, 20 mM HEPES - for 1000 mL)

- 1) 6.72 g NaCl
- 2) 133 mg CaCl₂
- 3) 114 mg MgCl₂
- 4) 418 mg K₂HPO₄
- 5) 4.77 g HEPES
- 6) Bring to 900 mL with ddH₂O
- 7) Adjust the pH to 7.4 with either HCl or NaOH
- 8) Bring the final volume to 1000 mL with ddH₂O
- 9) Filter solution through a Durapore-Stericup 0.45 um filter

[Solution is light sensitive and should be wrapped in foil and stored at 4 °C]

60% D-Sucrose Solution (for 500 mL)

- 1) 300.00 g D-Sucrose
- 2) Bring to 400 mL with ddH₂O and mix on stir plate with low heat.
- 3) Bring the final volume to 500 mL with ddH₂O
- 4) Filter solution through a Durapore-Stericup 0.45 um filter

[Store solution at 4 °C]

[Make diluting sucrose concentrations with the 60% D-Sucrose solution]

Sucrose Gradient Buffers (for 50 mL)

	New % of Sucrose	60% Sucrose (mL)	ddH ₂ O (mL)
	42.8	35.67	14.33



	42.3	35.25	14.75
	41.8	34.83	15.17
	41.0	34.17	15.83
	39.0	32.50	17.50
	37.0	30.83	19.17

Digestion Buffer (100 mM NH_4HCO_3 , 10% Sodium Deoxycholate, both made with ddH₂O - for 400 uL per sample)

- 1) 40 uL of 10% Sodium Deoxycholate (SDC)
- 2) 360 uL of 100 mM NH_4HCO_3

[The 10% SDC needs to be made fresh every time you are preparing a sample]

Reduction with 5 mM Dithiothreitol (DTT) in NH_4HCO_3 (50 mM NH_4HCO_3 , 1000 mM DTT - for 400 uL per sample)

- 1) 2 uL of 1000 mM DTT
- 2) 398 uL of 50 mM NH_4HCO_3

Alkylation with Iodoacetamide (IAA)

- 1) 92.48 g of IAA in 1000 mL makes 0.5 M IAA

[I would make a smaller 250 uL - 500 uL stock solution of 0.5 M IAA, depending on the number of samples]

[Adding 16 uL of 0.5 M IAA to one 400 uL sample makes the final concentration 20 mM]

[0.5 M IAA needs to be made fresh every time it is used]



[IAA is extremely light sensitive and needs to be stored in the dark at  4 °C]

Safety warnings

-  This protocol needs prior approval by the users' institutional review board (IRB) or equivalent ethics committee(s).

DAY 1

- 1 After all buffers are prepared, retrieve PDX/tumor samples or cells and thaw on ice. Add 1 tablet of **Roche cOmplete™ Protease Inhibitor Cocktail EASYpack** per 50 mL of homogenization buffer. Vortex to dissolve tablet.
 - For cells, aim to have roughly **100 million**.
 - For solid PDX tumor, aim to have **between 250 and 350 milligrams**.
- 2 Resuspend each sample in 10 mL of **HOMOGENIZATION BUFFER** in a 50 mL conical and use an electric TissueRuptor® II (from Qiagen) to dissociate tumor chunks into a fully fluid and homogenous mixture with no particulate or remaining solid tissue.
 - **Keep samples on ice throughout!**
 - Be sure to autoclave the TissueRuptor stems before use and change stems between samples.
- 3 Retrieve a second 50 mL conical and pass the sample through an 18-gauge needle into the second conical at least once.
 - If passaging is easy, move on to the next step. If syringe becomes clogged, further electric homogenization may be necessary.
- 4 Pass the sample through a 22-gauge needle between the first and second 50 mL conical 3 times.
 - If sample becomes clogged in the 22-gauge syringe, further electric homogenization may be necessary.
- 5 Transfer fully homogenized samples to 15 mL conicals and centrifuge at  1000 x g, 4°C for  00:15:00 .
- 6 Collect the supernatant, which is the **Post-Nuclear Cytosol (PNC)**, in a new 15 mL conical.
 - **Keep samples on ice!!!**
- 7 Retrieve Beckman Coulter® Ultra-Clear Centrifuge Tubes, rotor (SW-41-Ti), sample holders, and necessary equipment for centrifugation. Add 2 mL of 60% sucrose as a cushion to the ultra-clear tubes slowly with a P-1000. Layer the first 2 mL of PNC supernatant from your samples on top of the 60% sucrose cushion **VERY SLOWLY** and carefully so as not to disturb the 60% gradient layer (can use a serological pipette).
 - **Samples running opposite of each other on the rotor should be balanced in their respective holders to the 100th decimal with homogenization buffer!**
- 8 Centrifuge samples at  100000 x g, 4°C for  01:00:00 .

- 9 After centrifugation, use a serological pipette to carefully take out roughly 75% of the homogenization buffer supernatant making sure not to disturb the protein layer on top of the 60% cushion. Then using a P-200, harvest the **Crude Plasma Membrane (CPM)** fraction from the top of the 60% sucrose cushion and place in a new ultra-clear centrifuge tube.
- You can use a swirling technique with the P-200 to suck up the CPM fraction.
 - Try to **minimize the amount of 60% sucrose carry over** when transferring the CPM fraction to new ultra tubes.
- 10 Using a P-1000, set the volume to 750 uL and add 1.5 mL of each sucrose gradient on top of the sample and each other **VERY SLOWLY in a "tear drop" style fashion** starting with 42.8% sucrose, then 42.3%, then 41.8%, then 41.0%, then 39.0%, and lastly 37.0% sucrose at the top.
- You can fill the rest of the tubes (**SLOWLY**) and balance opposite facing samples to the 100th decimal with 37% sucrose.
 - It is **VERY IMPORTANT** to add your consecutive gradients slowly and steadily so as to have well-defined layer separation and prevent mixing of the gradients.
- 11 Centrifuge samples at  100000 x g, 4°C overnight for a minimum of  18:00:00 .
- I would recommend setting the ultra-centrifuge to "hold" during this spin in case you are doing other experiments.

DAY 2

- 12 The next day recover the **Enriched Plasma Membrane (EPM)** fraction from the top of the 37% sucrose cushion using a P-200 and place in a new ultra-clear centrifuge tube.
- You can use a swirling technique with the P-200 to suck up the EPM fraction.
 - Retrieve as much of the EPM fraction on top of the 37% cushion as possible, while also **limiting the amount of 37% sucrose carry-over**, and place into new ultra-clear tubes. The key is to not get greedy!
- 13 Fill up the ultra-tubes with **HEPES BUFFER** and balance opposing tubes to the 100th decimal before spinning.
- 14 Centrifuge samples at  150000 x g, 4°C for a minimum of  02:30:00 .
- Sometimes you will have to centrifuge your samples anywhere between  02:30:00 and  05:00:00 during this step to guarantee total pellet sedimentation of the available EPM.
 - **Would recommend starting with a 3-hour spin every time.**



- 15 Decant the HEPES supernatant out of the ultra-clear tubes being careful not to disturb the protein pellet. **Resuspend the pellet in 400 uL of DIGESTION BUFFER** (10% SDC + 100 mM NH_4HCO_3) and vortex.
- After resuspension with digestion buffer you can store your samples in  -80 °C , however, **it is ideal to complete the protocol up through the trypsin digestion.**
 - It would also be helpful to run a BCA Assay or other protein quantifications to determine the protein concentration per sample and determine how much trypsin to use for each sample.
 - If pellets are not solvating into the digestion buffer well, you can sonicate the samples on ice for  00:10:00 .
- 16 Retrieve Amicon® 0.5 mL Centrifugal Filter Units (for up to 10 KDa). Apply samples to the filter held in eppendorf tubes and spin down the samples between  8000 x g and  12000 x g for  00:10:00 at  4 °C to bind protein to the filter column.
- A small portion of liquid sample might remain at the bottom of the filter unit, but this is okay.
- 17 Decant the flow-through waste from the collection tube and wash the filters with 400 uL of ddH₂O by spinning samples down between  8000 x g and  12000 x g for  00:10:00 at  4 °C . **Repeat this step a second time.**
- Again, a small portion of H₂O might remain at the bottom of the filter unit, but this is okay.
- 18 Next, you want to perform a **REDUCTION** on your protein samples using a **5 mM Dithiothreitol (DTT)** solution made in **50 mM NH_4HCO_3** . Add 400 uL of the 5 mM DTT solution to each sample's filter unit and let **incubate at room temperature** for  01:00:00 for full reduction to occur. **DO NOT SPIN DOWN AFTER.**



- 19 You now want to cap/block all the reduced sulfide bonds on the proteins by **ALKYLATION with IODOACETAMIDE (IAA)** in order to prevent DTT from reverse reacting and reforming disulfide bonds. You want a ratio of **4 IAA equivalents for every 1 DTT equivalent** for full deactivation. Therefore, you will need a minimum of **20 mM of IAA to combat the 5 mM of DTT per sample**. Using a **stock solution of 0.5 M IAA** (made with ddH₂O), you will need roughly **16.67 uL of IAA** per each 400 uL sample. Add directly to the filter units with a P-20 and **INCUBATE IN THE DARK FOR**  00:30:00 **AT ROOM TEMPERATURE**.
- 0.5 M of IAA should be made fresh with ddH₂O for every new experiment.
- 20 After the incubation, spin down the samples at  12000 x g, 4°C for  00:05:00 and decant out the waste from the collection tube.
- You may have to spin the samples for up to  00:15:00 in order for the waste to fully pass through the column into the collection tube.
- 21 After quantifying your total protein, make a 400 uL solution of **50 mM NH₄HCO₃** that includes the appropriate amount of **TRYPSIN** for each sample. A **ratio of 1 ug of trypsin to 20 ug of protein** is our standard ratio. Add the respective trypsin solutions to each sample and **let the reaction sit overnight at room temperature for full peptide digestion**.
- It would also be useful to test the pH of the samples after trypsin is added to make sure the solution is at a  8 . This is ideal for optimal enzymatic activity.
 - Some reagents require you to add **TEAB** to the trypsin before creating the 50 mM NH₄HCO₃ solutions for each sample. The trypsin used in the materials section does NOT require this.

DAY 3

- 22 Using a P-200, lightly resuspend the solution in the filter to make sure your peptides are mixed and detached from the column. Retrieve fresh centrifuge tubes and carefully tilt the sample in the filters upside-down into the new tubes. Centrifuge the upside-down filters in the new collection tubes at  12000 x g, 4°C for  00:02:00 . This ensures that the total amount of peptides fall into the new collection tube.
- 23 In order to **promote peptide binding to the stage tips**, use **LC/MS-Grade TRIFLUOROACETIC ACID (TFA)** to acidify your samples. Start by adding 3 uL - 5 uL of TFA to the samples inside a fume hood. You want to make sure the pH of your sample is  2 . Let the samples **incubate on the bench at room temperature** for  00:30:00 .
- TFA is highly corrosive and will burn your skin. Handle with care IN A FUME HOOD.



- TFA also deactivates the trypsin by quenching the reaction as well as pulling the SDC out of the sample solution as a precipitate.
- While the samples sit for 30 minutes, you can begin **STAGE TIPPING with CARBON-18 RESIN** for your samples.

- 24 Spin down the samples after the incubation at **maximum speed** for  00:15:00 at  4 °C . While this occurs you can continue to **STAGE TIP with C-18 RESIN**. About 4-6 punches of the resin per sample are optimal to pack a P-200 pipette tip. Using pipette tip holders, or **ADAPTERS**, you can place the tips in fresh eppendorf tubes.
- Punches are easily made using P-1000 pipette tips that have the tip-ends snipped around half an inch.
 - The C-18 resin helps to cleanse the sample of excess salt and contaminants.
- 25 Wash the stage tips with **100 uL of 100% Acetonitrile** and spin them down at  2000 x g for  00:02:00 . Decant the waste.
- 26 Wash the stage tips with **100 uL of 0.1% TFA** (diluted with ddH₂O) and spin them down at  2000 x g for  00:02:00 . Decant the waste.
- 27 Add the sample to each of their own stage tips and spin them down at  2000 x g for  00:01:00 so they bind to the resin. Decant the waste.
- 28 Wash the samples with **100 uL of 0.1% TFA** and spin them down at  2000 x g for  00:02:00 . Decant the waste.
- 29 Finally, elute your samples into fresh eppendorf tubes by using **100 uL of 70% Acetonitrile** and centrifuging them at  1500 x g for  00:02:00 .
- The samples will be put into the speed vacuum next, so it is okay if they start to dry out.
- 30 Place your samples in the speed vacuum to dry down the samples. This can take anywhere between 2-6 hours.
- Freeze back the speed vacuumed samples in  -20 °C until ready to run on the mass spectrometer.