

Apr 07, 2023

Synthesis and preparation of DOPA pheomelanin and eumelanin

DOI

<https://dx.doi.org/10.17504/protocols.io.4r3l27j4qg1y/v1>

Kazumasa Wakamatsu¹

¹Institute for Melanin Chemistry, Fujita Health University, Toyoake, Japan



Jackson Schumacher

Massachusetts General Hospital

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.4r3l27j4qg1y/v1>

Protocol Citation: Kazumasa Wakamatsu 2023. Synthesis and preparation of DOPA pheomelanin and eumelanin. **protocols.io** <https://dx.doi.org/10.17504/protocols.io.4r3l27j4qg1y/v1>

License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: April 07, 2023

Last Modified: December 13, 2024

Protocol Integer ID: 80143

Keywords: ASAPCRN, preparation of dopa pheomelanin, preparing dopa pheomelanin, dopa pheomelanin, dopa eumelanin, using tyrosinase, pheomelanin, eumelanin, synthesis, preparation

Funders Acknowledgements:

From cancer associations to altered immunity in the pathogenesis of Parkinson's disease

Grant ID: ASAP-000312

Abstract

This is the protocol for synthesizing and preparing DOPA pheomelanin and DOPA eumelanin using tyrosinase.

Attachments



pcmr12121-sup-0001-

a...

3.5MB

- 1 As precursor(s) of melanin, 1) l-dopa, 2) DHI, 3) a mixture of DHI and DHICA, 4) DHICA, 5) a mixture of l-dopa and l-cysteine, and 6) 5SCD in the presence of l-dopa is used.
- 2 To a solution of the precursor(s) (1 mmol) in 0.1 M sodium phosphate buffer, pH 6.8 (98 ml), add a solution of mushroom tyrosinase in the same buffer (2 ml) at 25°C.
- 2.1 The amount of mushroom tyrosinase (Sigma-Aldrich, 4270 U/mg) is either 50,000 U or 100,000 U.
- 3 Carry out oxidation with vigorous (maximal, but avoiding splashing) stirring in a 500 ml Erlenmeyer flask.
- 4 Stop oxidation (1-4) at 4 h by adding 6 M HCl (2 ml) to make pH 1 and keep the mixture at 4°C overnight to ensure complete precipitation of the pigment.
- 5 Collect the precipitate by centrifugation at 3,000 rpm for 10 min and wash with 0.1 M HCl (40 ml x 3).
- 6 For the preparation of pheomelanins (5 and 6), acidify the oxidation mixture with acetic acid (3 ml) to make pH 3, and wash the precipitate with 1% acetic acid (40 ml x 3).
- 7 Dry the melanin powder by lyophilization and equilibrate with moisture in a desiccator with a saturated CaCl_2 solution.
- 8 Prior to treatment, suspend the synthetic DOPA pheomelanin (DOPA:Cys = 1:1) and DOPA eumelanin (DOPA:Cys = 1:0) in sterile PBS and sonicate overnight.

Protocol references

d'Ischia, M., Wakamatsu, K., Napolitano, A., Briganti, S., Garcia-Borron, J. C., Kovacs, D., ... & Ito, S. (2013). Melanins and melanogenesis: methods, standards, protocols. *Pigment cell & melanoma research*, 26(5), 616-633.