



Oct 07, 2025

Synthetic peptide-based indirect ELISA using S1c (of the sequence CPVAIHADQLTPTWRVYSTC) as the antigen, performed on regular polystyrene plates

Forked from a deleted protocol

DOI

<https://dx.doi.org/10.17504/protocols.io.6qpvrw753lmk/v1>

Brian Andrich Pollo^{1,2}, Ruby Anne King^{3,1}, Fresthel Monica M. Climacosa^{4,1}, Salvador Eugenio C. Caoili¹

¹Biomedical Innovations Research for Translational Health Science (BIRTHS) Laboratory, Department of Biochemistry and Molecular Biology, College of Medicine, University of the Philippines Manila;

²School of Medicine, The Manila Times College of Subic;

³Department of Science and Technology - Philippine Council for Health Research and Development (DOST-PCHRD);

⁴Department of Medical Microbiology, College of Public Health, University of the Philippines Manila

Brian Andrich Pollo: corresponding author;

Coronavirus Method De...



Brian Andrich L. Pollo

University of the Philippines Manila

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.6qpvrw753lmk/v1>

Protocol Citation: Brian Andrich Pollo, Ruby Anne King, Fresthel Monica M. Climacosa, Salvador Eugenio C. Caoili 2025. Synthetic peptide-based indirect ELISA using S1c (of the sequence CPVAIHADQLTPTWRVYSTC) as the antigen, performed on regular polystyrene plates. **protocols.io** <https://dx.doi.org/10.17504/protocols.io.6qpvrw753lmk/v1>

License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: October 06, 2025

Last Modified: October 07, 2025

Protocol Integer ID: 229158

Keywords: SARS-CoV-2, COVID-19, Spike protein, Synthetic peptides, Enzyme-Linked Immunosorbent Assay, ELISA, Antibody detection, Serologic tests, Protein A, Peroxidase, TMB substrate, Peptide antigens, Immunoassay methods, Polystyrene microplates, Carbonate-bicarbonate buffer, Tween 20, Skimmed milk, Blocking agents, Phosphate-Citrate Buffer, Absorbance spectrophotometry, Diagnostic assay development, Immunodiagnostics, Seroprevalence studies, based indirect elisa, indirect elisa, antibody detection in serum sample, immunosorbent assay, linked immunosorbent assay, primary antibody incubation, standard elisa plate, immobilized antigen, antibody detection, antigen, serum sample, synthetic peptide, milk in pbst, based indirect enzyme, elisa, cpvaihadqltptwrvystc, peptide s1c, enzyme, indirect enzyme, sequence cpvaihadqltptwrvystc, regular polystyrene plates this protocol, polystyrene plate, assay

Funders Acknowledgements:

Dissertation grant

IDC 211 grant

Abacov project

NIH faculty grant



Disclaimer

DISCLAIMER – FOR INFORMATIONAL PURPOSES ONLY; USE AT YOUR OWN RISK

The protocol content here is for informational purposes only and does not constitute legal, medical, clinical, or safety advice, or otherwise; content added to **protocols.io** is not peer reviewed and may not have undergone a formal approval of any kind. Information presented in this protocol should not substitute for independent professional judgment, advice, diagnosis, or treatment. Any action you take or refrain from taking using or relying upon the information presented here is strictly at your own risk. You agree that neither the Company nor any of the authors, contributors, administrators, or anyone else associated with **protocols.io**, can be held responsible for your use of the information contained in or linked to this protocol or any of our Sites/Apps and Services.

Abstract

This protocol describes a synthetic peptide-based indirect **enzyme-linked immunosorbent assay** (ELISA) using the peptide **S1c (sequence: CPVAIHADQLTPTWRVYSTC)** as the immobilized antigen. The assay is designed for antibody detection in serum samples and uses regular polystyrene microplates. Plates are coated overnight at 4 °C with 100 µL/well of S1c in carbonate–bicarbonate buffer (pH 9.6), then washed thrice with 0.05% Tween 20 in PBS (PBST). After a 15 min equilibration to room temperature, wells are blocked with 200 µL/well of 5% skimmed milk in PBST for 1 h at 37 °C and washed again three times. Primary antibody incubation is performed for 1 h at room temperature using 100 µL/well of sera diluted 1:100 in 0.05% skimmed milk–PBST, followed by washing. Detection is achieved with Protein A–peroxidase (0.5 µg/mL) diluted 1:2000 in PBST, incubated for 1 h at room temperature. After a final wash, the reaction is developed with 0.1 mg/mL TMB in phosphate–citrate buffer (pH 6.0) for 1 h at room temperature, stopped with 50 µL of 1 M H₂SO₄, and absorbance is measured at 450 nm. This optimized workflow enables reliable detection of anti-S1c antibodies using accessible reagents and standard ELISA plates.



Coat

- 1 overnight incubation at 4°C
100µL/well, Carbonate-bicarbonate buffer in water, pH 9.6

Wash

- 2 3x at 5 min interval
0.05% Tween 20 in 1X PBS (PBST)
freeze

Block

- 3 *leave plates at RT for 15min before blocking*
incubation at 37°C for 1hr
200µL/well, 5% skimmed milk in PBST

Wash

- 4 3x at 5 min interval
0.05% Tween 20 in 1X PBS

Primary antibody

- 5 incubation at RT for 1hr
100µL/well, 1:100 sera in dilution buffer (0.05% skimmed milk in PBST)

Wash

- 6 3x at 5 min interval
0.05% Tween 20 in 1X PBS

Secondary antibody

- 7 incubation at RT for 1hr
50µL/well, 1:2000 Protein A-peroxidase (0.5ug/ml) in dilution buffer

Wash



- 8 3x at 5 min interval
0.05% Tween 20 in 1X PBS

Substrate

- 9 incubation at RT for 1hr
50 μ L/well, 0.1mg/mL TMB in Phosphate Citrate Buffer in water, pH 6.0

Stop

- 10 50 μ L/well, 1M H₂SO₄
read at 450nm