

Sep 09, 2020

T4H protocol

DOI

dx.doi.org/10.17504/protocols.io.bk5iky4e

Ophir Auslaender¹, kalle.levon²

¹NYU Grossman School of Medicine; ²NYU Tandon School of Engineering

Tech4Health



kalle.levon

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.bk5iky4e>

Protocol Citation: Ophir Auslaender, kalle.levon 2020. T4H protocol. **protocols.io**
<https://dx.doi.org/10.17504/protocols.io.bk5iky4e>

License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: September 08, 2020

Last Modified: September 09, 2020

Protocol Integer ID: 41866

Keywords: t4h protocol, protocol

Troubleshooting

- 1 Clean Au chip: insert in piranha solution (75% H_2SO_4 , 25% H_2O_2), wash with DI water, dry with nitrogen gas.
- 2 Mix the solution for imprinting in 19/1 volume ratio: 19 for the analyte, 1 for the monolayer former at self-assembled monolayer (SAM) concentration.
- 3 Place Au chip in the imprinting solution.
Incubate overnight RT.
Wash multiple times with 3M NaCl solution.
Leave overnight in DI water.
- 4 At this point the chip can be sealed for future use.
Prior to measurement chip is inserted in PBS for 2 hours.
- 5 The chip is connected as the working electrode in a potentiostat circuit.
A reference electrode is added to the reaction vessel and recording of the open-circuit potential (OCP) is started.
When the voltage stabilizes, the analyte is added.
- 6 Adding the analyte shifts the OCP if the analyte is recognized by the imprinted working electrode.

In the calibration phase, controlled serial dilutions of analyte are added, and resulting OCP values are recorded. The relationship between analyte concentration and OCP is established from these calibration measurements, and is used for the detection of unknown analytes.