May 17, 2018

## Tail-flick test

DOI

[dx.doi.org/10.17504/protocols.io.p72drqe](https://dx.doi.org/10.17504/protocols.io.p72drqe)

Sunamita Lima<sup>1</sup>

<sup>1</sup>Unifesp

Sunamita Lima

### Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS

DOI: <https://dx.doi.org/10.17504/protocols.io.p72drqe>

**Protocol Citation:** Sunamita Lima 2018. Tail-flick test. **protocols.io** <https://dx.doi.org/10.17504/protocols.io.p72drqe>

**Manuscript citation:**

***Pharmacological characterization of crotamine effects on mice hind limb paralysis employing both ex vivo and in vivo assays: insights into the involvement of voltage-gated ion channels in the crotamine action on skeletal muscles***

Sunamita de Carvalho Lima<sup>1</sup>, Lucas de Carvalho Porta<sup>1</sup>, Álvaro da Costa Lima<sup>1</sup>, Joana D'Arc Campeiro<sup>1</sup>, Ywlliane Meurer<sup>2</sup>, Nathália Bernardes Teixeira<sup>3</sup>, Thiago Duarte<sup>1</sup>, Gisele Picolo<sup>3</sup>, Rosely Oliveira Godinho<sup>1</sup>, Regina Helena Silva<sup>1</sup>, Mirian Akemi Furuie Hayashi<sup>1, \*</sup>

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**License:** This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

**Protocol status:** Working

**We use this protocol and it's working**

**Created:** May 17, 2018

**Last Modified:** May 17, 2018

**Protocol Integer ID:** 12250

**Keywords:** flick test, tail

## Troubleshooting

- 1 Mice are allowed to habituate for 30 min to the testing environment and the injection occur immediately the start of the test.
- 2 Each mouse is lightly restrained in a denim pocket while the distal half of it's tail is exposed into a thermostatically controlled temperature
- 3 Latency to respond to the heat stimulus by vigorous flexion of the tail is measured with a stopwatch.
- 4 Each mouse is given 3-5 trials with an inter-trial interval of 10 s, and a maximum trial duration of 8 s.
- 5 To prevent tissue injury each mouse's tail is promptly removed from the hot water after 30 s.
- 6 Latency performance of the last 3 trials are recorded.