

Jan 28, 2026

## Target enrichment of 10X Genomics scRNA-seq gene expression libraries

DOI

<https://dx.doi.org/10.17504/protocols.io.5qpvo1qpbg4o/v1>

Boxun Li<sup>1</sup>, Charles Gersbach<sup>1</sup>

<sup>1</sup>Duke University

Gersbach Lab



Andrea R Daniel

Duke University

### Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



**DOI:** <https://dx.doi.org/10.17504/protocols.io.5qpvo1qpbg4o/v1>

**Protocol Citation:** Boxun Li, Charles Gersbach 2026. Target enrichment of 10X Genomics scRNA-seq gene expression libraries. protocols.io <https://dx.doi.org/10.17504/protocols.io.5qpvo1qpbg4o/v1>

**Manuscript citation:**



**License:** This is an open access protocol distributed under the terms of the **[Creative Commons Attribution License](#)**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

**Protocol status:** Working

**We use this protocol and it's working**

**Created:** January 26, 2026

**Last Modified:** January 28, 2026

**Protocol Integer ID:** 241588

**Keywords:** 10X Genomics, scRNA-seq, neuron, astrocyte, target enrichment, target enrichment of 10x genomics scrna, enrichment of 10x genomics scrna, seq gene expression libraries of neuron mono, 10x genomics scrna, seq gene expression library, seq gene expression libraries this protocol, 10x genomic, mouse primary astrocyte, primary astrocyte, neuron mono

**Funders Acknowledgements:**

NIH

Grant ID: HG012053

## Disclaimer

This protocol is adapted from the work of Boxun Li in the Gersbach lab at the Duke University.

## Abstract

This protocol describes target enrichment of 10X Genomics scRNA-seq gene expression libraries of neuron mono-culture or co-culture with mouse primary astrocytes.

## Target enrichment of 10X Genomics scRNA-seq gene expression libraries

- 1 Enrichment DNA probe panel for target enrichment for scRNA-seq gene expression libraries using a hybridization-based method (Twist Bioscience, PN 101001, 104445, 100856, and 101262) was designed by Twist Bioscience for the 5' UTR and first ~400bp of the coding sequence of all transcripts annotated in Gencode v41 of the 222 transcription factor genes of interest along with 4 positive control genes (TFRC, UBQLN2, GRN, and CDH2) as well as 7 housekeeping genes (DPM1, KRIT1, BAD, ARF5, POLDIP2, HCCS, TSR3).
- 2 Target enrichment is performed on the 10X Genomics 5' HT v2 gene expression sequencing libraries of the putative regulatory element (pRE) Perturb-seq screens, according to the manufacturer's recommended protocol Standard Hybridization v2 Protocol (Twist Bioscience, DOC-001273 REV 3.0)
- 3 The enriched libraries were amplified using the 10X Genomics Amp Mix (10X Genomics, PN 2000047) and purified using Twist Purification Beads.
- 4 The target-enriched gene expression libraries were sequenced using Illumina Nextseq 2000, NovaSeq 6000 or NovaSeq X Plus sequencing to target depth of  $\sim 5 \times 10^3$  reads/cell.