

Sep 19, 2019

Testing selective agents for the Ichthyosporeans *Abeoforma whisleri*, *Pirum gemata* and *Spaheroforma artica*

DOI

dx.doi.org/10.17504/protocols.io.z5nf85e

Elena Casacuberta¹, Cristina Aresté²

¹Institute of Evolutionary Biology; ²IIBB

Protist Research to Opti...

Multicellgenomelab

1 more workspace



Elena Casacuberta

Institute of Evolutionary Biology

OPEN  ACCESS



DOI: dx.doi.org/10.17504/protocols.io.z5nf85e

Protocol Citation: Elena Casacuberta, Cristina Aresté 2019. Testing selective agents for the Ichthyosporeans *Abeoforma whisleri*, *Pirum gemata* and *Spaheroforma artica*. **protocols.io** <https://dx.doi.org/10.17504/protocols.io.z5nf85e>

Manuscript citation:

License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: April 17, 2019

Last Modified: September 19, 2019

Protocol Integer ID: 22414


Keywords: Selective agents, *Abeoforma whisleri*

Abstract

Minimal protocol for testing selective agents for *Abeoforma whisleri*, *Sphaeroforma arctica*, *Pirum gemmata*

Materials

MATERIALS

 Benomyl

 Phleomycin

 100 mg Nourseothricin Sulfate **G-Biosciences Catalog #RC-188**

 1 g Puromycin Dihydrochloride **G-Biosciences Catalog #RC-270**

 Marine Broth 2216 **Becton Dickinson (BD) Catalog #279110**

Carboxin, Fluorotic Acid

- 1 Seed cells on 24 well/plate. Density depends on the organism but should allow to screen for growth and death easily
Seed the same density to test four or five different dilutions of each drug
Seed enough wells to have 3X replicates for each condition
Seeded cells for *Abeoforma whisleri*, *Pirum gemmata* and *Sphaeroforma artica* 100 cells/well 500 microliters of medium.
- 2 Add to each well the corresponding amount of drug. Following the amounts tested in our case for *Abeoforma whisleri*, *Pirum gemmata*, *Sphaeroforma artica*

Nourseothricin	10-50 µgr/ml
Phleomycin	10-300 µgr/ml
Fluorotic Acid	250-30ugr/ml
Benomyl Carboxine Puromycin	1-20 µgr/ml 20-300µgr/ml 100-500 µgr/ml

- 3 Spin down cells every 24 hours and refresh media with new antibiotic.
Observe for up to a week.
If effects on growth are suspected repeat experiment with the appropriate concentration of drug and count cells every 24 hours.
- 4 Summary of our results

<i>Abeoforma whisleri</i>	Nourseothricin	10-50 µgr/ml (not sensitive)
	Phleomycin	10-300 µgr/ml (not sensitive)
	Fluorotic Acid	250-30 µgr/ml (not sensitive)
	Benomyl Carboxine P uromycin	1-20 (500) µgr/ml 20-300 µgr/ml 100-500 µgr/ml (cytostatic)
<i>Pirum gemmata</i>	Nourseothricin	10-50 µgr/ml (not sensitive)
	Phleomycin	10-300 µgr/ml
	Fluorotic Acid	250-30ugr/ml (not sensitive)
	Benomyl Carboxine P uromycin	1-20 µgr/ml 20-300µgr/ml 100-500 µgr/ml (cytostatic)



	<i>Sphaeroforma arctica</i>	Nourseothricin	10-50 µgr/ml (not sensitive)
		Phleomycin	10-300 µgr/ml
		Fluorotic Acid	250-30 µgr/ml (not sensitive)
		Benomyl	1-20 µgr/ml (cytostatic)
		Carboxine	20-300 µgr/ml
		Puromycin	100-500 µgr/ml (cytostatic)