

Apr 21, 2019

The Illumina libraries preparation for the Scapharca broughtonii

 [GigaScience](#)

 In 1 collection

DOI

<https://dx.doi.org/10.17504/protocols.io.z75f9q6>

Chang-Ming Bai¹

¹Yellow Sea Fisheries Research Institute, Chinese Academy of Fishery Sciences

GigaScience Press



Chang-Ming CM Bai

Yellow Sea Fisheries Research Institute, Chinese Academy of ...



Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.z75f9q6>

External link: <https://doi.org/10.1093/gigascience/giz067>



Protocol Citation: Chang-Ming Bai 2019. The Illumina libraries preparation for the *Scapharca broughtonii*. **protocols.io**
<https://dx.doi.org/10.17504/protocols.io.z75f9q6>

Manuscript citation:

Chang-Ming Bai, Lu-Sheng Xin, Umberto Rosani, Biao Wu, Qing-Chen Wang, Xiao-Ke Duan, Zhi-Hong Liu, Chong-Ming Wang, Chromosomal-level assembly of the blood clam, *Scapharca (Anadara) broughtonii*, using long sequence reads and Hi-C, *GigaScience*, Volume 8, Issue 7, July 2019, giz067, <https://doi.org/10.1093/gigascience/giz067>

License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: April 21, 2019

Last Modified: April 21, 2019

Protocol Integer ID: 22493

Keywords: preparation for the scapharca broughtonii genome, scapharca broughtonii genome, preparation for the scapharca broughtonii, scapharca broughtonii, genome, process of illumina, illumina, preparation

Abstract

This protocol is used to detail the process of Illumina libraries preparation for the *Scapharca broughtonii* genome.

- 1 The extracted DNA was sheared into fragments about 350 bp in size using a Bioruptor Pico Sonication System (Diagenode, Seraing, Belgium), and verified by agarose gel electrophoresis.

Note

The sizes of the main fragments should be between 200 bp and 500 bp.

- 2 Fragments with size > 300 bp were purified using VAHTSTM DNA Clean Beads (Vazyme Biotech Co., Ltd, Nanjing, China).
- 3 Repaired using NEBNext® End Prep Enzyme Mix (NEB, E6091) and NEBNext® End Repair Reaction Buffer (NEB, B6052) kept at 20°C, 30min; 65°C, 30min with the end-repaired fragments to obtain blunt ends which were then 3'-adenylated to create sticky ends.
- 4 These DNA fragments were ligated at both ends to T-tailed adapters and amplified.
- 5 PCR was performed using NEBNext® Ultra™ II Q5® Master Mix (NEB, M0544), and the Index i7 and Universal i5 primers. The temperature profile was 30 sec. at 98 °C followed by 10 cycles of 10 sec. at 98 °C, 75 sec. at 65 °C, and more 5 min. at 65 °C for further elongation.

Note

INDEX i7 PRIMER: 5'-CAA GCA GAA GAC GGC ATA CGA GAT ATGACGTC GTG ACT GGA GTT CAG ACG TGT GCT CTT CCG ATC*T-3'

NEBNext Universal i5 PCR Primer : 5'-AAT GAT ACG GCG ACC ACC GAG ATC TAC ACT CTT TCC CTA CAC GAC GCT CTT CCG ATC*T-3'

- 6 VAHTSTM DNA Clean Beads (Vazyme Biotech Co., Ltd, Nanjing, China) was used to purify the PCR production.