



May 14, 2018

Thermocoagulation induced focal ischemic model

 [PLOS One](#)

DOI

<https://dx.doi.org/10.17504/protocols.io.pasdiee>

Silva HR¹, Nucci MP², Mamani JB¹, Mendez-Otero R³, Nucci LP⁴, Tannus A⁵, Gamarra LF¹

¹Hospital Israelita Albert Einstein, São Paulo, Brazil;

²Hospital das Clínicas HCFMUSP, Universidade de São Paulo, São Paulo, SP – Brazil;

³Instituto de Biofísica Carlos Chagas Filho, Universidade Federal do Rio de Janeiro, Rio de Janeiro, Brazil;

⁴Universidade Federal de São Paulo, São Paulo, Brazil;

⁵CIERMag-Instituto de Física de São Carlos, Universidade de São Paulo, São Paulo, Brazil

Nanobiotechnology



Lionel Gamarra

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

[Create free account](#)

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.pasdiee>

External link: <https://doi.org/10.1371/journal.pone.0200135>

Protocol Citation: Silva HR, Nucci MP, Mamani JB, Mendez-Otero R, Nucci LP, Tannus A, Gamarra LF 2018. Thermocoagulation induced focal ischemic model. protocols.io <https://dx.doi.org/10.17504/protocols.io.pasdiee>

Manuscript citation:

Silva Hd, Nucci MP, Mamani JB, Mendez-Otero R, Nucci LP, Tannus A, Gamarra LF (2018) Evaluation of temperature induction in focal ischemic thermocoagulation model. PLoS ONE 13(7): e0200135. doi: [10.1371/journal.pone.0200135](https://doi.org/10.1371/journal.pone.0200135)

License: This is an open access protocol distributed under the terms of the [Creative Commons Attribution License](#), which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: April 07, 2018

Last Modified: May 14, 2018

Protocol Integer ID: 11314

Keywords: thermocoagulation, stroke, real-time blood perfusion, MRI, thermal dissipation, stroke model, permanent ischemic stroke model, focal ischemic model the thermocoagulation model, focal cerebral ischemia, type of focal cerebral ischemia, thermocoagulation model, focal ischemic model, permanent ischemic brain lesion, permanent ischemic brain lesions with good reproducibility, craniectomy

Abstract

The thermocoagulation model is a type of focal cerebral ischemia with craniectomy. It represents an opportunity to study permanent ischemic brain lesions with good reproducibility and low mortality

Attachments



Protocol-
Thermocoagu...

12KB

- 1 A focal ischemic lesion is induced by thermocoagulation of blood in the pial blood vessels of primary motor and somatosensory cortex
- 2 The animals are anesthetized with ketamine hydrochloride (90 mg/kg, i.p.) and xylazine hydrochloride (12 mg/kg, i.p.) and place in a stereotaxic apparatus
- 3 A craniectomy is made to expose the left somatosensory cortex, region of the left middle cerebral artery (+ 2 mm to -6 mm in anterior-posterior and +2 mm in medial-lateral axis from Bregma), according to the atlas of Paxinos and Watson.
- 4 Basal blood perfusion image is acquired in the exposed brain area with a Perfusion Speckle Imager before induction.
- 5 Thermocoagulation is performed in three different animal groups at temperatures of 200 °C, 300 °C, or 400°C, respectively, providing a better dynamic evaluation of the lesion in the pial blood vessels of the somatosensory cortex according to the temperature used.
- 6 Superficial blood vessels of the left sensorimotor brain area are thermocoagulated transdurally by a hot probe into the dura matter (~2 mm) for 30 minutes.
- 7 After 30 min of TCI at 400 °C, vessel lesions are evaluated macroscopically through the color changes in the region targeted, from light red to dark red to indicate complete thermocoagulation of blood. Thus TCI is performed for 30 min at the three temperatures. A decrease in blood perfusion by 75% of the initial measure (before induction) after 30 min of TCI is considered as a parameter for complete ischemic lesion.
- 8 The procedure is completed with suturing of the incision tissue and topical application of lidocaine and IV administration of tramadol (5 mg/kg) every 12 h for six days.
- 9 Throughout anesthesia, the rats are placed on a heating pad to maintain the rectal temperature at $37.0\text{ }^{\circ}\text{C} \pm 0.5\text{ }^{\circ}\text{C}$.
- 10 Animals from the sham group are underwent the whole procedure up to, but not including the TCI. The probe is positioned turned off for 30 min and thereafter, the incision tissue is sutured.