



Jul 11, 2024

Tile Scan Imaging and Cell Counting

DOI

<https://dx.doi.org/10.17504/protocols.io.bp2l6231kgqe/v1>

Shiyi Wang¹

¹Duke University



Shiyi Wang

Duke University

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.bp2l6231kgqe/v1>

Protocol Citation: Shiyi Wang 2024. Tile Scan Imaging and Cell Counting. **protocols.io**
<https://dx.doi.org/10.17504/protocols.io.bp2l6231kgqe/v1>

License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: July 11, 2024

Last Modified: July 11, 2024

Protocol Integer ID: 103200

Keywords: ASAPCRN, cell counting tile scan imaging, cell counting, tile, imaging

Funders Acknowledgements:

Aligning Science Across Parkinson's (ASAP) initiative

Grant ID: ASAP-020607

Disclaimer

DISCLAIMER – FOR INFORMATIONAL PURPOSES ONLY; USE AT YOUR OWN RISK

The protocol content here is for informational purposes only and does not constitute legal, medical, clinical, or safety advice, or otherwise; content added to **protocols.io** is not peer reviewed and may not have undergone a formal approval of any kind. Information presented in this protocol should not substitute for independent professional judgment, advice, diagnosis, or treatment. Any action you take or refrain from taking using or relying upon the information presented here is strictly at your own risk. You agree that neither the Company nor any of the authors, contributors, administrators, or anyone else associated with **protocols.io**, can be held responsible for your use of the information contained in or linked to this protocol or any of our Sites/Apps and Services.

Abstract

Tile Scan Imaging and Cell Counting

1 ****Sample Preparation****

- 1.1 - Prepare 30 μm thick coronal sections containing the anterior cingulate cortex (ACC) and primary motor cortex (MOp) from P21 WT and LRRK2 G2019Ski/ki; Aldh1l1-eGFP mice.

2 ****Confocal Microscopy****

- 2.1 - Acquire tile scan images using a confocal Leica SP8 STED microscope equipped with a galvo scanner and a 20x objective.
- 2.2 - Ensure consistent settings for image acquisition across all samples.

3 ****Double Positive Cell Identification****

- 3.1 - Identify cells labeled with ALDH1L1-EGFP (astrocyte marker) and SOX9 (astrocyte marker) in the ACC and MOp regions.
- 3.2 - Use FIJI software with the cell counter tool for manual cell counting.

4 ****Data Collection****

- 4.1 - Count ALDH1L1-EGFP and SOX9 double positive cells manually in 2-4 sections per brain.
- 4.2 - Analyze sections from 3 sex- and age-matched mice per genotype.