

Jun 13, 2023

Tissue and blood collection

DOI

<https://dx.doi.org/10.17504/protocols.io.bp2l69wmklqe/v1>

michela.deleidi^{1,2}, Bianca Marchetti^{3,4}, Federico Bertoli^{1,2}, Carmela Giachino³

¹Mitochondria and Inflammation in Neurodegenerative Diseases, DZNE, Tübingen-Germany;

²Hertie Institute for Clinical Brain Research, University of Tübingen;

³Neuropharmacology Laboratory, Oasi Research Institute-IRCCS, Troina, Italy;

⁴Biomedical and Biotechnological Sciences, Pharmacology Section, University of Catania-Italy



carmela giachino

institut imagine

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.bp2l69wmklqe/v1>

Protocol Citation: michela.deleidi , Bianca Marchetti, Federico Bertoli, Carmela Giachino 2023. Tissue and blood collection. protocols.io <https://dx.doi.org/10.17504/protocols.io.bp2l69wmklqe/v1>

License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited



Protocol status: Working

We use this protocol and it's working

Created: March 31, 2023

Last Modified: June 13, 2023

Protocol Integer ID: 79800

Keywords: Tissue collection, tissue collection, collection from mice, blood collection this protocol detail, blood collection, tissue, mice, blood, collection, protocol, protocol detail

Funders Acknowledgements:

ASAP-Aligning Science Across Parkinson's

Grant ID: ASAP-000420

Abstract

This protocol details tissue collection from mice.

Attachments



[k4dmb48if.docx](#)

13KB

Materials

Materials

- 0.9% NaCl
- 4% paraformaldehyde
- phosphate buffer
- EDTA-treated tubes
- 15% sucrose solution



Tissue and blood collection

- 1 For immunohistochemistry, mice (5-6 mice/genotype/age-group/treatment) were anesthetized and transcardially perfused with 0.9% NaCl, followed by 4 % paraformaldehyde in phosphate buffer (pH 7.2 at 4°C), the brains and peripheral tissues carefully removed and post-fixed for 2-4 hrs, in 4% paraformaldehyde in phosphate buffer (pH 7.2) and placed in 15% sucrose in the solution of phosphate buffer overnight at 4°C.
- 2 Blood was collected into EDTA-treated tubes and serum was separated by centrifugation and stored at -20C.
- 3 For gene expression and protein analysis, animals were sacrificed by cervical dislocation, the brains and peripheral tissues were quickly isolated and then stored at -80°C until assayed, the tails collected from each mouse were genotyped through PCR followed by electrophoretic analysis, as described in the manufacturer's instructions (Jackson Laboratory).

