

May 09, 2020

Version 4

TissueCyte Specimen Embedding Acrylamide Coembedding V.4

DOI

<https://dx.doi.org/10.17504/protocols.io.bf65jrg6>

Allen Institute for Brain Science¹

¹Allen Institute

BICCN / BICAN

Allen Institute for Brain S...



Allen Institute

Allen Institute

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.bf65jrg6>

Protocol Citation: Allen Institute for Brain Science 2020. TissueCyte Specimen Embedding Acrylamide Coembedding. protocols.io <https://dx.doi.org/10.17504/protocols.io.bf65jrg6>

License: This is an open access protocol distributed under the terms of the [Creative Commons Attribution License](#), which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: May 09, 2020

Last Modified: March 17, 2021

Protocol Integer ID: 36797

Keywords: Embedding, Coembedding, TissueCyte, Co-Embedding, PF0315, agarose block for blockface imaging, fixed mouse brain specimen, embedding acrylamide, mouse brain specimen, tissuecyte specimen, blockface imaging, agarose block, tissuecyte 1000 system, tissuecyte, imaging,

Abstract

This protocol describes the preparation of a fixed mouse brain specimen bonded to an agarose block for blockface imaging and sectioning using the TissueCyte 1000 system.

Note: Research reported in this publication was supported by the National Institute Of Mental Health of the National Institutes of Health under Award Number U19MH114830. The content is solely the responsibility of the authors and does not necessarily represent the official views of the National Institutes of Health.

Attachments



PF0315_TissueCyte_Sp..

↓

321KB

