



Feb 10, 2026

Toxicity Testing of Compounds in *Galleria mellonella* Larvae

DOI

<https://dx.doi.org/10.17504/protocols.io.ewov1k7wkgr2/v1>

Juan Jose Quispe Haro¹

¹University of Helsinki

Galleria protocols



Juan Jose Quispe Haro

University of Helsinki

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.ewov1k7wkgr2/v1>

Protocol Citation: Juan Jose Quispe Haro 2026. Toxicity Testing of Compounds in *Galleria mellonella* Larvae. **protocols.io**
<https://dx.doi.org/10.17504/protocols.io.ewov1k7wkgr2/v1>

License: This is an open access protocol distributed under the terms of the **[Creative Commons Attribution License](#)**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: In development

We are still developing and optimizing this protocol

Created: February 10, 2026

Last Modified: February 10, 2026

Protocol Integer ID: 242969

Keywords: Galleria mellonella, toxicity testing, invertebrate model, drug screening, larval injection, acute toxicity, survival assay, galleria mellonella larvae this protocol, galleria mellonella larvae, using galleria mellonella, toxicity testing of compound, preliminary toxicity screening, method for toxicity screening, cost preliminary toxicity screening in an invertebrate model, toxicity screening, galleria mellonella, larvae per treatment group, larvae, test compound, microbial metabolite, assay, 7th instar larvae, test compounds via the last proleg, chemical compound, compound, nanoparticle

Disclaimer

DISCLAIMER – FOR INFORMATIONAL PURPOSES ONLY; USE AT YOUR OWN RISK

The protocol content here is for informational purposes only and does not constitute legal, medical, clinical, or safety advice, or otherwise; content added to protocols.io is not peer reviewed and may not have undergone a formal approval of any kind. Information presented in this protocol should not substitute for independent professional judgment, advice, diagnosis, or treatment. Any action you take or refrain from taking using or relying upon the information presented here is strictly at your own risk. You agree that neither the Company nor any of the authors, contributors, administrators, or anyone else associated with protocols.io, can be held responsible for your use of the information contained in or linked to this protocol or any of our Sites/Apps and Services.

Abstract

This protocol describes a method for toxicity screening of chemical compounds, nanoparticles, or microbial metabolites using *Galleria mellonella* larvae. The procedure utilizes 7th instar larvae injected with test compounds via the last proleg, followed by survival monitoring over 72 hours. With a sample size of 10+ larvae per treatment group, this assay allows for statistically strong, rapid, low-cost preliminary toxicity screening in an invertebrate model that shares key physiological similarities with mammals.

Guidelines

- **Dose volume:** Do not exceed 10–20 µl per larva to avoid mechanical damage.
- **Vehicle toxicity:** Always include vehicle control; DMSO >10% can be toxic.
- **Larvae quality:** Use only healthy, active larvae of similar size.
- **Sterility:** Work aseptically to avoid infection-related death.
- **Replicates:** Repeat experiment at least 3 times independently for statistical power (total n=30 per group across replicates).



Materials

Galleria mellonella larvae: 7th instar larvae (≈ 300 mg each, cream-colored, active). Minimum 10 larvae per treatment group + 10% extra for selection.

Hamilton syringe (10 μ l) with removable needle.

6-well cell culture plates (sterile, non-treated).

Optional: Dissecting microscope or magnifying lamp.

Safety warnings

- ⚠ Follow institutional guidelines for invertebrate use. Euthanize larvae by freezing at -20°C for 2 h after experiment.

Before starting

- 1 Acclimate larvae at experimental temperature for 24 h before injection.
 30 °C or experimental temp.
- 2 Prepare compound solutions freshly on the day of injection. Filter-sterilize if possible (0.22 µm).
- 3 Plan groups (n=10-16 larvae per group):

Group 0: Optional handling control (to test if larvae survive handling/physical trauma).

Group 1: Vehicle control (same solvent used to prepare the compound solutions).

Group 2: Positive control (substance known to be toxic with similar mode of action of test compound)

Groups 3-X: Test compounds at desired concentrations (e.g., low, medium, high dose, or serial dilutions).

- 4 Assign random 7th instar larvae to each group to avoid health/size bias.

Injection Procedure

- 5 Surface sterilize each larvae before injection by wiping each larva gently with 70% ethanol and allowing to dry on filter paper.
- 6 Prepare a Hamilton syringe with 10 µl of the desired solution.
- 7 Grab a larva by its sides with the thumb and the index fingers and flip it upside down to expose its abdomen.
- 8 Localize the hind-left proleg of the larva and carefully insert the needle of the syringe about 5 mm into the body of the larva.
- 9 Inject the contents of the syringe slowly to prevent any leaking, and remove the syringe.



- 10 Place each individual larva in a well of a 6 or 12-well dish and repeat for all the larvae and conditions.

Incubation and Monitoring

- 11 Transfer the dishes to an incubator and incubate the larvae for 72+ hours at desired temperature with 12-12 hours illumination cycle. Do not provide food during experiment.
- 12 Monitor survival at 0, 2, 4, 6, 8, 12, 24, 48, 72 hours post-injection (adjust based on kinetics).
- 13 Monitor health using the Health Score Index points system every 24 hours.
Modified from DOI: 10.1080/21505594.2021.1950269

	A	B	C
	Category	Description	Score
	Activity	no movement	0
		movement only after stimulation	1
		letargic movement	2
		whole body movement	3
	Cocoon formation	no cocoon	0
		partial cocoon	1
		full cocoon	2
	Melanization	black larvae	0
		black spots on brown larvae	1
		≥3 spots on beige larvae	2



	A	B	C
		<3 spots on beige larvae	3
		no melanization	4

Data analysis

- 14 Calculate percent survival per group over time and plot Kaplan–Meier survival curves. This can be done easily in GraphPad Prism.
- 15 Compare groups using log-rank test or Cox proportional hazards (if multiple doses).
- 16 Plot the average health score vs time. Analyse the data by performing a Repeated Measures ANOVA, and post-hoc tests (Bonferroni or Tukey's HSD) to determine if the differences are significant.
- 17 Determine the LD_{50} using the following formula:
$$LD_{50} = \text{Log}_{10}(x_k) + (\text{Log}_{10}(d) * (0.5 - \sum p_i))$$
where:
 x_k = Lowest concentration that produces 100% mortality
 d = Log-interval between doses (10-fold dilutions)
 $\sum p_i$ = Sum of the proportions of mortality for all doses up to but not including x_k

Protocol references

Larvae Health Index Scoring System, modified from **Establishing an invertebrate *Galleria mellonella* greater wax moth larval model of *Neisseria gonorrhoeae* infection**, *Virulence*, 2021 Jul 25;12(1):1900–1920. doi: [10.1080/21505594.2021.1950269](https://doi.org/10.1080/21505594.2021.1950269)