Feb 13, 2019 Version 1

# Tranfection of sgRNA using SpCas9 containing plasmids to generate cell lines with Cas9 V.1

DOI

dx.doi.org/10.17504/protocols.io.x37fqrn

Binnypreet Kaur<sup>1</sup>, 2<sup>2</sup>, Drahomíra Faktorová<sup>1</sup>, 2<sup>2</sup>, and Julius Lukeš<sup>1</sup>, <sup>2</sup>, <sup>3</sup>

<sup>1</sup>Institute of Parasitology, Czech Academy of Sciences;

<sup>2</sup>Faculty of Sciences, University of South Bohemia, České Budějovice, Czech Republic; <sup>3</sup>Canadian Institute for Advanced Research, Toronto, Canada

#### Binnypreet Kaur

1 Institute of Parasitology, Biology Centre, Czech Academy o...





#### DOI: dx.doi.org/10.17504/protocols.io.x37fqrn

**Protocol Citation:** Binnypreet Kaur, 2, Drahomíra Faktorová, 2, and Julius Lukeš, , 2019. Tranfection of sgRNA using SpCas9 containing plasmids to generate cell lines with Cas9. **protocols.io** <u>https://dx.doi.org/10.17504/protocols.io.x37fqrn</u>

License: This is an open access protocol distributed under the terms of the <u>Creative Commons Attribution License</u>, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: In development We are still developing and optimizing this protocol

Created: February 13, 2019

Last Modified: February 13, 2019

Protocol Integer ID: 20319

## Abstract

This protocol describes transfection of ribonucleoprotein (RNP) complexes that consist of purified Cas9 nuclease duplexed with synthetic guide RNA in cultured cells using the AMAXA Nucleo AMAXA Nucleofector II Electroporation system using Program X-001.

RNPs complex have only transient expression inside the cell, this allows for the highest levels of editing efficiency and greatly reduces the chances of possible off target and toxic effects of integration of foreignDNA in host genome

### Guidelines

#### **Pre-Electroporation**

Subculture cells 2-3 days before electroporation and seed cells in the appropriately sized vessel so that they are 70-80% confluent on the day of transfection. Each transfection reaction will require approximately  $5 \times 10^7$  cells.

## Materials

- 1. Synthetic guide RNA (sgRNA or crRNA:tracrRNA)
- 2. 2NLS-Cas9 nuclease
- 3. AMAXA Nucleoporator
- 4. AMAXA T- cell solution
- 5. AMAXA T- supplement
- 6. Cell counter
- 7. Normal growth medium
- 8. Tissue culture plates

- 1 Step1: Assemble RNP complexes with (30pmol/µl) of gRNA and (20pmol/µl) of Cas9 in the ratio of (Cas9:gRNA) 1:9 in resuspension buffer before electroporation and keep it at room temperature for 10 min and then transfer it to ice untill use.
- 2 Step2: Follow the protocol 7 for transfection