

Feb 13, 2019

Version 3

Tranfection of sgRNA using SpCas9 containing plasmids to generate cell lines with Cas9 V.3

DOI

<https://dx.doi.org/10.17504/protocols.io.x39fqr6>

Binnypreet Kaur¹, Drahomíra aktorová¹, Julius Lukeš^{2, 3, 4}

¹Institute of Parasitology, Czech Academy of Sciences and Faculty of Sciences, University of South Bohemia, České Budějovice, Czech Republic;

²Institute of Parasitology, Czech Academy of Sciences;

³Faculty of Sciences, University of South Bohemia, České Budějovice, Czech Republic;

⁴Canadian Institute for Advanced Research, Toronto, Canada

Protist Research to Opti...

Julius Lukes



Binnypreet Kaur

1 Institute of Parasitology, Biology Centre, Czech Academy o...

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.x39fqr6>



Protocol Citation: Binnypreet Kaur, Drahomíra aktorová, Julius Lukeš, , 2019. Transfection of sgRNA using SpCas9 containing plasmids to generate cell lines with Cas9. **protocols.io** <https://dx.doi.org/10.17504/protocols.io.x39fqr6>

License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: In development

We are still developing and optimizing this protocol

Created: February 13, 2019

Last Modified: February 13, 2019

Protocol Integer ID: 20321

Keywords: synthetic guide rna in cultured cell, synthetic guide rna, purified cas9 nuclease, foreigndna in host genome, cell lines with cas9, using spcas9, host genome, ribonucleoprotein, genome, transfection, foreigndna, containing plasmid, spcas9, cas9, amaxa nucleo amaxa nucleofector ii electroporation system, cultured cell, rnps complex

Abstract

This protocol describes transfection of ribonucleoprotein (RNP) complexes that consist of purified Cas9 nuclease duplexed with synthetic guide RNA in cultured cells using the AMAXA Nucleo AMAXA Nucleofector II Electroporation system using Program X-001.

RNPs complex have only transient expression inside the cell, this allows for the highest levels of editing efficiency and greatly reduces the chances of possible off target and toxic effects of integration of foreignDNA in host genome

Guidelines

Pre-Electroporation

Subculture cells 2-3 days before electroporation and seed cells in the appropriately sized vessel so that they are 70-80% confluent on the day of transfection. Each transfection reaction will require approximately 5×10^7 cells.

Materials

1. Synthetic guide RNA (sgRNA or crRNA:tracrRNA)
2. 2NLS-Cas9 nuclease
3. AMAXA Nucleoprotector
4. AMAXA T- cell solution
5. AMAXA T- supplement
6. Cell counter
7. Normal growth medium
8. Tissue culture plates



- 1 Step1: Assemble RNP complexes with (30pmol/ μ l) of gRNA and (20pmol/ μ l) of Cas9 in the ratio of (Cas9:gRNA) 1:9 in resuspension buffer before electroporation and keep it at room temperature for 10 min and then transfer it to ice until use.
- 2 Step2: Follow the protocol 7 for transfection