

Aug 30, 2018

Transformation of *Acanthamoeba castellanii* with Qiagen SuperFect reagent

DOI

<https://dx.doi.org/10.17504/protocols.io.s4regv6>

Morgan Colp¹

¹Dalhousie University

Protist Research to Opti...



Morgan Colp

Dalhousie University

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.s4regv6>

External link: <https://www.sciencedirect.com/science/article/pii/S0167488904002095>

Protocol Citation: Morgan Colp 2018. Transformation of *Acanthamoeba castellanii* with Qiagen SuperFect reagent. **protocols.io** <https://dx.doi.org/10.17504/protocols.io.s4regv6>

License: This is an open access protocol distributed under the terms of the [Creative Commons Attribution License](#), which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited



Protocol status: Working

We use this protocol and it's working

Created: August 30, 2018

Last Modified: August 30, 2018

Protocol Integer ID: 15217

Keywords: transformation of *acanthamoeba castellanii*, plasmid dna, *acanthamoeba castellanii*, qiagen superfect reagent transformation, qiagen superfect, dna, plasmid, reagent transformation

Abstract

Transformation of *Acanthamoeba castellanii* Neff (ATCC-30010) with plasmid DNA can be achieved with a modified version of the protocol described in Peng, Omaruddin, & Bateman (2005) using constructs from Bateman (2010).

Before start

Materials needed:

- PYG medium with additives (0.75% yeast extract, 0.75% proteose peptone, 2 mM KH_2PO_4 , 1 mM MgSO_4 , 1.5% glucose, 0.1 mM ferric citrate, 0.05 mM CaCl_2 , 1 $\mu\text{g/mL}$ thiamine, 0.2 $\mu\text{g/mL}$ D-biotin, and 1 ng/mL vitamin B12.
- Plasmid DNA
- Qiagen SuperFect
- Encystment medium (20 mM Tris-HCl [pH 8.8], 100 mM KCl, 8 mM MgSO_4 , 0.4 mM CaCl_2 , 1 mM NaHCO_3)

Strain and Culture Conditions

- 1 *A. castellanii* Neff (ATCC-30010) cells were grown axenically in PYG medium at room temperature in 75 cm³ vented tissue culture flasks with no shaking.

Plasmids

- 2 **pGAPDH-EGFP:** This plasmid contains a bacterial ampicillin resistance gene for propagation, a neomycin resistance gene (*neoR*; an aminoglycoside phosphotransferase) is expressed from the TBP promoter, and an EGFP gene is expressed from the GAPDH promoter.

pTPBF-EGFP: This plasmid contains a bacterial ampicillin resistance gene for propagation, a neomycin resistance gene (*neoR*; an aminoglycoside phosphotransferase) is expressed from the TBP promoter, and an EGFP gene is expressed from the TPBF promoter.

These plasmids were provided by the Andrew Roger lab at Dalhousie University who in turn acquired them from the Erik Bateman lab at the University of Vermont.

Bateman, E. (2010). Expression plasmids and production of EGFP in stably transfected *Acanthamoeba*. *Protein Expression and Purification*, 70(1), 95. doi: 10.1016/j.pep.2009.10.008

Cell Preparation

- 3 Per reaction: 5×10^5 cells are pelleted by centrifugation at 1000 *g* for 5 minutes at 4 °C, and resuspended in 500 microlitres of encystment medium (20 mM Tris-HCl [pH 8.8], 100 mM KCl, 8 mM MgSO₄, 0.4 mM CaCl₂, 1 mM NaHCO₃).
- 4 The cell suspension is added to one well of a six well culture plate.

DNA Preparation

- 5 4 µg of plasmid DNA in 3 to 6 µL is diluted in 100µL of encystment medium.



- 6 20 μ L of Qiagen SuperFect reagent are added to the DNA solution desired for transfection (in this case pGAPDH-EGFP), and mixed by pipetting 5 times.
- 7 The DNA/reagent mixture is incubated for 10 minutes at room temperature.
- 8 600 μ L of encystment medium are added to the DNA/reagent mixture and mixed.

Transfection Reaction

- 9 The whole aliquot of DNA/reagent/medium mixture is added to the culture dish well that contains the cells.
- 10 The cells are incubated with DNA and the SuperFect reagent for 3 hours at room temperature.

Recovery and selection

- 11 The mixture can be pipetted off and replaced with 4 mL of growth (PYG) medium.
- 12 Prior to applying selection, the cells should be allowed to recover for 24 hours in their typical growth conditions.
- 13 Selection at ~20% of the final antibiotic concentration should be applied until growth is observed. This was 10 μ g/mL Geneticin (G418) for this experiment. It typically took about a week for sufficient growth to occur, and visually this corresponds to a cell density roughly half that of the peak density achieved during regular culturing. This can be verified with cell counts but visual confirmation seems to be fine.
- 14 Subsequent transfers can make use of the full concentration of antibiotic selection for indefinite maintenance, which in this case is 50 μ g/mL Geneticin (G418). If the cells are transferred diligently in 2-3 week intervals (or whichever interval involves transferring at peak density before a lot of encystation occurs), their growth should mirror that of wild-type cells even under continued selection, although sometimes they are slow to take off after a culture transfer.