



Mar 24, 2025

🌐 Transient Transfection of Cultured Mammalian Cells for Live 3D Fluorescence Microscopy

DOI

<https://dx.doi.org/10.17504/protocols.io.ewov1dr82vr2/v1>

Eve V. Kakudji¹, Samantha C Lewis¹

¹UC Berkeley



Samantha C Lewis

UC Berkeley

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.ewov1dr82vr2/v1>

Protocol Citation: Eve V. Kakudji, Samantha C Lewis 2025. Transient Transfection of Cultured Mammalian Cells for Live 3D Fluorescence Microscopy. **protocols.io** <https://dx.doi.org/10.17504/protocols.io.ewov1dr82vr2/v1>

License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working



Created: February 20, 2025

Last Modified: March 24, 2025

Protocol Integer ID: 123321

Keywords: Transfection, Fluorescence microscopy, Plasmids, mitochondria, live imaging, mitochondrial proteins into cultured mammalian cell, transient transfection of cultured mammalian cell, tagged mitochondrial protein, live 3d fluorescence microscopy, mitochondrial morphometric, cultured mammalian cell, various features of mitochondrial morphometric, fluorescence microscopy, sample of live cell, transient transfection, cell, plasmid, live cell, protocol for transient transfection

Funders Acknowledgements:

Michael J. Fox Foundation

Grant ID: 059455

Abstract

This is a protocol for transient transfection of plasmids encoding tagged mitochondrial proteins into cultured mammalian cells. The result is a sample of live cells in which various features of mitochondrial morphometrics can be observed by fluorescence microscopy.

Guidelines

Timeline

- **D0:** Seed U2OS cells (typically at ~ 20-30% confluency, so that they are at ~ 80% confluency on the day of the experiment).
- **D1:** Perform transfection – Replace with fresh media 5 hours post-transfection.
- **D2:** Perform experiment 24 hours post-transfection.



Materials

Reagents:

- Complete DMEM media

1. 500mL  DMEM high glucose no glutamine **Thermo Fisher Scientific Catalog #11960044**

2. 55mL

 Fetal Bovine Serum, heat inactivated, qualified, One Shot[®], United States **Thermo Fisher Catalog #A3840101**

3. 5.5 mL  L-Glutamine (200 mM) **Thermo Fisher Catalog #25030081**

4. 5.5mL  Penicillin-Streptomycin **Gibco - Thermo Fisher Scientific Catalog #15140122**

-  Opti-MEM[™] I Reduced Serum Medium **Thermo Fisher Scientific Catalog #31985070**

-  Lipofectamine[™] 2000 Transfection Reagent **Thermo Fisher Scientific Catalog #11668030**

- Plasmids

1.  mCherry-TOMM20-N-10 **addgene Catalog #55146**

2.  ATP5F1B (NM_001686) Human Tagged ORF Clone **OriGene Catalog #RG201638**

- 35-nm glass bottom dishes (Mattek, P35GC-1.5-14-C)

- Pipettes

- Pipette tips

- 1.5mL or 2mL microtubes

Equipment:

- Incubator

- Biosafety cabinet

- Fluorescence microscope



Procedure

1d 5h 15m

- 1 Seed U2OS cells in DMEM complete medium onto 35-nm glass bottom imaging dishes.
- 2 Incubate the cells at 37 °C , 5% CO₂ for ~ 24:00:00 . 1d
- 3 Thaw plasmid(s) and measure your desired concentration for transfection (500 ng for mCherry-TOMM20-N-10 and 250 ng for ATP5F1B-tGFP).
- 4 Take an aliquot of Opti-MEM and let it warm at Room temperature in the dark.
- 5 Take a microfuge tube: add 100 µL of Opti-MEM and 4 µL of lipofectamine 2000 per transfection (example: if you are doing two transfections, you will need 200 µL Opti-MEM and 8 µL lipofectamine 2000). Mix gently.
- 6 Take microfuge tubes. In each tube, add 100 µL of Opti-MEM and your desired concentration of plasmid (example: 500 ng for mCherry-TOMM20-N-10 / 250 ng for ATP5F1B-GFP). Mix gently.
- 7 Add 104 µL of the Opti-MEM/Lipofectamine 2000 mix in each of the tubes prepared in step 6. Mix gently.
- 8 Let the tubes incubate at Room temperature for 00:15:00 . 15m
- 9 Take the cells out of the incubator and add the transfection mixture dropwise to cells.
- 10 Incubate the cells at 37 °C , 5% CO₂ for 05:00:00 . 5h



- 11 Replace media with fresh DMEM complete media.
- 12 Perform experiment 24 hours post-transfection.

Acknowledgements

We thank the Lewis Lab for helpful discussion.