

Aug 05, 2020

Trypsin digestion

DOI

dx.doi.org/10.17504/protocols.io.bjdpki5n

avinash.kale¹

¹domnic colvin

UM-DAE Centre for Exce...



avinash.kale

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.bjdpki5n>

Document Citation: avinash.kale 2020. Trypsin digestion. **protocols.io** <https://dx.doi.org/10.17504/protocols.io.bjdpki5n>

License: This is an open access document distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Created: August 05, 2020

Last Modified: August 05, 2020

Document Integer ID: 40079

Troubleshooting



Trypsin digestion

In solution trypsin digestion of the TRIzol extracted protein sample was carried out by using Trypsin (Promega). Briefly, 15 μL of Digestion Buffer (50 mM Ammonium bicarbonate) and 1.5 μL of Reducing Buffer (100mM Dithiothreitol) were added to 10 μL of the protein solution to the micro-centrifuge tube and the final volume was adjusted to 27 μL with ultrapure water. The samples were incubated at 56 °C for 45 minutes and allowed to cool at room temperature. 3 μL of freshly prepared alkylation buffer (100mM Iodoacetamide) was then added to the tube and kept for incubation in the dark at room temperature for 20 minutes. Finally, 10 μL of Trypsin (20 $\mu\text{g}/\text{ml}$) solution was added in the tube and kept for incubation at 37°C for 15 hours (overnight)(25).