

Apr 05, 2024

Two-action sequence reinforcement

 Forked from [Single action sequence reinforcement](#)

DOI

<https://dx.doi.org/10.17504/protocols.io.3byl49248go5/v1>

Jonathan Tang^{1,2,3}

¹Department of Neuroscience, Zuckerman Mind Brain Behavior Institute, Columbia University, New York, NY 10027, USA;

²Seattle Children's Research Institute, Center for Integrative Brain Research, Seattle, WA, USA;

³University of Washington School of Medicine, Department of Pediatrics, Seattle, WA, USA



Christine Weber-Schmidt

Allen Institute

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.3byl49248go5/v1>

Protocol Citation: Jonathan Tang 2024. Two-action sequence reinforcement . protocols.io

<https://dx.doi.org/10.17504/protocols.io.3byl49248go5/v1>



Manuscript citation:

Tang, J.C.Y., Paixao, V., Carvalho, F. *et al.* Dynamic behaviour restructuring mediates dopamine-dependent credit assignment. *Nature* (2023). <https://doi.org/10.1038/s41586-023-06941-5>

License: This is an open access protocol distributed under the terms of the [Creative Commons Attribution License](#), which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: April 05, 2024

Last Modified: May 31, 2024

Protocol Integer ID: 97817

Keywords: ASAPCRN, mouse, behavior, brain, action sequence reinforcement protocol for mouse study, action sequence reinforcement protocol, action sequence reinforcement, mouse study, mouse, protocol

Funders Acknowledgements:

Life Sciences Research Fellowship

NINDS K99/R00 Award

Grant ID: 1K99NS112575

NIH

Grant ID: 5U19NS104649

Aligning Science Across Parkinson's (ASAP)

Grant ID: ASAP-020551

Disclaimer

The [protocols.io](#) team notes that research involving animals and humans must be conducted according to internationally-accepted standards and should always have prior approval from an Institutional Ethics Committee or Board.

Abstract

Two-action sequence reinforcement protocol for mouse studies from Tang et al 2023.

Guidelines

Individual mice were subjected to a single session of protocol each day, with sessions following each other on consecutive days.

Materials

The open-field box is made of 410 × 400 mm grey opaque acrylic walls and a 410 × 400 mm white matte acrylic base.



First reinforcement session+ Baseline

- 1 Mice were placed in a grey open-field behaviour recording for this protocol. To acquire baseline behavior, individual mice were allowed to behave freely inside the box for 30 minutes when the laser stimulation was not available for reinforcement.
- 2 After initial behavior acquisition, 60 minutes of laser availability was given. During the reinforcement periods, when the closed-loop system detects performance of the proximal action (T1) of interest, the algorithm enters a state in which the laser is triggered on performance of the distal action (T2), regardless of the amount of time that has elapsed between the latest T1 and T2.

Remaining Sessions

- 3 90 minutes of closed loop reinforcement were made available for individual mice during subsequent sessions. During the reinforcement periods, when the closed-loop system detects performance of the proximal action (T1) of interest, the algorithm enters a state in which the laser is triggered on performance of the distal action (T2), regardless of the amount of time that has elapsed between the latest T1 and T2.

Extinction protocol

- 4 An extinction protocol was carried out comprising of 20-minute maintenance of reinforced behavior with laser availability, followed by 60 minutes of extinction of reinforced behavior without laser availability, followed by 20-minute re-acquisition of reinforced behavior with laser availability.