



Sep 13, 2019

Version 2

UC Davis - Atheroma Quantification V.2

DOI

dx.doi.org/10.17504/protocols.io.7bwhipe



Amparo Villablanca¹

¹University of California, Davis

Mouse Metabolic Phenotyping Centers
Tech. support email: info@mmpc.org



Lili Liang

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.7bwhipe>

External link: <https://mmpc.org/shared/document.aspx?id=109&docType=Protocol>



Protocol Citation: Amparo Villablanca 2019. UC Davis - Atheroma Quantification. **protocols.io**

<https://dx.doi.org/10.17504/protocols.io.7bwhipe>

License: This is an open access protocol distributed under the terms of the **[Creative Commons Attribution License](#)**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: September 13, 2019

Last Modified: September 13, 2019

Protocol Integer ID: 27734

Keywords: Atheroma Quantification, neutral lipid accumulation, atherosclerotic lesion development, atherosclerotic cardiovascular disease, lipid accumulation in the artery wall, atheroma quantification summary, hallmarks of atherosclerotic cardiovascular disease, atheroma quantification, neutral lipid accumulation in the vessel wall, increased lipid accumulation, artery wall, aortic segment

Abstract

Summary:

Increased lipid accumulation in the artery wall and atherosclerotic lesion development are two hallmarks of atherosclerotic cardiovascular disease. Atheroma quantification and characterization assesses neutral lipid accumulation in the vessel wall, percentage of lesions per aortic segment (quantitative) and lesion morphology, complexity, and severity (qualitative).

Modified from: Villablanca AC, *et.al.* J Cardiovasc Transl Res. 2009 Sep;2(3):289-99.



Materials

MATERIALS

⊗ DPBS Invitrogen - Thermo Fisher Catalog #14190

⊗ formaldehyde Fisher Scientific Catalog #F79

⊗ Tissue-Tek OCT Fisher Scientific Catalog #14-373-65

⊗ Oil-red-O Merck MilliporeSigma (Sigma-Aldrich) Catalog #O0625

⊗ Gills hematoxylin Merck MilliporeSigma (Sigma-Aldrich)

Reagent Preparation:

Reagent 1: 4% & 10% paraformaldehyde

Formaldehyde (Fisher) is diluted to 4% or 10% in DPBS (Invitrogen)

Note:

Fisher Scientific, RRID:SCR_008452

Sigma-Aldrich, RRID:SCR_008988

Troubleshooting

Safety warnings



WARNING:

Formalin is, toxic, flammable and considered a carcinogen

All blood components and biological materials should be handled as potentially hazardous. Follow universal precautions established by CDC when handling and disposing of infectious agents.

- 1 Mice are anesthetized with an intraperitoneal injection with 50 mg pentobarbital/kg weight.
- 2 Immediately after sacrifice, aortas are perfused with phosphate-buffered saline to remove blood after venting and nicking the heart.
- 3 After careful removal of periadventitial fat, aortas are excised and fixed (4% followed by 10% paraformaldehyde overnight), and cut into three segments*
 - a. Proximal segment including all of the aortic arch
 - b. Midsegment consisting of the thoracic and suprarenal aorta
 - c. Distal segment consisting of the infrarenal aorta proximal to the iliac bifurcation.

* **Note:** aortas can also be analysed for a desired segment (for example: coronary sinus only, or proximal segment between aortic arch and the aortic valve only)
- 4 Aortic segments are placed side-by-side in cryomolds with optimum cutting temperature (OCT) compound frozen on dry ice in a cryomold, and then stored at -80°C.
- 5 Tissues are later sectioned with a microtome cryostat (HM 500 M) to 10 µm thick taken at -30°C, discarding every other section. (This allows sampling of all three segments of the aorta, permitting direct comparisons of each.)
 - a. Four sections per slide with a total of 9 slides
 - b. 36 sections per segment and 108 cryosections per aorta
 - c. Total of 360 µm from each segment and 1,000 µm from each aorta.

Note: Atheroma quantification can be completed on frozen tissue sections provided by the investigator

- 6 Quantitative morphometric analyses of aortic lesions is performed by staining cryosections with oil-red-O and Gill's hematoxylin
 - a. oil-red-O: define and characterize neutral lipid in the vessel wall
 - b. Gill's hematoxylin: provides histological assessment of lesions in artery segments
- 7 Sections are analyzed for lesion location and area by direct imaging (Olympus BX-40 microscope with a DP11 camera with smart media chip).
- 8 Atheromas are identified by a single observer blinded to experimental parameters, and mean lesion area ($\mu\text{m}^2 \pm \text{SE}$) determined for each atheroma by morphometric analysis with a computer-assisted imaging system (Image Pro Plus, version 4.1).

a. The quantification data is highly reproducible having an intraobserver error of ~6%.

- 9 Atheroma distribution is expressed as percentage of lesions per segment.
- 10 Qualitative morphologic assessment of lesion complexity is performed with light microscopy for the following parameters: absence or presence of foam cells, wall thickness, vessel wall architecture integrity, nuclear proliferation, and luminal extension.
- 11 Lesion morphology, complexity, and severity are used to classify the stage of lesions per the classification of Stary.
- 12 Lesion complexity can also be performed (e.g., for foam cells and macrophages)

Note: the above method is for quantitative microscopy. It is also possible to provide en face analysis of aortic atheroma by Sudan IV staining.