

May 14, 2019

UC Davis - Echocardiography

DOI

dx.doi.org/10.17504/protocols.io.yiafuae



Anil Singapuri¹

¹University of California, Davis

Mouse Metabolic Phenotyping Centers
Tech. support email: info@mmpc.org



Lili Liang

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.yiafuae>

External link: <https://mmpc.org/shared/document.aspx?id=115&docType=Protocol>

Protocol Citation: Anil Singapuri 2019. UC Davis - Echocardiography. **protocols.io**
<https://dx.doi.org/10.17504/protocols.io.yiafuae>



License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: February 25, 2019

Last Modified: May 14, 2019

Protocol Integer ID: 20770

Keywords: Cardiac hypertrophy, heart failure, Echocardiography, cardiac hypertrophy, development of cardiac hypertrophy, noninvasive imaging with echocardiogram, echocardiogram, common causes of heart failure, echocardiography summary, using noninvasive imaging, heart

Abstract

Summary:

Cardiac hypertrophy is one of the most common causes of heart failure. The development of cardiac hypertrophy and failure can be monitored using noninvasive imaging with echocardiogram in conscious animals.

Modified from: Timofeyev V. et al. J Mol Cell Cardiol. 2006 Jul;41(1):170-81.

Troubleshooting

- 1 The animals' chest will be shaved and then be placed in a supine position on a warmed (37°C) towel covered surgical table.
- 2 The mice are restrained in a 50 ml conical tube with an opening for the nose and chest, a standard procedure for imaging mice for this kind of study.
- 3 The imaging generally lasts for only 2-5 min, we were trained in this method by CLAS personnel and in our experience works very well.
- 4 M-mode and two-dimensional measurements are used to assess systolic function.
- 5 An average of six selected cardiac cycles from at least two separate scans performed in randomblindfashion with papillary muscles used as a point of reference.
- 6 End diastole was defined as the maximal left ventricular (LV) diastolic dimension and end systole was defined as the peak of posterior wall motion.
- 7 Fractional shortening (FS), a surrogate of systolic function, was calculated from LV dimensions as follows:
$$FS = ((EDD - ESD) / EDD) \times 100\%$$
where EDD and ESD are LV end diastolic and end systolic dimension, respectively.