

Aug 08, 2024

Version 2

URM C TriState SenNet Mouse Lung Digestion V.2

 Version 1 is forked from [Mouse Ischemia Experiment](#)

DOI

<https://dx.doi.org/10.17504/protocols.io.5qpvok9z9l4o/v2>

Gagandeep Kaur¹, Irfan Rahman²

¹Department of Environmental Medicine, University of Rochester Medical Center, Rochester, NY;

²University of Rochester

TriState SenNet



Gagandeep Kaur

University of Rochester Medical Center

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.5qpvok9z9l4o/v2>

Protocol Citation: Gagandeep Kaur, Irfan Rahman 2024. URM C TriState SenNet Mouse Lung Digestion. **protocols.io** <https://dx.doi.org/10.17504/protocols.io.5qpvok9z9l4o/v2> Version created by [Gagandeep Kaur](#)



License: This is an open access protocol distributed under the terms of the **[Creative Commons Attribution License](#)**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: August 08, 2024

Last Modified: August 08, 2024

Protocol Integer ID: 104916

Keywords: cellular senescence, lung tissue, Dissociation, urmc tristate sennet mouse lung digestion the objective, urmc tristate sennet mouse lung digestion, tristate sennet tmc bioanalyses core, tristate sennet tmc bioanalyses core at the university, scrna analysis, cellular senescence, part of the cellular senescence network program, mouse lung, cellular senescence network program, minimum damage to the cell

Funders Acknowledgements:

TriState SenNet

Grant ID: U54AG075931

Disclaimer

The authors have no conflict of interest to declare.

Abstract

The objective of this document is to share the material and kits used and steps involved in digesting the mouse lungs to perform scRNA analyses. This study was performed for the TriState SenNet TMC Bioanalyses Core at the University of Rochester, as part of the Cellular Senescence Network Program (SenNet). This protocol ensures minimum damage to the cells in suspension and ensures maximum viability to conduct downstream analyses.



Materials

Materials, Equipment & Reagents including Formulations

1. Liberase: (Cat# 5401127001; Roche)
2. DNase I (Cat# 4380-096-06; R&D)
3. Serum free DMEM

4. Disposables:
 - i) C-tubes for GentleMACS (Cat# 130-093-237, Miltenyi Biotec)
 - ii) MACS SmartStrainers 70µm (Cat# 130-098-462, Miltenyi Biotec)
 - iii) Sterile serological pipets; 50, 25,10, 5, 2, 1 mL size
 - iv) Sterile pipet tips; 1000, 200 and 20 µl
 - v) Sterile 15mL or 50mL conical polypropylene tubes
 - vi) Test tube racks for 15mL or 50mL conical tubes
 - vii) LD columns (Cat# 130-042-901; Miltenyi Biotec)

5. Equipment:
 - i) GentleMACS Tissue Octo Dissociator (Miltenyi Biotec Inc San Diego, CA)
 - ii) MACSmix tube rotator
 - iii) Table Top Centrifuge
 - iv) Biological Safety Cabinet, Class II
 - v) Dissecting forceps and scalpel



Lung Tissue Dissociation

- 1 Mince tissue finely, place into 50ml GentleMACS C-tube with liberase enzymatic cocktail .
- 1.1 For 50mg tissue use 0.5 ml liberase + 2ml DMEM + 2ul 5 unit/ul DNase I.
- 2 Immediately transfer the tubes to MACS Tissue Dissociator and run user-defined program named m_lung_01_02 for mouse lungs.
- 3 Thereafter, place 50ml conical with digestion into an incubating rocker for 30 minutes at  37 °C .

NOTE: If you still see chunks of lung tissue in the tubes after this step then transfer to the MACS Tissue Dissociator and run user-defined program named m_lung_02_01 for mouse lungs for 10-15 sec.
- 4 Remove conical from rocker and strain the cell suspension through 70 micron into 50mL tube
- 5 Add 3 ml DMEM (10% FBS) through the strainer to collect the remaining cell suspension
- 6 After straining, centrifuge at 300g for 5 minutes at  4 °C .

 
- 7 Remove supernatant, add 500 µl of red blood cell lysis to cell pellet. Incubate the cell suspension  On ice for 5 min.


- 8 After incubation add 4.5 mL of DMEM (10% FBS) to the suspension and centrifuge at 300g for 5 minutes at  4 °C .

 
- 9 Remove supernatant, re-suspend pellet in 2 mL DMEM (10%FBS)



10 Count cells and viability using AO/PI method, confirm reading under microscope