


Dec 26, 2023

## User Guide for CytoFLEX S Analyzer

DOI

<https://dx.doi.org/10.17504/protocols.io.yxmvm3br6l3p/v1>

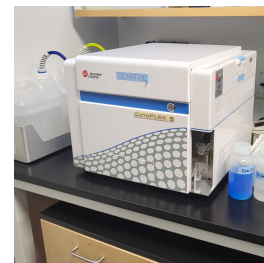

Jamie Tijerina<sup>1</sup>

<sup>1</sup>California Institute of Technology



Jamie C Tijerina

Universidade de Lisboa, California Institute of Technology

### Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



**DOI:** <https://dx.doi.org/10.17504/protocols.io.yxmvm3br6l3p/v1>

**Protocol Citation:** Jamie Tijerina 2023. User Guide for CytoFLEX S Analyzer. **protocols.io**  
<https://dx.doi.org/10.17504/protocols.io.yxmvm3br6l3p/v1>

**License:** This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited



**Protocol status:** Working

**We use this protocol and it's working**

**Created:** December 22, 2023

**Last Modified:** December 26, 2023

**Protocol Integer ID:** 92684

**Keywords:** caltech flow cytometry facility, instrument in both tube mode, instructions for troubleshooting, tube mode, basic instruction, user guide, instruction, guide

## Abstract

This is the Caltech Flow Cytometry Facility's User Guide for the CytoFLEX S Analyzer. This guide is given to all investigators who are trained to run the CytoFLEX S in the facility. A copy is kept in a binder next to the instrument at all times along with all other CytoFLEX S related protocols. This user guide includes basic instructions for operating the instrument in both tube mode and plate mode. It also contains instructions for troubleshooting.

## Before start

Make sure that you have checked the configuration of your instrument before purchasing reagents and preparing your samples. Before switching on the instrument, wipe down the area around the instrument with 70% ethanol or CaviWipes. Refill the sheath tank and empty the waste tank.



1

 Cytoflex User Guide Rev Dec 2022.... 849KB

Download the full user guide here.