

Mar 29, 2024

Validating Responsiveness of AAV-DIO-hM3Dq-DREADD

 [eLife](#)

DOI

dx.doi.org/10.17504/protocols.io.6qpvr3yrpvmk/v1

Katerina Rademacher¹, Ken Nakamura¹

¹Gladstone Institute of Neurological Disease



Haru Yamamoto

UCSF

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.6qpvr3yrpvmk/v1>

External link: <https://doi.org/10.7554/eLife.98775>

Protocol Citation: Katerina Rademacher, Ken Nakamura 2024. Validating Responsiveness of AAV-DIO-hM3Dq-DREADD. [protocols.io](https://dx.doi.org/10.17504/protocols.io.6qpvr3yrpvmk/v1) <https://dx.doi.org/10.17504/protocols.io.6qpvr3yrpvmk/v1>

Manuscript citation:

Rademacher K, Doric Z, Haddad D, Mamaligas A, Liao S, Creed R, Kano K, Chatterton Z, Fu Y, Garcia JH, Vance VM, Sei YJ, Kreitzer A, Halliday G, Nelson AB, Margolis E, Nakamura K Chronic hyperactivation of midbrain dopamine neurons causes preferential dopamine neuron degeneration. eLife 13(). doi: [10.7554/eLife.98775](https://doi.org/10.7554/eLife.98775)



License: This is an open access protocol distributed under the terms of the **[Creative Commons Attribution License](#)**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: January 07, 2024

Last Modified: May 31, 2024

Protocol Integer ID: 93034

Keywords: ASAPCRN, dopamine neurons with cno, activation of dreadd receptor, dreadd receptor, dopamine neuron, dreadd in datirescre mice, datirescre mice, responsiveness of aav, neuron, hm3dq, robust increase in locomotion, locomotion, dreadd, mice

Funders Acknowledgements:

ASAP

Grant ID: 020529

Abstract

To confirm responsiveness of AAV-DIO-hM3Dq-DREADD in DAT^{IRESCre} mice. Activation of DREADD receptors in dopamine neurons with CNO should induce a robust increase in locomotion.

Materials

- DAT^{IRESCre} mice bilaterally injected in the midbrain with AAV-DIO-hM3Dq-DREADD
- Med Associates wireless activity wheels (ENV-047)
- Med Associates Hub (DIG-807)
- Med Associates software (SOF-860)
- Clozapine-n-oxide (CNO; Tocris 4936-MTO)
- Sterile saline

Troubleshooting



Preparation

- 1 Mice should be at least two-weeks post-surgery.
- 2 Habituate mice to wireless running wheels in individually-housed cages.
 - 2.1 See Victoria Vance, Katerina Rademacher, Ken Nakamura 2023. Behavior Tracking with Running Wheels. protocols.io <https://protocols.io/view/behavior-tracking-with-running-wheels-cymrxu56>
- 3 Prepare CNO for IP injection.
 - 3.1 0.5mg/kg CNO dissolved in sterile saline, 500µL volume per mouse.
 - 3.2 Prepared fresh the day of injection.

Procedure

- 4 Note: All injections should be performed at the same time of day, at least 2–3 hours before the dark cycle.
- 5 Day 1: IP inject mice with 500µL of sterile saline.
- 6 Day 2: IP inject mice with 500µL CNO.
- 7 Day 3 or 4: IP inject mice again with 500µL CNO.
- 8 Responders increase their running wheel activity in the 10-60min following IP injection of CNO. Exclude mice that do not respond to IP CNO from further analysis.