


May 16, 2019

Vandy - Exercise Stress Test

DOI

<https://dx.doi.org/10.17504/protocols.io.yxcfxiw>

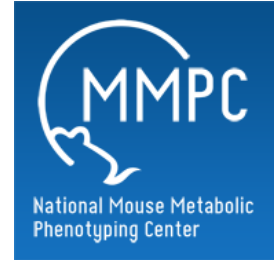

Louise Lantier¹

¹Vanderbilt University

Mouse Metabolic Phenotyping Centers
Tech. support email: info@mmpc.org



Lili Liang

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.yxcfxiw>

External link: <https://mmpc.org/shared/document.aspx?id=237&docType=Protocol>

Protocol Citation: Louise Lantier 2019. Vandy - Exercise Stress Test. **protocols.io**
<https://dx.doi.org/10.17504/protocols.io.yxcfxiw>



License: This is an open access protocol distributed under the terms of the [Creative Commons Attribution License](#), which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: March 07, 2019

Last Modified: May 16, 2019

Protocol Integer ID: 21188

Keywords: exercise test, mice, exercise stress test summary, exercise stress test, maximal exercise test, exhaustion as the treadmill speed, vo2 max, carbon dioxide output by the mouse, treadmill speed, maximum amount of oxygen, oxygen consumption, direct measurements of the oxygen consumption, stress, exhaustion, vandy, oxygen, test, maximal exercise

Abstract

Summary:

The exercise stress test is a maximal exercise test used to calculate VO2 max. Mice are run to complete exhaustion as the treadmill speed is increased every three minutes (See table 1). During the test, direct measurements of the oxygen consumption and carbon dioxide output by the mouse are made, which are then used to calculate VO2 max, VCO2, and RER. VO2 max refers to the maximum amount of oxygen that the mouse can utilize during maximal exercise.

Materials

MATERIALS

 Treadmill **Columbus Instruments**

 Oxymax **Columbus Instruments**

- 1 Mice are familiarized with the treadmill two days prior to the experiment by exercising them for 10 minutes at 10m/min.
- 2 On the day of the experiment, the mouse is placed in the stopped treadmill for 45 min to acclimatize and the Oxymax system is started. The treadmill is equipped with a shock pad. Shocks are set at 1.5mA, 200ms pulses, 4Hz.
- 3 After 45 min, basal measurements are made for 15 min in the stopped treadmill.
- 4 Treadmill is started at 10m/min.
- 5 After 3 min, speed is increased by 4m/min.
- 6 Every three minutes thereafter, speed is increased by 4m/min until mouse has reached exhaustion (table 1). Exhaustion is defined as the point at which the mouse remains on the shock pad for 5 continuous seconds.

Time (minutes)	Speed (m/min)
0	10
3	14
6	18
9	22
12	26
15	30
18	34
21	38
24	42
27	46

Table 1

- 7 Once exhaustion is reached, the mouse is left 30 min in the stopped treadmill to recover.