

May 16, 2019

Vandy - Mouse Echocardiography

DOI

dx.doi.org/10.17504/protocols.io.yyifxue



Chee Lim¹

¹Vanderbilt University

Mouse Metabolic Phenotyping Centers
Tech. support email: info@mmpc.org



Lili Liang

OPEN  ACCESS



DOI: dx.doi.org/10.17504/protocols.io.yyifxue

External link: <https://mmpc.org/shared/document.aspx?id=225&docType=Protocol>

Protocol Citation: Chee Lim 2019. Vandy - Mouse Echocardiography. **protocols.io**
<https://dx.doi.org/10.17504/protocols.io.yyifxue>

License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: March 09, 2019

Last Modified: May 16, 2019

Protocol Integer ID: 21226

Keywords: mouse echocardiography



Abstract

Summary:

Transthoracic mouse echocardiography is used to provide noninvasive imaging of the heart and allows for quantification of myocardial wall and chamber dimensions and systolic and diastolic performance.

Materials

MATERIALS



Isoflurane



Coupling gel

Vevo2100 Imaging System **VisualSonics Catalog #Vevo2100**

- 1 The chest hair is first shaved and a topical depilatory agent (e.g. Nair) is used to remove any remaining body hair.
- 2 The conscious mouse is held in the prone position, decreasing vagal reflexes and associated abnormalities of heartrate or AV conduction.
If anesthesia is required, 3% isoflurane is used for sedation and the mouse is placed in a supine position on a heated platform (to maintain body temperature) with embedded ECG leads (Visual Sonics). A nose cone is placed over the snout and isoflurane (1%) is delivered to maintain sedation throughout the procedure.
- 3 Ultrasound coupling gel heated to 34C is applied to the chest area and a linear array transducer (18-23 MHz) is positioned to obtain two-dimensional B-mode parasternal long and short axis views at the mid-ventricular level (Vevo 2100, VisualSonics).
- 4 One-dimensional M-mode images are obtained for measurement in the short axis view to obtain cardiac wall and chamber dimensions.
- 5 Digital images are permanently archived. Depending on the type of measurement, the imaging procedure can last from one to several minutes.
- 6 Left ventricular (LV) chamber size and wall thickness are measured off-line in the M-mode from at least three consecutive beats and averaged. LV wall thickness: intraventricular septum (IVS) and posterior wall (PW) at systole and diastole; and LV internal dimensions (LVID) during systole and diastole are measured. LV percent fractional shortening (FS) and ejection fraction (EF) are calculated from the M-mode measurements.