

Mar 19, 2025

Version 1

Viral extraction from Sediment V.1

DOI

<https://dx.doi.org/10.17504/protocols.io.6qpvr97zbvmk/v1>

Aditi K. Narayanan¹, Alon Philosof¹

¹California Institute of Technology

Aditi K. Narayanan: ORCID: 0000-0003-0627-1859

Alon Philosof: ORCID: 0000-0003-2684-8678

Orphan Lab



Victoria J Orphan

California Institute of Technology

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.6qpvr97zbvmk/v1>

Protocol Citation: Aditi K. Narayanan, Alon Philosof 2025. Viral extraction from Sediment. **protocols.io**
<https://dx.doi.org/10.17504/protocols.io.6qpvr97zbvmk/v1>

License: This is an open access protocol distributed under the terms of the **[Creative Commons Attribution License](#)**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: January 16, 2025

Last Modified: March 19, 2025

Protocol Integer ID: 118534

Keywords: viruses from sediment particle, sediment virus, viral extraction, detaching virus, sediment particle, subsurface sediment, extraction, density separation

Funders Acknowledgements:

NOMIS Foundation

U.S. Department of Energy

Grant ID: DE-SC0020373; DE-SC0022991

Disclaimer

DISCLAIMER – FOR INFORMATIONAL PURPOSES ONLY; USE AT YOUR OWN RISK

The protocol content here is for informational purposes only and does not constitute legal, medical, clinical, or safety advice, or otherwise; content added to [protocols.io](#) is not peer reviewed and may not have undergone a formal approval of any kind. Information presented in this protocol should not substitute for independent professional judgment, advice, diagnosis, or treatment. Any action you take or refrain from taking using or relying upon the information presented here is strictly at your own risk. You agree that neither the Company nor any of the authors, contributors, administrators, or anyone else associated with [protocols.io](#), can be held responsible for your use of the information contained in or linked to this protocol or any of our Sites/Apps and Services.

Abstract

Viruses are often undercounted in subsurface sediments due to their propensity to stick to surfaces. This protocol outlines a method for detaching viruses from sediment particles via a combination of chemical and physical agitation followed by density separation.

Aditi K. Narayanan and Alon Philosof

- 1 Based on *An Improved Method for Extracting Viruses From Sediment: Detection of Far More Viruses in the Subseafloor Than Previously Reported*

<https://doi.org/10.3389/fmicb.2019.00878> (Pan et al 2019, Frontiers in Microbiology)

Materials

- 2
 1. 20mM tetrasodium pyrophosphate solution (0.532g in water for 100mL final volume), CAS #7722-88-5, Millipore-Sigma Cat. #P8010
 2. 10% NaCl solution (10g in water for 100mL final volume), CAS #7647-14-5
 3. 30% Nycodenz solution (4.5g in water for 15mL final volume). Nycodenz is also called Histodenz or Iohexol), CAS #66108-95-0, Millipore-Sigma Cat. #D2158
 4. 50% Nycodenz solution (7.5g in water for 15mL final volume)
 5. 15mL Falcon tubes or similar (e.g. Becton Dickinson Cat. #4-959-49B), capable of withstanding ~4000xg
 6. Sonication wand (e.g. QSonica CL-188)
 7. Centrifuge capable of holding 15mL Falcon Tubes (e.g. Beckman Coulter Allegra X-15R Centrifuge with SX4570 swinging bucket rotor)
 8. Syringes, various volumes
 9. Hypodermic needles, between preferably 3-4 inches in length if possible (otherwise 1.5 inches will do)
 10. Glutaraldehyde solution (optional, for fixation)
 11. Liquid nitrogen (optional, for storage after fixation)

Protocol

- 3 Take 0.1-0.33cm³ of sample (or 1mL of slurry) and mix enough NaCl solution and tetrasodium pyrophosphate solution to get to a final concentration of 2.5% NaCl and 5mM tetrasodium pyrophosphate in a 6mL volume. Your stock solutions under the Materials section should be 4X concentrations. Get rid of any clumps by shaking. Works best in a 15mL Falcon tube.

For example, to each 1mL of slurry, add 1.5mL of 4X tetrasodium pyrophosphate, 1.5mL of 4X NaCl, and 2mL of virus-free water.

- 4 Sonicate on ice at 25 amps for 10 seconds. Rest for another 10 seconds. Repeat 2 more times for a total of 30 seconds of sonication. Our sonicator is a QSonica CL-188.
- 5 Create Nycodenz 30% and 50% gradient in a 15mL falcon tube using the underlay method:



1. Fill a syringe with 1mL of 30% solution. Attach a needle to the end.
 2. Tap out any air bubbles in the syringe
 3. Dispense the 30% solution into the very bottom of the Falcon tube
 4. Repeat the above steps 1-2 with the 50% solution
 5. Push the needle of the syringe containing the 50% solution all the way to the bottom of the Falcon tube
 6. CAREFULLY dispense the 50% solution under the 30% solution, allowing the 30% layer to rise above it
- 6 Layer the sonicated sample over the gradient and centrifuge at 2900xg for 30min in a swinging bucket rotor.
- 7 Use a needle and syringe to remove the topwater, 30% fraction, and liquid portion of the 50% fraction. These fractions will contain your viruses, though you will need to stain with SYBR Gold to determine which of your fractions carries the largest number.
- 8 Filter the viral fraction through a 0.2µm membrane and store. If you choose to fix before storing, mix enough glutaraldehyde with the sample for a final concentration of 0.1% glutaraldehyde. Allow the samples to sit at 4°C for 30 minutes, then flash freeze and store at -80°C.
- 9 If you choose, you can repeat the pyrophosphate, sonication, and gradient steps on the sediment pellet to ensure maximum extraction efficiency.
1. Add 5-6mL of 1:1 solution (final concentration 5mM tetrasodium pyrophosphate and 2.5% NaCl)
 2. Repeat steps 2-6

Protocol references

<https://doi.org/10.3389/fmicb.2019.00878> (Pan et al 2019, Frontiers in Microbiology)