

Jul 12, 2023

Water Sampling and Plating Generations (Antimicrobial Group)

DOI


<https://dx.doi.org/10.17504/protocols.io.ewov1yd87vr2/v1>

Melissa B Duhaime¹, Aubrey Reed¹, Nicole Sunshine¹, Renae Lyons¹

¹University of Michigan



Nicole Sunshine

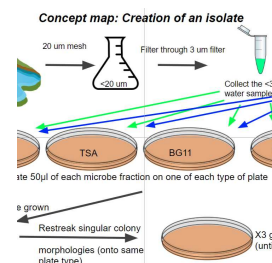

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.ewov1yd87vr2/v1>

Protocol Citation: Melissa B Duhaime, Aubrey Reed, Nicole Sunshine, Renae Lyons 2023. Water Sampling and Plating Generations (Antimicrobial Group). **protocols.io** <https://dx.doi.org/10.17504/protocols.io.ewov1yd87vr2/v1>

License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: December 09, 2018

Last Modified: July 12, 2023

Protocol Integer ID: 18493

Keywords: microbes from freshwater lake, advanced environmental microbiology lab, antimicrobial group class project, microbiology, environmental reservoirs of antibiotic resistance, antimicrobial group, microbe, antibiotic production in nature, microbes in the wild, antibiotic production, antibiotic resistance, antimicrobial compound, freshwater lake, antimicrobial compounds in order, michigan biological station, environmental reservoir, protocol for class project, plating generation, sampling, class project

Abstract

Protocol developed by the Antimicrobial Group class project. The goal was to isolate microbes from freshwater lakes that produce antimicrobial compounds in order to study antibiotic production in nature and environmental reservoirs of antibiotic resistance.

Protocol for class project in **Microbes in the Wild: Advanced Environmental Microbiology Lab (EEB 447)** course at University of Michigan Biological Station and on campus.

- 1 Collect whole water at six sites using a Van Dorn water collector for deep (~122.5-123.7 m) samples and 2 L containers for surface samples (~15 cm below the surface).
- 2 Set site names and classifications

Lake	Location Type	Location Name	Latitude	Longitude	Depth
Lake Michigan	Pristine	Off Shore A) Shallow B) Deep	45.28287	-85.25048	A) 15cm B) 123.7M
Lake Michigan	Pristine	Near Shore A) Shallow B) Deep	45.21355	-85.16976	A) 15cm B) 122.5M
Lake Michigan	Disturbed	Clad 2	45.2778	-85.135415	15cm
Douglas Lake	Disturbed	Septic	45.555901	-84.665300	15cm

Table 1: Sampling site specifics. We took samples from both Lake Michigan and Douglas Lake (inland). We took note of whether the site was anthropologically disturbed or pristine, gave it a specific location title, noted latitude and longitude of the site, and noted the depth at which the sample was taken.

Sites samples and description


3. Prefilter whole water using 20µm mesh.
 4. Set up peristaltic pump with 3 µm filter
 5. Filter approximately 500mL of sample water.
 6. Collect the filter in an eppendorf with 1mL of PBS as the particle-associated large fragment
- 3 Collect the <3µm filtered water as the free-living (FL) smaller fraction.
 - 4 For the particle-associated (PA) fraction, vortex tube, then partially remove filter with an edge under the closed cap
 - 5 Centrifuge at 1500 RPM for 1 minute
 - 6 Fully remove filter and resuspend pellet

- 7 Pipette 10µL of both PA and FL onto separate plates of 4 agar types (TSA, LB, BG11, and LAKE) and spread.

This is the parent generation.

- 8 PROPAGATE NEXT GENERATIONS:

- 8.1 From the lawns initially plated, by sterile loop pick up a single colony.
- 8.2 Select a plate of the same media.
- 8.3 Streak for isolation using a three phase streak pattern as shown below, sterilizing the loop in between each quadrant.

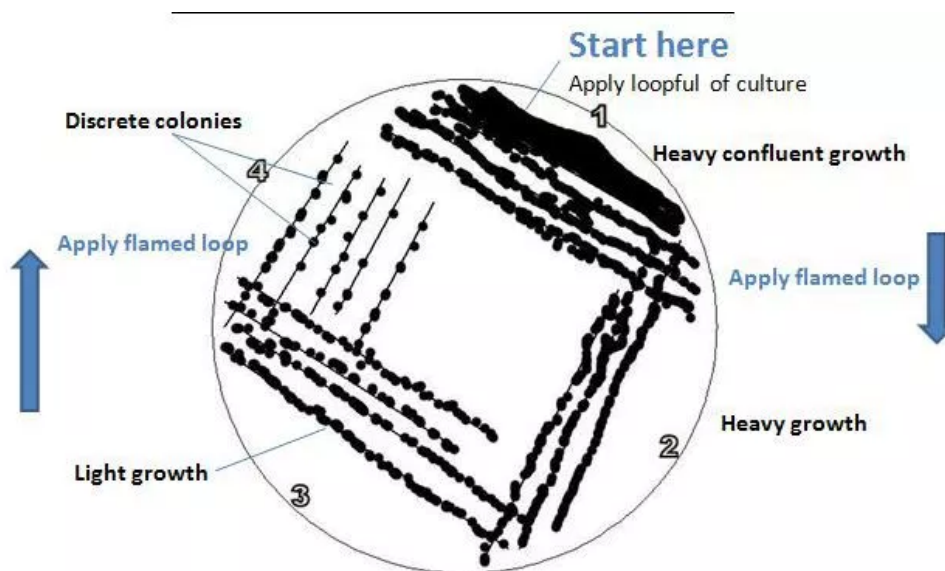

4. Label these plates "generation 1"
5. Continue until generation 3 is obtained.
6. Use generation 3 for testing.

