



Apr 03, 2023

Western Blot Analysis

 In 1 collection

DOI

<https://dx.doi.org/10.17504/protocols.io.j8nlkwp515r/v1>

michela.deleidi¹, Federico Bertoli¹

¹German Center for Neurodegenerative Diseases (DZNE), Tübingen, 72076 Germany



Federico Bertoli

ASAP

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.j8nlkwp515r/v1>

Protocol Citation: michela.deleidi , Federico Bertoli 2023. Western Blot Analysis. **protocols.io**

<https://dx.doi.org/10.17504/protocols.io.j8nlkwp515r/v1>

License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working



Created: April 01, 2023

Last Modified: May 31, 2024

Protocol Integer ID: 79873

Keywords: ASAPCRN, western blot analysis, western blot, protocol

Abstract

Western blot protocol

- 1 Wash cells 1X in PBS.
- 2 Detach cells using Accutase for 5 minutes at 37°C and collect them.
- 3 Spin cells in a centrifuge at 250g for 5 minutes at room temperature.
- 4 Remove the supernatant.
- 5 Lyse cells in 1% TBS + 0.5% NP40 + PI/PHI (Pierce #A32959).
- 6 Determine protein concentration was determined using the Pierce BCA Protein Assay Kit (Thermo Fisher Scientific).
- 7 Dilute samples in 6x Laemmli buffer containing 12.5% β -mercaptoethanol.
- 8 Boil samples 5 minutes at 95°C in a thermoblocker.
- 9 Load protein in a self-casted 7.5-15% acrylamide gel or in precast NuPage 4-12% Bis-Tris Protein gel.
- 10 Run gel at 100V for 30min-2h.
- 11 Transfer proteins to PDVF membrane via wet transfer (20% methanol) overnight at 20V at 4°C.
- 12 Block membrane in TBS 0.1% Tween containing 5% milk or 5% bovine serum albumin.



- 13 Incubate primary or secondary antibodies in 5% Roche Block solution or 5% BSA in TBS-T supplemented with 0.04% sodium azide.