

Nov 10, 2022

Western Blot

DOI

dx.doi.org/10.17504/protocols.io.14egn22emg5d/v1

Addgene The Nonprofit Plasmid Repository¹

¹Addgene



Addgene The Nonprofit Plasmid Repository

Addgene

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.14egn22emg5d/v1>

Protocol Citation: Addgene The Nonprofit Plasmid Repository 2022. Western Blot. **protocols.io**
<https://dx.doi.org/10.17504/protocols.io.14egn22emg5d/v1>

License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: November 04, 2022



Last Modified: November 10, 2022

Protocol Integer ID: 72299

Keywords: western blot western blot, western blot, total protein concentration in the lysate, lysing cell, basic steps for lysing cell, antibodies specific to the protein, lysate, using antibody, antibody, determining total protein concentration, proteins by size, protein

Disclaimer

We have listed the specific equipment, reagents, and methods that we use in our lab at Addgene. Equipment and reagents from other vendors should produce similar outcomes when using these protocols. However, please be aware that your protocol may need to be adjusted to accommodate slight differences between products. Addgene does not endorse or recommend specific products or equipment. Inclusion of this information is solely for transparency intended to support reproducibility in science.

Abstract

Western blot is a technique used to separate proteins by size followed by detection using antibodies specific to the protein of interest. This protocol describes the basic steps for lysing cells, determining total protein concentration in the lysate, running a precast SDS-PAGE gel, and immunoblotting.



Guidelines

General Considerations

Acrylamide percentage varies in SDS-PAGE gels. Use lower percentage acrylamide when immunoblotting high molecular weight proteins and higher percentage acrylamide when immunoblotting low molecular weight proteins.

This protocol uses a dry transfer device but can be adapted for wet and semi-dry transfer approaches.

Workflow Timeline

Day 1: Prepare lysates, run SDS-PAGE, transfer, block, incubate with primary antibody

Day 2: Incubate with secondary antibody

Tips and Troubleshooting:

The optimal lysis buffer will vary depending on the sample type and cellular location of the protein of interest. You may need to try a variety of lysis buffers to find the best for your target.

The optimal concentration and blocking buffer will vary between antibodies. Review the instructions before starting your experiment and consider titrating your antibody to determine the optimal dose.

To ensure that your antibody is both functioning as expected and is specific, include a positive control sample that you know expresses the protein, such as cells transfected with the protein of interest, and a negative control sample that does not express the protein of interest.

Materials

Equipment:

- Microcentrifuge
- 0.5–10 µL single channel pipette
- 2–20 µL single channel pipette
- 20–200 µL single channel pipette
- 200–1000 µL single channel pipette
- Pipette controller
- Pipette tips and pipettes
- Spectrophotometer
- Heat block
- Mini gel tank chamber
- Power supply
- iBlot 2 Gel Transfer Device
- Roller
- Spatula
- Platform shaker
- Cold room
- Gel imager
- -80 °C freezer

Reagents and Consumables:

- 1X PBS
- Lysis buffer e.g., RIPA lysis buffer
- Microcentrifuge tubes
- BCA Assay, Thermo Fisher 23227
- β-mercaptoethanol
- 4X protein loading buffer
- Precast SDS-PAGE gel
- SDS-PAGE running buffer
- Prestained protein ladder
- Ethanol
- iBlot 2 PVDF Mini Stack, Thermo Fisher IB24002
- 20X TBS
- Tween-20
- Nonfat milk powder
- 96-well microtiter plate
- Chemiluminescence substrate
- Plastic wrap
- Primary antibody



- Secondary antibody
- Deionized water

Troubleshooting







Before start

Refer to the manufacturer's instructions for additional information specific to your antibody, such as ideal blocking buffer and optimal antibody concentrations. Consider titrating your antibody to determine the optimal dose.

Secondary antibodies must match the host species of the primary antibody. For example, use an anti-mouse secondary antibody for primary antibodies raised in a mouse.







Lyse cells


- 1 Centrifuge  0 μL for  00:05:00 at  100 x g . 5m
- 2 Gently remove supernatant.
- 3 Resuspend cell pellet in  1 mL of 1X PBS and transfer to a microcentrifuge tube.
- 4 Centrifuge for  00:05:00 at  100 x g . 5m
- 5 Carefully remove supernatant.
- 6 Resuspend cell pellet in appropriate volume of cold lysis buffer.

Note

The volume of lysis buffer will vary depending on the size of your cell pellet but will generally be between 250–1000 μL .

Note

The ideal lysis buffer will vary depending on the cellular location of the protein of interest. RIPA buffer is suitable for most proteins but more stringent buffers and a sonication step may be required for hard to extract proteins such as those in the nucleus.
- 7 Incubate on ice for  00:30:00 . 30m
- 8 Centrifuge lysate for  00:15:00 at  14.000 x g at  4 $^{\circ}\text{C}$. 15m
- 9 Transfer supernatant to a clean microcentrifuge tube.



10 Use the lysate immediately or store at  -80 °C until ready to use.

Determine the total protein concentration and prepare the lysate for SDS-PAGE

11 Determine the protein concentration using a Pierce BCA Assay Kit or other preferred method for protein determination.

11.1 Prepare a 50:1 Reagent A to Reagent B dilution of the BCA assay.

11.2 Prepare serial dilutions of the BSA standard that range from 0–2000 µg/mL.

11.3 In duplicate, dilute  10 µL of standard, blank, and lysate samples into  200 µL of BCA reagent in a 96-well microtiter plate.

11.4 Incubate for  00:30:00 at  37 °C .

30m

11.5 Determine the absorbance at 590 nm.

11.6 Calculate the average absorbance of the duplicate samples on the plate.

11.7 Subtract the average absorbance of the blank from all of the samples.

11.8 Plot a standard curve of BSA standard concentration versus absorbance.

11.9 Extrapolate the total protein concentration of the sample from the standard curve.

12 Determine the volume of sample required to load an equivalent amount of total protein for each sample.

**Note**

The ideal total protein loaded will vary between samples and target proteins but is typically between 10–50 µg. If the protein is in low abundance in the sample you will need to load a greater amount of total protein.

13 Prepare sample for loading as follows:

13.1 Add 10% v/v β-mercaptoethanol to the 4X protein loading buffer.

13.2 Dilute 4X protein loading buffer in the sample to 1X.

14 Boil the samples for  00:10:00 at  100 °C .

10m

SDS-PAGE

15 Prepare the precast gel as follows:

15.1 Remove the gel from the plastic packaging.

15.2 Remove tape and plastic comb.



15.3 Rinse the wells with deionized water 3x.

15.4 Shake the gel gently between washes to remove residual water.

15.5 Load the gel into the chamber of the SDS-PAGE gel tank.

**Note**



Different SDS-PAGE gel tanks will have different gel orientations. Refer to the instructions for your specific tank.

- 15.6 Close the clamp.
- 16 Prepare 1X running buffer as follows:
- 16.1 Dilute  25 mL of 20X running buffer to  500 mL with deionized water. Mix well.
- 17 Fill the chamber of the SDS-PAGE gel tank with the 1X running buffer.
- 18 Load the samples into the gel.
- 19 Load 5–10 μL of the prestained protein ladder.
- 20 Place the lid on the tank and plug the electrode cords into the power supply.
- 21 Run the gel at **100 V** for 10–15 min or until the samples have moved out of the wells and into the gel.
- 22 Increase the voltage to **150 V** and continue running the gel until you have obtained the desired separation.
- 23 Gently open the gel with the spatula.

Dry transfer

- 24 Soak the gel for  00:15:00 in 20% ethanol in deionized water.

15m

- 24.1 To prepare 20% ethanol, dilute  2 mL of ethanol into  8 mL of deionized water and mix well.
- 25 Assemble the transfer sandwich as follows:
- 25.1 Unseal the iBlot 2 PVDF Mini transfer stack.
- 25.2 Set the Top Stack to one side and discard the white separator.
- 25.3 Keep the Bottom Stack in the plastic tray.
- 25.4 Place the Bottom Stack on the blotting surface.
- 25.5 Align the electrical contacts on the blotting surface of the iBlot 2 Gel Transfer Device.
- 25.6 Wet the pre-run gel in deionized water and place it on the transfer membrane of the bottom stack.
- 25.7 Soak a piece of iBlot Filter Paper in deionized water.
- 25.8 Place the pre-soaked iBlot Filter Paper on the gel and remove the air bubbles with the roller.
- 25.9 Place the Top Stack over the pre-soaked filter paper.
- 25.10 Remove air bubbles using the roller.
- 25.11 Place the Absorbent Pad on top of the Top Stack such that the electrical contacts are aligned with the corresponding electrical contacts on the blotting surface of the iBlot 2



Gel Transfer Device.

25.12 Close the lid of the device.

25.13 Select the desired method and make sure the parameters are correct.

25.14 Transfer for 5–6 min for proteins <30 kDa.


25.15 Transfer for 8–10 min for proteins >150 kDa.


25.16 Select Start Run.


26 When the run is complete, select Done.

Blocking

27 Prepare 1X TBST as follows:

27.1  25 mL of 20X TBS



27.2  2.5 mL of Tween-20

27.3  472.5 mL of deionized water

27.4 Mix well

28 Prepare blocking buffer as follows:





- 28.1 Dilute 5% w/v non-fat milk into 1X TBST. (Example:  5 g of non-fat milk into  100 mL of 1X TBST.)



Note

The ideal blocking buffer will vary between antibodies. Refer to the manufacturer's information before running your experiment.

- 29 Remove the membrane from the transfer device and place it in a container with 1X TBST, protein side up.

- 30 Block the membrane in blocking buffer for  01:00:00 at  Room temperature on a shaking platform.

1h

- 31 Wash the membrane 3x for  00:05:00 in 1X TBST at  Room temperature on a shaking platform.

5m

Antibody incubation

1h 15m



- 32 Dilute the primary antibody to the desired concentration in blocking buffer.

Note



The optimal concentration will vary between antibodies but is typically between 1–10 µg/mL.

- 33 Incubate the membrane overnight in primary antibody at  4 °C on a rocking platform.

Note

Primary antibody incubation can be reduced to  02:00:00 at  Room temperature but may result in more non-specific binding.







34 Wash the membrane 3x for  00:05:00 in 1X TBST at  Room temperature on a shaking platform. 5m

35 Dilute the horseradish peroxidase-conjugated secondary antibody to the desired concentration in blocking buffer.



Note

The optimal concentration will vary between antibodies but is typically between 1–10 µg/mL.

36 Incubate the membrane with secondary antibody for  01:00:00 at  Room temperature on a shaking platform. 1h

37 Wash the membrane 3x for  00:05:00 in 1X TBST at  Room temperature on a shaking platform. 5m

38 Prepare the chemiluminescence substrate by mixing 1:1 reagent A to reagent B.

39 Gently incubate the membrane in the chemiluminescence reagent for  00:05:00 at  Room temperature . 5m

40 Cover the membrane with clear plastic and use a gel imager with chemiluminescence detection or a dark room to detect the bands.